

Agilent 325 Dual Wavelength UV-Vis Detector



Features, Technical Details, Specifications, and Ordering Details

The Agilent 325 Dual Wavelength UV-Vis Detector delivers superb sensitivity for detection in the UV and Visible light ranges. Dual path-length flow cell technology offers a dynamic absorbance range 40 times greater than conventional detectors, especially for preparative LC applications.

Features


- Secure compound identification Simultaneously handling two user-defined UV or visible wavelengths increases analyte information and enhances productivity.
- **Superb dynamic range** Dual path-length flow cell delivers a dynamic absorbance range 40 times greater than conventional detectors.
- Scale-up easily from analytical to preparative flow rates. Interchangeable, cartridge-style flow cells quickly switch flow rates from one microliter to hundreds of milliliters per minute.
- Agilent OpenLAB CDS ChemStation Edition Software or Agilent LC Responder Software provide full system control, data analysis and multiple reporting options plus tools for 21 CFR Part 11 compliance.
- Stable, precise wavelength settings Deuterium source is used for the ultraviolet range and a quartz halogen lamp for the visible range.
- **Tool-free maintenance** Routine maintenance is easy with front panel access. Simple thumbscrews secure the lamps and flow cells, no tools are needed.
- **Regulatory compliance** Agilent offers complete IQ, QQ, and PQ documentation packages and contract services to help meet regulatory requirements.



Technical Details

Dual pathlength flow cell


The dual path-length flow cell in the 325 Dual Wavelength UV-Vis Detector handles both low to high flow rates and low to high absorbance. The long path length measures low absorbance readings (<0.8 AUFS) and the shorter path-length measures high absorbance readings (>1.2 AUFS) combining to deliver a dynamic absorbance range from from 0.0005 AU up to 80 AU up to 40 times more than conventional detectors.

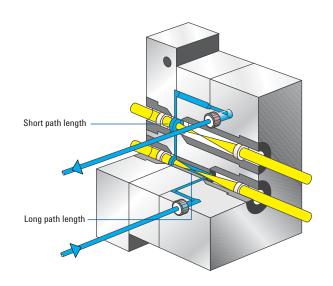

Achieve up to 80 AU when moving from analytical to preparative operation, without changing the dual path-length flow cell. Above shows dynamic range from 2 to 21 AU of 5-hydroxytryptophan sample.

Column: Agilent Pursuit XRs C18
Mobile phase: 85% water, 15% methanol
Detection: UV at 230 and 280 nm

Flow: 21 mL/min

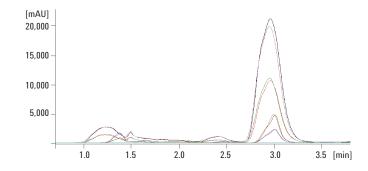

Specifications

Wavelength range	190 to 900 nm	
Light sources	Deuterium for UV and quartz halogen for visible range	
Spectral bandwidth	6 nm	
Wavelength precision	±0.1 nm	
Wavelength accuracy	±1 nm at 260 nm	
Noise, single wavelength	$\pm 5 \times 10^{-6}$ AU peak-to-peak at 254 nm, 2 seconds response time, 1 mL/min MeOH, 9-mm flow cell	
Noise, dual wavelength	$\pm 10 \times 10^{-6}$ AU peak-to-peak at 254/280 nm, 2 seconds response time, 1 mL/min MeOH, 9-mm flow cell	
Noise, dry cell	Typically, $\pm 3 \times 10^{-6}$ AU peak-to-peak at 254 nm, 2 seconds response time, 9 mm flow cell	
Drift:	Less than 1 mAU/hr	
Linearity, with acetone in water	Analytical 1% to 2 AU; preparative1% to 20 AU at 265 nm	
Absorbance range	Up to 80 AU with preparative flow cell	
Response time, selectable	0.05, 0.5, 1, and 2 seconds	
Data rate	20 Hz single channel	
Analog outputs	Two with selectable absorbance range of 0.001 to 100 AU full scale at 1 V	
External	Ethernet	
Contact closure outputs	Four time-programmable external event relays, one peak sense relay, three synchronization signals (ready out, start out, and fault out)	
Contact closure inputs	Lamp off, autozero, three synchronization signals (ready in, start in, and fault in)	
Software	Agilent OpenLab ChemStation	
Dual channel ratio calculation	Peak detection through OpenLab ChemStation	
Dual pathlength flow cells	Preparative flow cell: $9 \text{ mm} \times 1 \text{ mm}$; 4×0.25 Super preparative flow cell: $4 \text{ mm} \times 0.15 \text{ mm}$	
Maximum flow rate preparative flow cell	Preparative flow cell: 500 mL/min Super preparative flow cell: 10 L/min	
Pressure	1,000 psi maximum	

Ordering Details

Agilent 325 Dual Wavelength UV-Vis Detector

Description	
Agilent 325 UV Dual Wavelength UV-Vis Detector (190–900 nm)	
PWB Synchronization Interface 325/335	G9309A #183

Note: You must order the appropriate flow cell separately.

Flow Cells

Description	Part number
Flow cell 9 × 1 mm, dual path-length (Prep)	G9309A #180
Flow cell 4 × 0.25 mm dual path-length (Semi-prep)	G9309A #181
Flow cell 4 × 0.15 mm dual path-length (Super-prep)	G9309A #182

www.agilent.com/chem/ purification

© Agilent Technologies, Inc., 2013 Published in the USA, January 1, 2013 5991-1639EN

