




## Using the CapLok™ Tool to Undo and Tighten 1/4 inch Brass Long-Term Storage Caps

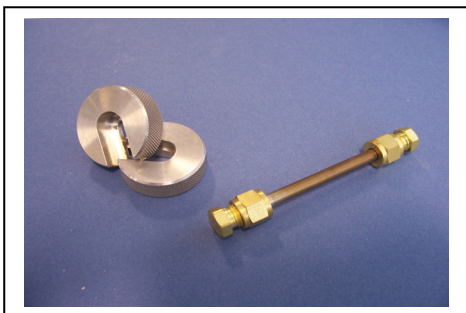

1. CapLok tool and sorbent tube capped with 1/4 inch brass SwageLok type long-term

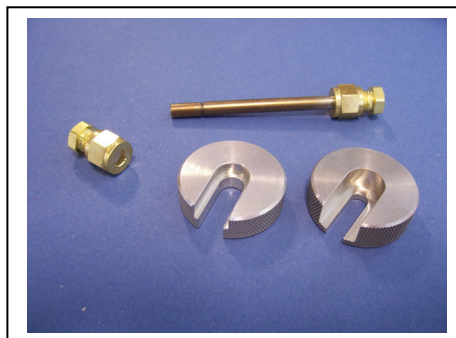

2. The two halves of the CapLok tool are slightly different. The disk etched with the Markes Logo should be used on the top of the cap and the other disk should be used over the tube on the nut end

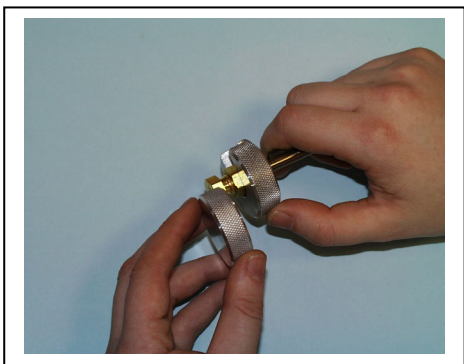

3. Loosen the brass cap using the CapLok tool as shown (by twisting the two discs in opposing directions). Then pull the nuts from the end of the tubes.



4. Remove the caps from the end of the tubes



5. To replace the caps, push them onto the end of the tube and tighten until finger tight. Use the CapLok tool (by twisting the two discs in opposite directions) to tighten the nut a further quarter to half turn.



6, Gently pull the brass caps at each end to ensure they are secure.