



Agilent 1260 Infinity Bio-inert Quaternary Pump

User Manual



Agilent Technologies

Notices

© Agilent Technologies, Inc. 2011-2012

No part of this manual may be reproduced in any form or by any means (including electronic storage and retrieval or translation into a foreign language) without prior agreement and written consent from Agilent Technologies, Inc. as governed by United States and international copyright laws.

Manual Part Number

G5611-90000 Rev. C

Edition

05/12

Printed in Germany

Agilent Technologies
Hewlett-Packard-Strasse 8
76337 Waldbronn

This product may be used as a component of an in vitro diagnostic system if the system is registered with the appropriate authorities and complies with the relevant regulations. Otherwise, it is intended only for general laboratory use.

Warranty

The material contained in this document is provided “as is,” and is subject to being changed, without notice, in future editions. Further, to the maximum extent permitted by applicable law, Agilent disclaims all warranties, either express or implied, with regard to this manual and any information contained herein, including but not limited to the implied warranties of merchantability and fitness for a particular purpose. Agilent shall not be liable for errors or for incidental or consequential damages in connection with the furnishing, use, or performance of this document or of any information contained herein. Should Agilent and the user have a separate written agreement with warranty terms covering the material in this document that conflict with these terms, the warranty terms in the separate agreement shall control.

Technology Licenses

The hardware and/or software described in this document are furnished under a license and may be used or copied only in accordance with the terms of such license.

Restricted Rights Legend

If software is for use in the performance of a U.S. Government prime contract or subcontract, Software is delivered and licensed as “Commercial computer software” as defined in DFAR 252.227-7014 (June 1995), or as a “commercial item” as defined in FAR 2.101(a) or as “Restricted computer software” as defined in FAR 52.227-19 (June 1987) or any equivalent agency regulation or contract clause. Use, duplication or disclosure of Software is subject to Agilent Technologies’ standard commercial license terms, and non-DOD Departments and Agencies of the U.S. Government will

receive no greater than Restricted Rights as defined in FAR 52.227-19(c)(1-2) (June 1987). U.S. Government users will receive no greater than Limited Rights as defined in FAR 52.227-14 (June 1987) or DFAR 252.227-7015 (b)(2) (November 1995), as applicable in any technical data.

Safety Notices

CAUTION

A **CAUTION** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a **CAUTION** notice until the indicated conditions are fully understood and met.

WARNING

A **WARNING** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a **WARNING** notice until the indicated conditions are fully understood and met.

In This Guide...

This manual covers the Agilent 1260 Infinity Bio-inert Quaternary Pump G5611A.

1 Introduction

This chapter gives an introduction to the module, instrument overview and internal connectors

2 Site Requirements and Specifications

This chapter provides information on environmental requirements, physical and performance specifications.

3 Installing the Pump

This chapter gives information about the preferred stack setup for your system and the installation of your module.

4 Using the Pump

This chapter explains the operational parameters of the module.

5 Optimizing Performance

This chapter gives hints on how to optimize the performance or use additional devices.

6 Troubleshooting and Diagnostics

This chapter gives an overview about the troubleshooting and diagnostic features and the different user interfaces.

7 Error Information

This chapter describes the meaning of error messages, and provides information on probable causes and suggested actions how to recover from error conditions.

In This Guide...

8 Test Functions and Calibration

This chapter describes the tests for the module.

9 Maintenance

This chapter describes the maintenance of the module.

10 Parts for Maintenance

This chapter provides information on parts for maintenance.

11 Identifying Cables

This chapter provides information on cables used with the Agilent 1200 Infinity Series modules.

12 Hardware Information

This chapter describes the pump in more detail on hardware and electronics.

13 Appendix

This chapter provides addition information on safety, legal and web.

Contents

1	Introduction	9
	Introduction to the Pump	10
	Overview of the Hydraulic Path	11
	Bio-inert Materials	17
	Early Maintenance Feedback	19
	Instrument Layout	20
2	Site Requirements and Specifications	21
	Site Requirements	22
	Physical Specifications	25
	Performance Specifications	26
3	Installing the Pump	27
	Unpacking the Pump	28
	Optimizing the Stack Configuration	30
	Installing the Pump	33
	Connecting Modules and Control Software	36
	Flow Connections of the Pump	38
	Priming the System	41
4	Using the Pump	47
	Hints for Successful Use of the Quaternary Pump	48
	Prevent Blocking of Solvent Filters	49
	Algae Growth in HPLC Systems	50
	Solvent Information	51
5	Optimizing Performance	55
	When to Use a Degasser	56
	Operational Hints for the Multi Channel Gradient Valve (MCGV)	57
	When to use the Seal Wash Option	58
	Choosing the Right Pump Seals	59
	Optimize the Compressibility Compensation Setting	60

6	Troubleshooting and Diagnostics	63
	Overview of the Module's Indicators and Test Functions	64
	Status Indicators	66
	User Interfaces	68
	Agilent Lab Advisor Software	69
7	Error Information	71
	What Are Error Messages	73
	General Error Messages	74
	Module Error Messages	80
8	Test Functions and Calibration	95
	Introduction	96
	System Pressure Test	97
	Leak Rate Test	102
9	Maintenance	107
	Introduction to Maintenance	108
	Warnings and Cautions	109
	Overview of Maintenance	111
	Cleaning the Module	112
	Checking and Replacing the Solvent Filter	113
	Exchanging the Active Inlet Valve (AIV) or its Cartridge	114
	Exchanging the Outlet Valve	117
	Exchanging the Purge Valve Frit or the Purge Valve	119
	Removing the Pump Head Assembly	122
	Seal Wear-in Procedure	125
	Maintenance of the Pump Head	126
	Reinstalling the Pump Head Assembly	129
	Exchanging the Multi-Channel Gradient Valve (MCGV)	131
	Exchanging the Optional Interface Board	134
	Replacing the Module's Firmware	136

10 Parts for Maintenance 137

Pump Head Assembly with Seal Wash Option	138
Outlet Valve	140
Purge Valve Assembly	141
Active Inlet Valve	142
Accessory Kit, Bio-inert (G5611-68755)	144
Starter-Kit Bio-inert	145
Solvent Cabinet	146
Bottle Head Assembly	148
Hydraulic Path	149
System Tool Kit	151

11 Identifying Cables 153

Cable Overview	154
Analog Cables	156
Remote Cables	158
BCD Cables	161
CAN/LAN Cables	163
External Contact Cable	164
Agilent Module to PC	165
Agilent 1200 Module to Printer	166

12 Hardware Information 167

Electrical Connections	168
Interfaces	170
Setting the 8-bit Configuration Switch (without On-board) LAN	176

13 Appendix 181

General Safety Information	182
The Waste Electrical and Electronic Equipment Directive	185
Batteries Information	186
Radio Interference	187
Sound Emission	188
Agilent Technologies on Internet	189

Contents



1 Introduction

Introduction to the Pump	10
Overview of the Hydraulic Path	11
Hydraulic Path	12
How Does the Pump Work?	13
How Does Compressibility Compensation Work?	15
How Does Variable Stroke Volume Work?	16
Bio-inert Materials	17
Early Maintenance Feedback	19
Instrument Layout	20

This chapter gives an introduction to the module, instrument overview and internal connectors.



Introduction to the Pump

The bio-inert quaternary pump comprises an optional solvent cabinet, a vacuum degasser and a four-channel gradient pump. The four-channel gradient pump comprises a high-speed proportioning valve and a pump assembly. It provides gradient generation by low pressure mixing. A solvent cabinet provides enough space for four one-liter bottles. The pump is optimized for the use of buffer solutions by the active seal wash function, which cleans pistons and seals.

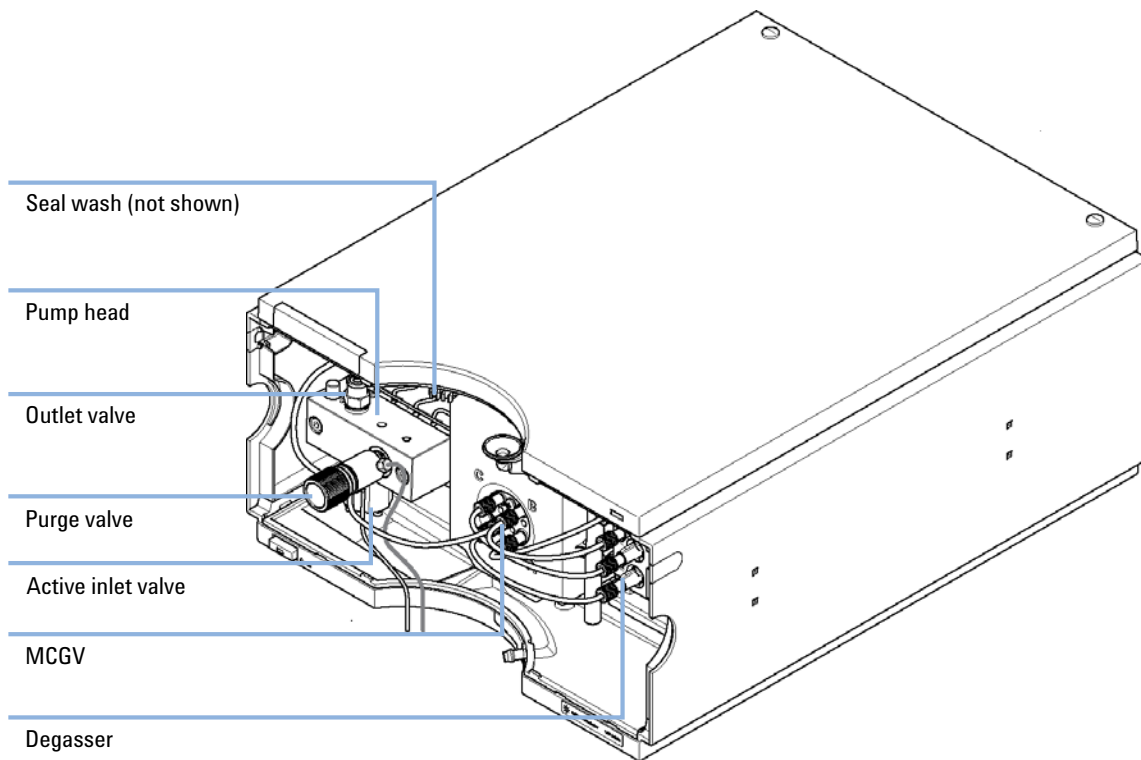


Figure 1 Overview of the bio-inert quaternary pump

Overview of the Hydraulic Path

The bio-inert quaternary pump is based on a two-channel, dual-piston in-series design which comprises all essential functions that a solvent delivery system has to fulfill. Metering of solvent and delivery to the high-pressure side are performed by one pump assembly which can generate pressure up to 600 bar.

Degassing of the solvents is done in a built-in vacuum degasser. Solvent compositions are generated on the low-pressure side by a high-speed proportioning valve (MCGV).

The pump assembly includes a pump head with an active inlet valve and an outlet valve. A damping unit is connected between the two piston chambers. A purge valve including a PTFE frit is fitted at the pump outlet for convenient priming of the pump head.

The pump includes an active seal wash function, which cleans pistons and seals for the use of buffer solutions.

Hydraulic Path

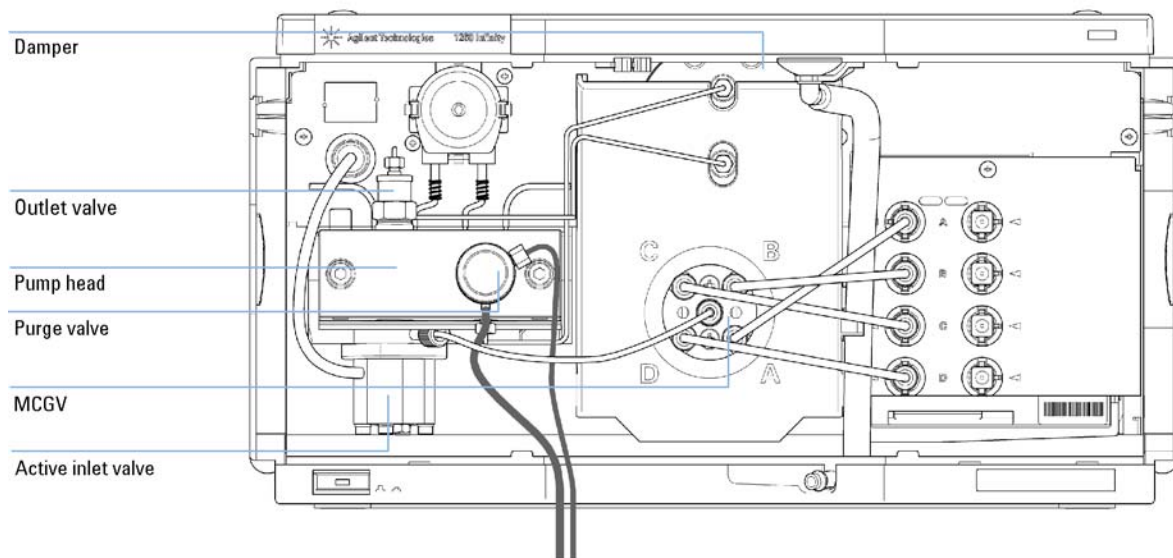


Figure 2 Hydraulic Path of the Bio-inert Quaternary Pump

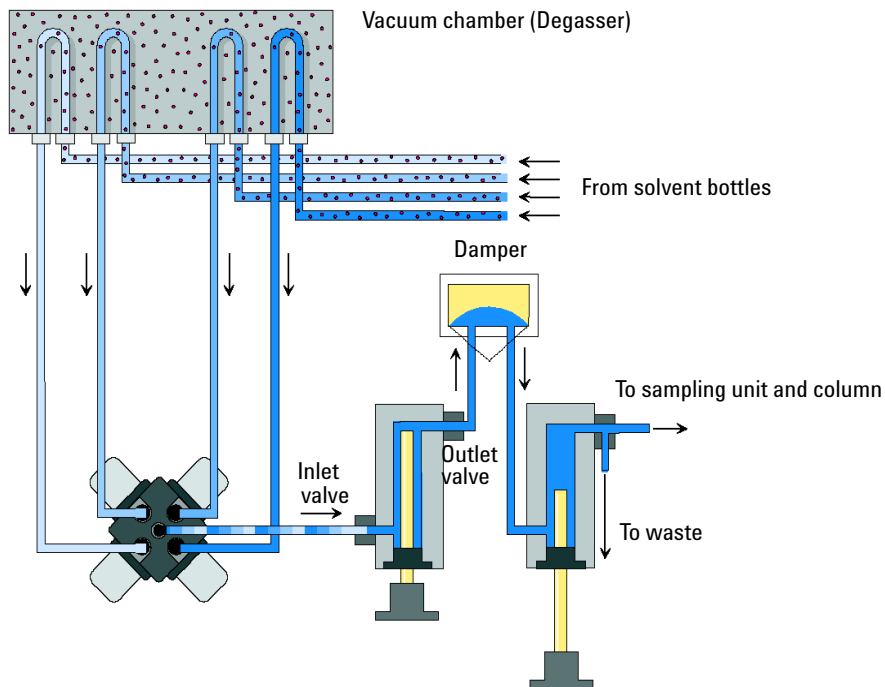


Figure 3 Hydraulic Path of the Quaternary Pump

How Does the Pump Work?

In the pump, the liquid runs from the solvent reservoir through the degasser to the MCGV and from there to the inlet valve.

The pump assembly comprises two substantially identical piston/chamber units. Both piston/chamber units comprise a ball-screw drive and a pump head with one sapphire piston for reciprocating movement in it.

A servo-controlled variable reluctance motor drives the two ball screw drives in opposite directions. The gears for the ball-screw drives have different circumferences (ratio 2:1) allowing the first piston to move at twice the speed of the second piston. The solvent enters the pump head close to the bottom limit and leaves it at its top. The outer diameter of the piston is smaller than the inner diameter of the pump head chamber allowing the solvent to fill the gap inbetween. The first piston has a stroke volume in the range of 20 – 100 μ l

1 Introduction

Overview of the Hydraulic Path

depending on the flow rate. The microprocessor controls all flow rates in a range of 1 $\mu\text{L}/\text{min}$ – 10 mL/min . The inlet of the first pumping unit is connected to the active inlet valve.

The outlet of the first piston/chamber unit is connected through the outlet valve and the damping unit to the inlet of the second piston/chamber unit. The outlet of the purge valve assembly is then connected to the following chromatographic system.

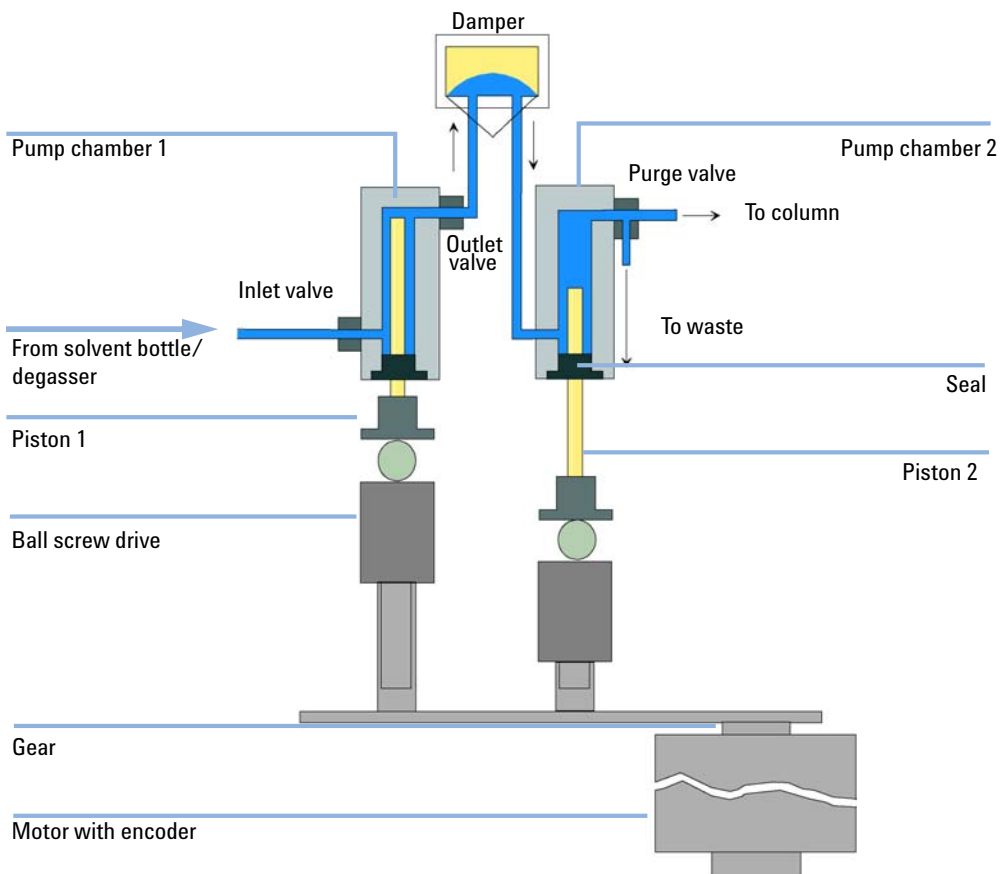


Figure 4 Principle of the Pump

When turned on, the pump runs through an initialization procedure to determine the upper dead position of the first piston. The first piston moves slowly upwards into the mechanical stop of the pump chamber and from there it moves back for a defined distance. The controller stores this piston position in memory. After this initialization the pump starts operation with the set parameters. The active inlet valve opens and the down-moving piston draws solvent into the first pump chamber. At the same time the second piston moves upwards delivering to the system. After a controller-defined stroke length that depends on the flow rate the drive motor is stopped and the active inlet valve closes. The motor direction is reversed and moves the first piston up until it reaches the stored upper limit and at the same time the second piston moves downwards. Then the sequence starts again moving the pistons up and down between the two limits. During the up movement of the first piston the solvent in the pump chamber is pressed through the outlet valve into the second pump chamber. The second piston draws in half of the volume displaced by the first piston and the remaining half volume is directly delivered to the system. During the drawing stroke of the first piston, the second piston delivers the drawn volume to the system.

For solvent compositions from the solvent bottles A, B, C, D the controller divides the length of the intake stroke in certain fractions in which the gradient valve connects the specified solvent channel to the pump input.

For specifications of the quaternary pump, see [“Performance Specifications”](#) on page 26.

How Does Compressibility Compensation Work?

The compressibility of the solvents in use will affect retention-time stability when the back pressure in the system changes (for example, ageing of column). In order to minimize this effect, the pump provides a compressibility compensation feature which optimizes the flow stability according to the solvent type. The compressibility compensation is set to a default value and can be changed through the user interface.

Without a compressibility compensation the following will happen during a stroke of the first piston. The pressure in the piston chamber increases and the volume in the chamber will be compressed depending on back pressure and solvent type. The volume displaced into the system will be reduced by the compressed volume.

With a compressibility value set the processor calculates a compensation volume, that depends on the back pressure of the system and the selected compressibility. This compensation volume will be added to the normal stroke volume and compensates the previously described *loss* of volume during the delivery stroke of the first piston.

How Does Variable Stroke Volume Work?

Due to the compression of the pump-chamber volume each piston stroke of the pump will generate a small pressure pulsation, influencing the flow stability of the pump. The amplitude of the pressure pulsation depends mainly on the stroke volume and the compressibility compensation for the solvent in use. Small stroke volumes generate pressure pulsations of smaller amplitude than higher stroke volumes at the same flow rate. In addition, the frequency of the pressure pulsations is higher. This decreases the influence of flow pulsations on quantitative results.

In gradient mode smaller stroke volumes result in a lower flow ripple improve composition ripple.

The module uses a processor-controlled spindle system for driving its pistons. The normal stroke volume is optimized for the selected flow rate. Small flow rates use a small stroke volume while higher flow rates use a higher stroke volume.

By default, the stroke volume for the pump is set to AUTO mode. This means that the stroke is optimized for the flow rate in use. A change to larger stroke volumes is possible but not recommended.

Bio-inert Materials

For the Agilent 1260 Infinity Bio-inert LC system, Agilent Technologies uses highest quality materials in the flow path (also referred to as wetted parts), which are widely accepted by life scientists, as they are known for optimum inertness towards biological samples and ensure best compatibility to common samples and solvents over a wide pH range. Explicitly, the complete flow path is free of stainless steel and free of other alloys containing metals like iron, nickel, cobalt, chromium, molybdenum or copper, which can interfere with biological samples. The flow downstream of the sample introduction does not contain metals at all.

Table 1 Bio-inert Materials

Module	Materials
Agilent 1260 Infinity Bio-inert Quaternary Pump (G5611A)	Titanium, gold, platinum-iridium, zirconium oxide, ruby, PTFE, PEEK
Agilent 1260 Infinity Bio-inert High-Performance Autosampler (G5667A)	Upstream of sample-introduction: <ul style="list-style-type: none">• Titanium, gold, PTFE, PEEK, zirconium oxide Downstream of sample-introduction: <ul style="list-style-type: none">• PEEK, zirconium oxide
Agilent 1260 Infinity Bio-inert Manual Injector (G5628A)	PEEK, zirconium oxide
Agilent 1260 Infinity Bio-inert Analytical Fraction Collector (G5664A)	PEEK, zirconium oxide, PTFE

Table 1 Bio-inert Materials

Module	Materials
<p>Bio-inert Flow Cells: Standard flow cell bio-inert, 10 mm, 13 µl, 120 bar (12 MPa) for MWD/DAD, includes Capillary Kit Flow Cells BIO (p/n G5615-68755) (p/n G5615-60022) (for Agilent 1260 Infinity Diode Array Detectors DAD G1315C/D)</p>	PEEK, zirconium oxide, sapphire, PTFE
<p>Max-Light Cartridge Cell Bio-inert (10 mm, V(σ) 1.0 µl) (p/n G5615-60018) and Max-Light Cartridge Cell Bio-inert (60 mm, V(σ) 4.0 µl) (p/n G5615-60017) (for Agilent 1290 Infinity Diode Array Detectors DAD G4212A/B)</p>	PEEK, fused silica
<p>Bio-inert flow cell, 8 µL, 20 bar (pH 1–12) includes Capillary Kit Flow Cells BIO (p/n G5615-68755) (p/n G5615-60005) (for Agilent 1260 Infinity Fluorescence Detector FLD G1321B)</p>	PEEK, fused silica, PTFE
<p>Bio-inert heat-exchanger G5616-60050 for 1290 Infinity Thermostatted Column Compartment (G1316C)</p>	PEEK (steel-cladded)
<p>Bio-inert Valve heads</p>	G4235A, G5631A, G5639A: PEEK, ceramic (Al ₂ O ₃ based)
<p>Bio-inert Connection capillaries</p>	<p>Upstream of sample-introduction:</p> <ul style="list-style-type: none"> • Titanium <p>Downstream of sample-introduction:</p> <ul style="list-style-type: none"> • Agilent uses stainless-steel-cladded PEEK capillaries, which keep the flow path free of steel and provide pressure stability to more than 600 bar.

For ensuring optimum bio-compatibility of your Agilent 1260 Infinity Bio-inert LC system, do not include non-inert standard modules or parts to the flow path. Do not use any parts that are not labeled as Agilent “Bio-inert”. For solvent compatibility of these materials, see [“Solvent information for parts of the 1260 Infinity Bio-inert LC system”](#) on page 51.

Early Maintenance Feedback

Maintenance requires the exchange of components which are subject to wear or stress. Ideally, the frequency at which components are exchanged should be based on the intensity of usage of the module and the analytical conditions, and not on a predefined time interval. The early maintenance feedback (**EMF**) feature monitors the usage of specific components in the instrument, and provides feedback when the user-selectable limits have been exceeded. The visual feedback in the user interface provides an indication that maintenance procedures should be scheduled.

EMF Counters

EMF counters increment with use and can be assigned a maximum limit which provides visual feedback in the user interface when the limit is exceeded. Some counters can be reset to zero after the required maintenance procedure.

Using the EMF Counters

The user-settable **EMF** limits for the **EMF Counters** enable the early maintenance feedback to be adapted to specific user requirements. The useful maintenance cycle is dependent on the requirements for use. Therefore, the definition of the maximum limits need to be determined based on the specific operating conditions of the instrument.

Setting the EMF Limits

The setting of the **EMF** limits must be optimized over one or two maintenance cycles. Initially the default **EMF** limits should be set. When instrument performance indicates maintenance is necessary, take note of the values displayed by the **EMF counters**. Enter these values (or values slightly less than the displayed values) as **EMF** limits, and then reset the **EMF counters** to zero. The next time the **EMF counters** exceed the new **EMF** limits, the **EMF** flag will be displayed, providing a reminder that maintenance needs to be scheduled.

Instrument Layout

The industrial design of the module incorporates several innovative features. It uses Agilent's E-PAC concept for the packaging of electronics and mechanical assemblies. This concept is based upon the use of expanded polypropylene (EPP) layers of foam plastic spacers in which the mechanical and electronic boards components of the module are placed. This pack is then housed in a metal inner cabinet which is enclosed by a plastic external cabinet. The advantages of this packaging technology are:

- virtual elimination of fixing screws, bolts or ties, reducing the number of components and increasing the speed of assembly/disassembly,
- the plastic layers have air channels molded into them so that cooling air can be guided exactly to the required locations,
- the plastic layers help cushion the electronic and mechanical parts from physical shock, and
- the metal inner cabinet shields the internal electronics from electromagnetic interference and also helps to reduce or eliminate radio frequency emissions from the instrument itself.



2 Site Requirements and Specifications

Site Requirements	22
Physical Specifications	25
Performance Specifications	26

This chapter provides information on environmental requirements, physical and performance specifications.



Site Requirements

A suitable environment is important to ensure optimal performance of the instrument.

Power Considerations

The module power supply has wide ranging capability. It accepts any line voltage in the range described in [Table 2](#) on page 25. Consequently there is no voltage selector in the rear of the module. There are also no externally accessible fuses, because automatic electronic fuses are implemented in the power supply.

WARNING

Hazard of electrical shock or damage of your instrumentation can result, if the devices are connected to a line voltage higher than specified.

→ Connect your instrument to the specified line voltage only.

WARNING

The module is partially energized when switched off, as long as the power cord is plugged in.

Repair work at the module can lead to personal injuries, e.g. electrical shock, when the cover is opened and the module is connected to power.

→ Always unplug the power cable before opening the cover.

→ Do not connect the power cable to the instrument while the covers are removed.

CAUTION

Inaccessible power plug.

In case of emergency it must be possible to disconnect the instrument from the power line at any time.

- Make sure the power connector of the instrument can be easily reached and unplugged.
 - Provide sufficient space behind the power socket of the instrument to unplug the cable.
-

Power Cords

Different power cords are offered as options with the module. The female end of all power cords is identical. It plugs into the power-input socket at the rear. The male end of each power cord is different and designed to match the wall socket of a particular country or region.

WARNING

Absence of ground connection or use of unspecified power cord

The absence of ground connection or the use of unspecified power cord can lead to electric shock or short circuit.

- Never operate your instrumentation from a power outlet that has no ground connection.
 - Never use a power cord other than the Agilent Technologies power cord designed for your region.
-

WARNING

Use of unsupplied cables

Using cables not supplied by Agilent Technologies can lead to damage of the electronic components or personal injury.

- Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.
-

WARNING

Unintended use of supplied power cords

Using power cords for unintended purposes can lead to personal injury or damage of electronic equipment.

- Never use the power cords that Agilent Technologies supplies with this instrument for any other equipment.
-

Bench Space

The module dimensions and weight (see [Table 2](#) on page 25) allow you to place the module on almost any desk or laboratory bench. It needs an additional 2.5 cm (1.0 inches) of space on either side and approximately 8 cm (3.1 inches) in the rear for air circulation and electric connections.

If the bench shall carry a complete HPLC system, make sure that the bench is designed to bear the weight of all modules.

The module should be operated in a horizontal position.

Condensation

CAUTION

Condensation within the module

Condensation will damage the system electronics.

- Do not store, ship or use your module under conditions where temperature fluctuations could cause condensation within the module.
 - If your module was shipped in cold weather, leave it in its box and allow it to warm slowly to room temperature to avoid condensation.
-

Physical Specifications

Table 2 Physical Specifications


Type	Specification	Comments
Weight	14.5 kg (32 lbs)	
Dimensions (height × width × depth)	180 x 345 x 435 mm (7.0 x 13.5 x 17 inches)	
Line voltage	100 – 240 VAC, ± 10%	Wide-ranging capability
Line frequency	50 or 60 Hz, ± 5%	
Power consumption	180 VA, 110W / 375 BTU	
Ambient operating temperature	4–55 °C (41–131 °F)	
Ambient non-operating temperature	-40–70 °C (-4–158 °F)	
Humidity	< 95%, at 25–40 °C (77–104 °F)	Non-condensing
Operating altitude	Up to 2000 m (6562 ft)	
Non-operating altitude	Up to 4600 m (15091 ft)	For storing the module
Safety standards: IEC, CSA, UL	Installation category II, Pollution degree 2	For indoor use only.

Performance Specifications

Table 3 Specifications - Agilent 1260 Infinity Bio-inert Quaternary Pump (G5611A)

Type	Specifications
Hydraulic system	Dual piston in series pump with proprietary servo-controlled variable stroke drive, floating pistons and active inlet valve, integrated 4-channel degassing unit
Setable flow range	0.001 – 10 mL/min, in 0.001 mL/min increments
Recommended flow range	0.2 – 10 mL/min
Flow precision	< 0.07 % RSD, or < 0.02 min SD whatever is greater, based on retention time at constant room temperature
Flow accuracy	± 1 % or 10 µL/min whatever is greater
Pressure	Operating range up to 60 MPa (600 bar, 8700 psi) up to 5 mL/min Operating range up to 20 MPa (200 bar, 2950 psi) up to 10 mL/min
Pressure pulsation	< 2 % amplitude (typically < 1.3 %), at 1 mL/min isopropanol, at all pressures > 1 MPa (10 bar)
Compressibility compensation	User-selectable, based on mobile phase compressibility
Recommended pH range	1 – 13 , short term 14 ¹
Gradient formation	Low pressure quaternary mixing/gradient capability using proprietary high-speed proportioning valve. Delay volume 600 – 900 µL, dependent on back pressure.
Composition range	0 – 95 % or 5 – 100 %, user selectable
Composition precision	< 0.2 % RSD, at 0.2 and 1 mL/min
Communications	Controller-area network (CAN), RS-232C, APG Remote: ready, start, stop and shut-down signals, LAN optional
Materials in flow path	Titanium, Gold, Platin-Iridium, Sapphire, PEEK, PTFE
Active seal wash	Included

¹ For solvent compatibility, refer to section "Solvent information for parts of the 1260 Infinity Bio-inert LC system" in the manual

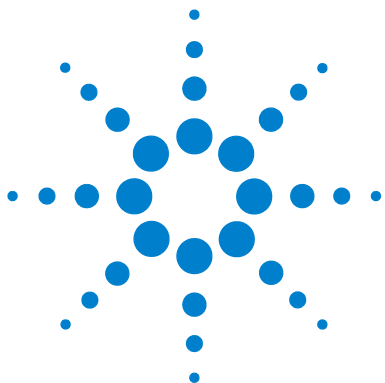

3 Installing the Pump

Unpacking the Pump	28
Delivery Checklist	28
Accessory Kit	29
Optimizing the Stack Configuration	30
One Stack Configuration	30
Installing the Pump	33
Connecting Modules and Control Software	36
Connecting Modules	36
Connecting Control Software and/or G4208 A Instant Pilot	37
Flow Connections of the Pump	38
Priming the System	41
Initial Priming	41
Regular Priming	43
Changing Solvents	44

This chapter gives information about the preferred stack setup for your system and the installation of your module.



Unpacking the Pump

If the delivery packaging shows signs of external damage, please call your Agilent Technologies sales and service office immediately. Inform your service representative that the instrument may have been damaged during shipment.

CAUTION

"Defective on arrival" problems

If there are signs of damage, please do not attempt to install the module. Inspection by Agilent is required to evaluate if the instrument is in good condition or damaged.

- Notify your Agilent sales and service office about the damage.
- An Agilent service representative will inspect the instrument at your site and initiate appropriate actions.

Delivery Checklist

Ensure all parts and materials have been delivered with the pump. The delivery checklist is shown below. To aid in parts identification, please refer to chapter Parts and Materials for Maintenance. Please report missing or damaged parts to your local Agilent Technologies sales and service office.

Delivery Checklist

p/n	Description
G4800-64500	Agilent LC Hardware Documentation DVD
G1311-60003 (4x)	Bottle-head assembly
M85550	Lab Advisor CD
5065-9981	Solvent cabinet, including all plastic parts
G4203-68708	HPLC System Tool Kit
G5611-68707	Starter-Kit Bio-inert HPLC
959961-902	Column Eclipse Plus C18, 4.6 x 100 mm, 3.5 µm (optional)

p/n	Description
699975-902	Column Poroshell 120 EC-C18, 4.6 x 50 mm, 2.7 µm (optional)
883975-902	Column SB-C18, 4.6 x 150 mm, 5 µm (optional)
G1369-60002	LAN Communication Card
	Power cord

Accessory Kit

Accessory Kit (Bio-inert) (p/n G5611-68755)

p/n	Description
5062-2461	Waste tube, 5 m (reorder pack)
5063-6527	Tubing assembly, i.d. 6 mm, o.d. 9 mm, 1.2 m (to waste)
5181-1519	CAN cable, Agilent module to module, 1 m
G5611-60502	Capillary 900 x 0.17 mm, titanium (Bio-inert) pump to thermostatted autosampler
5042-9954	Tubing clip (2x), re-order 4/pk
G5611-605003	Capillary 400 x 0.17 mm, titanium (Bio-inert) pump to injector

Optimizing the Stack Configuration

If your module is part of a complete Agilent 1260 Infinity Liquid Chromatograph, you can ensure optimum performance by installing the following configurations. These configurations optimize the system flow path, ensuring minimum delay volume.

One Stack Configuration

Ensure optimum performance by installing the modules of the Agilent 1260 Infinity LC System in the following configuration (see [Figure 5](#) on page 31 and [Figure 6](#) on page 32). This configuration optimizes the flow path for minimum delay volume and minimizes the bench space required.

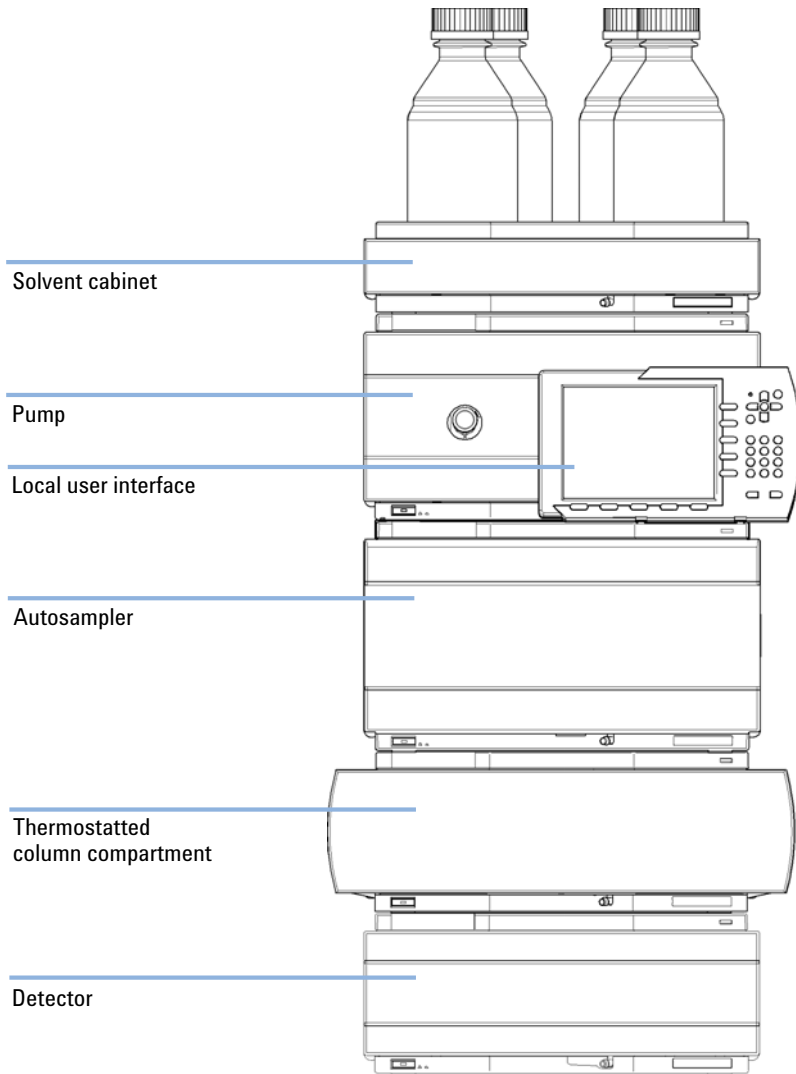


Figure 5 Recommended Stack Configuration (Front View)

3 Installing the Pump

Optimizing the Stack Configuration

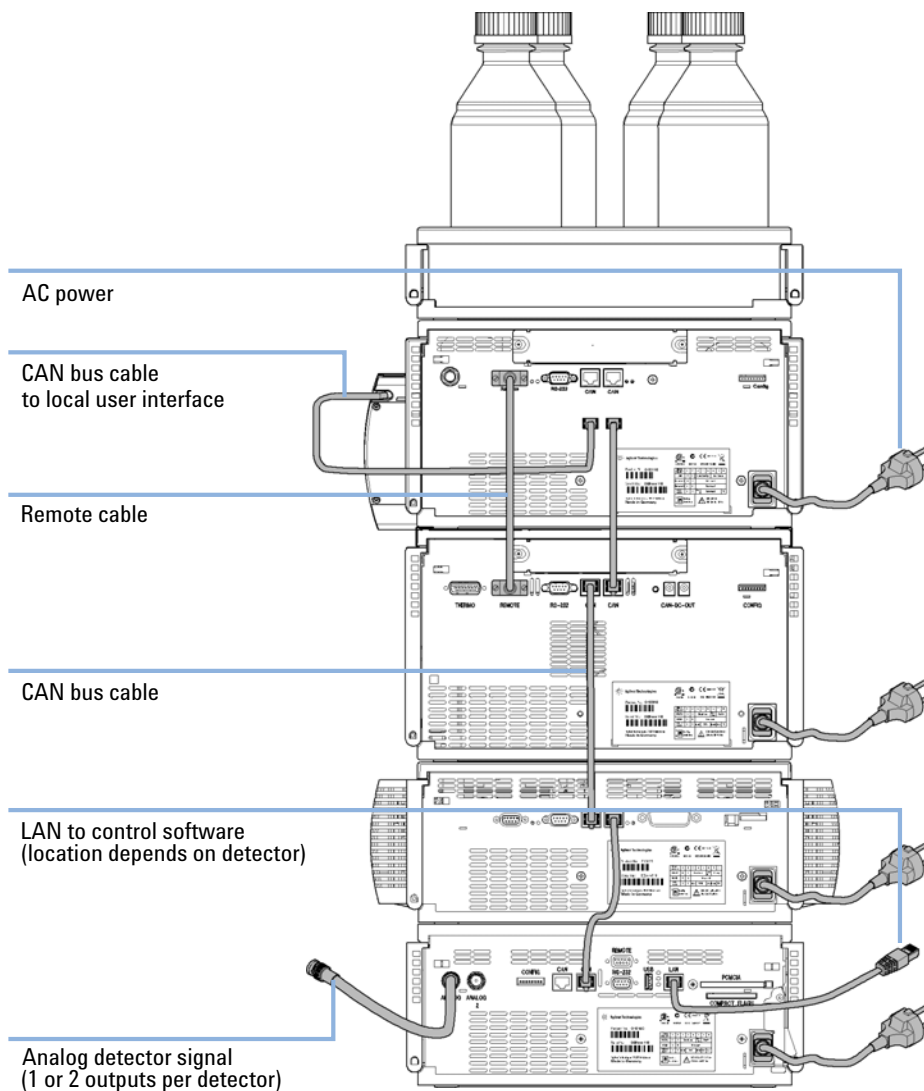


Figure 6 Recommended Stack Configuration (Rear View)

Installing the Pump

Parts required	#	p/n	Description
	1		Pump
	1		Data System
	1	G4208A	Instant Pilot
	1		Power cord

For other cables see text below and “[Cable Overview](#)” on page 154.

- Preparations**
- Locate bench space.
 - Provide power connections.
 - Unpack the module.

WARNING

Module is partially energized when switched off, as long as the power cord is plugged in.

Repair work at the module can lead to personal injuries, e.g. shock hazard, when the cover is opened and the module is connected to power.

- Make sure that it is always possible to access the power plug.
- Remove the power cable from the instrument before opening the cover.
- Do not connect the power cable to the Instrument while the covers are removed.

CAUTION

"Defective on arrival" problems

If there are signs of damage, please do not attempt to install the module. Inspection by Agilent is required to evaluate if the instrument is in good condition or damaged.

- Notify your Agilent sales and service office about the damage.
- An Agilent service representative will inspect the instrument at your site and initiate appropriate actions.

3 Installing the Pump

Installing the Pump

- 1 Place the module on the bench in a horizontal position.
- 2 Ensure the power switch on the front of the module is OFF (switch stands out).


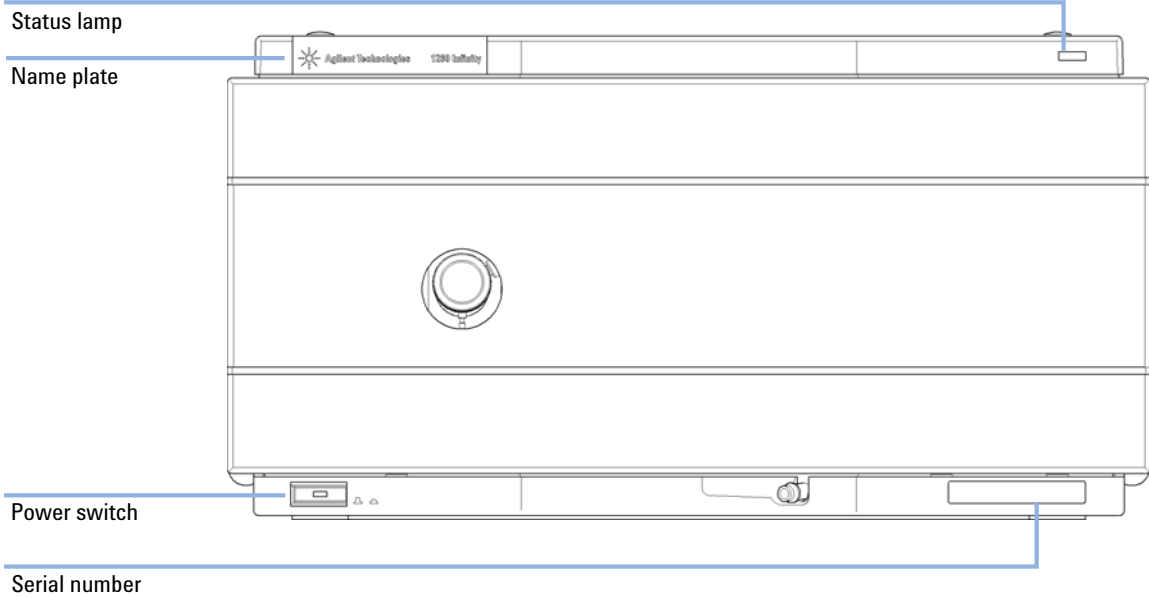


Figure 7 Front of Pump

- 3 Connect the power cable to the power connector at the rear of the module.

- 4 Connect the required interface cables to the quaternary pump, see “Connecting Modules and Control Software” on page 36.

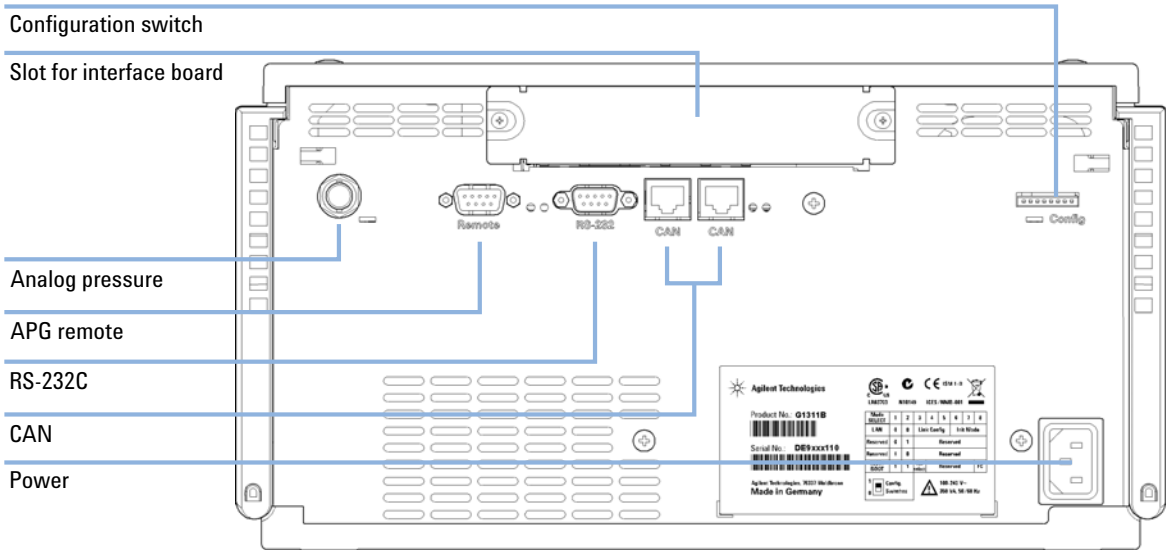

- 5 Connect all capillaries, solvent tubes and waste tubing (see “Flow Connections of the Pump” on page 38).
- 6 Press the power switch to turn on the module.

NOTE

The power switch stays pressed in and a green indicator lamp in the power switch is on when the module is turned on. When the line power switch stands out and the green light is off, the module is turned off.

- 7 Purge the quaternary pump (see “Initial Priming” on page 41).

NOTE

The pump was shipped with default configuration settings. To change these settings, see “Setting the 8-bit Configuration Switch (without On-board) LAN” on page 176.

Connecting Modules and Control Software

WARNING

Use of unsupplied cables

Using cables not supplied by Agilent Technologies can lead to damage of the electronic components or personal injury.

- Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.
-

Connecting Modules

- 1 Place the individual modules in a stack configuration as shown in [Figure 5](#) on page 31.
- 2 Ensure the power switches on the front of the modules are OFF (switches stand out).
- 3 Plug a CAN cable into the CAN connector at the rear of the respective module (except vacuum degasser).
- 4 Connect the CAN cable to the CAN connector of the next module, see [Figure 6](#) on page 32.
- 5 Press in the power switches to turn on the modules.

Connecting Control Software and/or G4208 A Instant Pilot

NOTE

With the introduction of the Agilent 1260 Infinity, all GPIB interfaces have been removed. The preferred communication is LAN.

NOTE

Usually the detector is producing the most data in the stack, followed by the pump, and it is therefore highly recommended to use either of these modules for the LAN connection.

- 1 Ensure the power switches on the front of the modules in the stack are OFF (switches stand out).
- 2 If there are no other 1260 with LAN port in the HPLC stack, install a G1369B LAN board into the extension slot of the pump.
- 3 Connect the LAN enabled module with a LAN cable to the data system.
- 4 Plug the CAN connector of the Instant Pilot into any available CAN port of the 1260 system.
- 5 Plug a CAN cable into the CAN connector of the Instant Pilot.
- 6 Connect the CAN cable to the CAN connector of one of the modules.
- 7 Press in the power switches to turn on the modules.

NOTE

The Agilent control software can also be connected to the system through a LAN cable, which requires the installation of a LAN-board. For more information about connecting the Instant Pilot or Agilent control software refer to the respective user manual. For connecting the Agilent 1260 Infinity equipment to non-Agilent 1260 Infinity equipment, see [“Introduction to the Pump”](#) on page 10.

Flow Connections of the Pump

Parts required	#	Description
	1	Other modules
	1	Parts from accessory kit
	2	wrenches 1/4 - 5/16 inch for capillary connections

Preparations Pump is installed in the LC system


WARNING

Toxic, flammable and hazardous solvents, samples and reagents

The handling of solvents, samples and reagents can hold health and safety risks.

- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor, and follow good laboratory practice.
 - The volume of substances should be reduced to the minimum required for the analysis.
 - Do not operate the instrument in an explosive atmosphere.
-

- 1 Remove the front cover by pressing the snap fasteners on both sides.

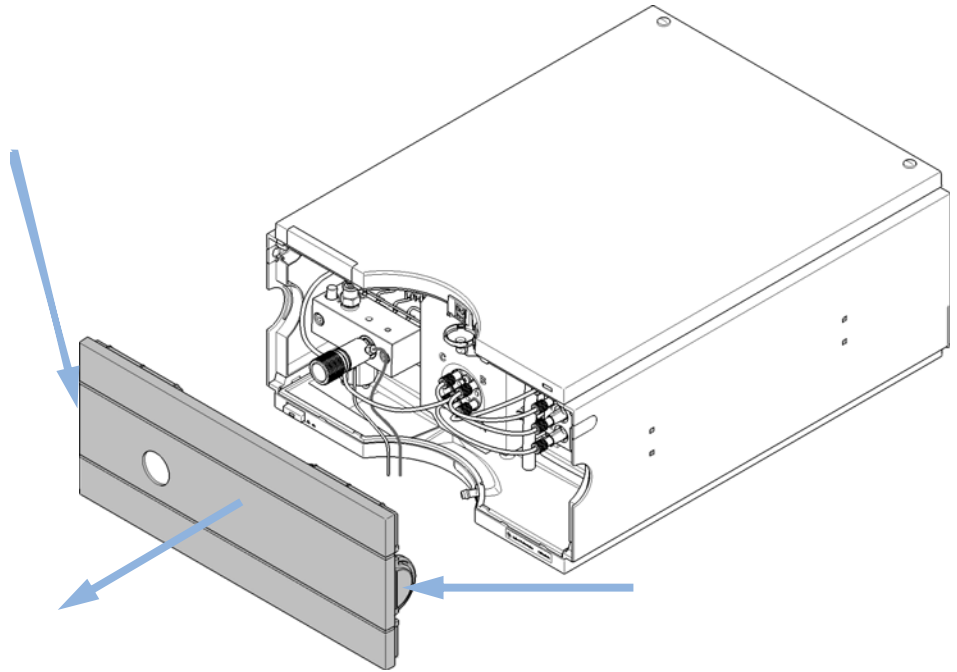

- 2 Place the solvent cabinet on top of the quaternary pump.
- 3 Put the bottle-head assemblies into solvent reservoirs containing your mobile phase and place the bottle in the solvent cabinet.
- 4 Connect the inlet tubes from the bottle-head assemblies to the inlet connectors A to D (typically the left connection of the channel) of the vacuum degasser. Fix the tubes in the tube clips of the pump.
- 5 Connect the solvent tubes to the outlet connectors (typically right connection of the channel) of the vacuum degasser.
- 6 Using a piece of sanding paper connect the waste tubing to the purge valve and place it into your waste system.
- 7 If the pump is not part of an Agilent 1260 Infinity system stack or placed on the bottom of a stack, connect the corrugated waste tube to the waste outlet of the pump leak handling system.
- 8 Connect the pump outlet capillary (pump to injection device) to the outlet of the purge valve.

3 Installing the Pump

Flow Connections of the Pump

9 Prime your system before first use (see “Initial Priming” on page 41).

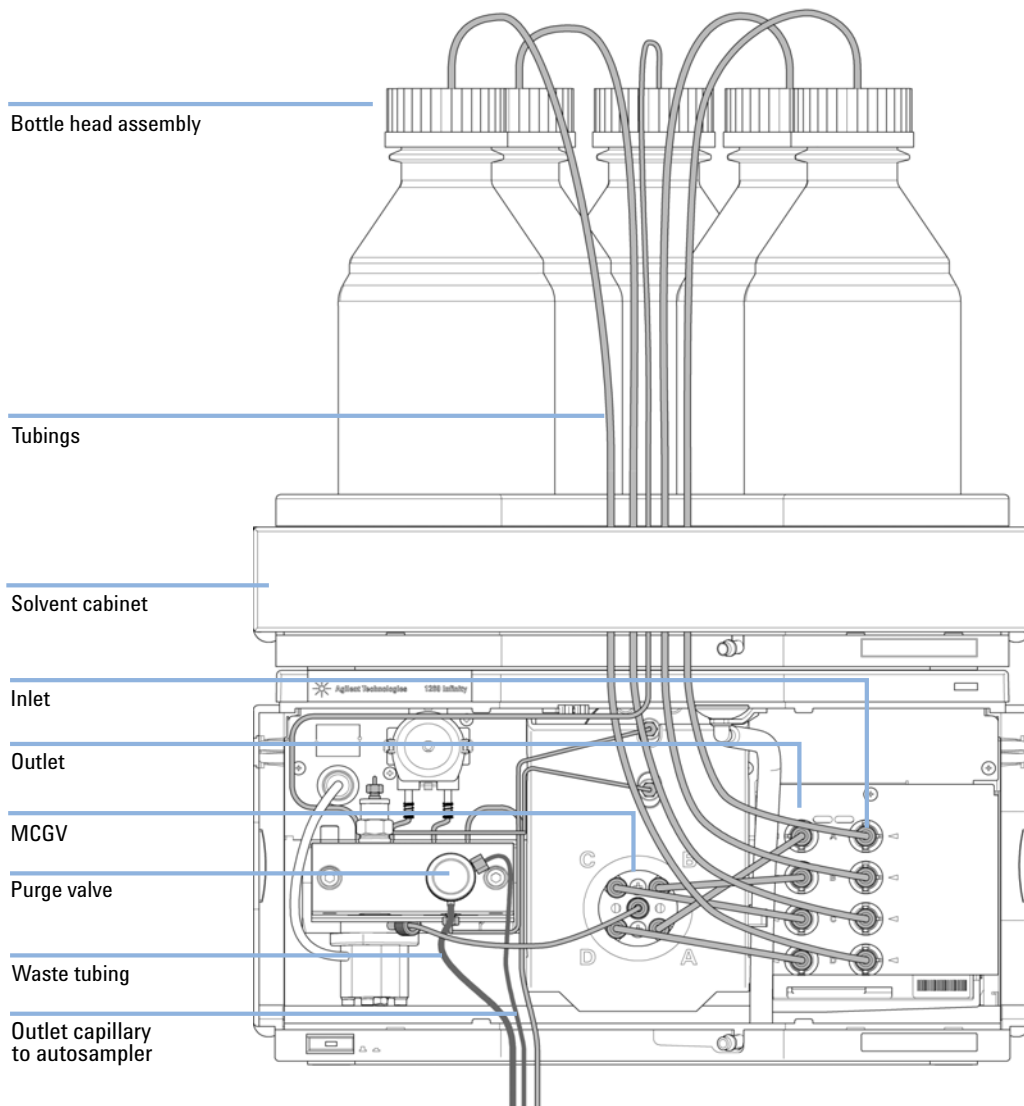


Figure 8 Flow Connections of the Pump

Priming the System

Initial Priming

When Before a new degasser or new solvent tubing can be used, it is necessary to prime the system. Isopropanol (IPA) is recommended as priming solvent due to its miscibility with nearly all HPLC solvents and its excellent wetting properties.

Parts required

#	Description
1	Isopropanol

Preparations Connect all modules hydraulically as described in the respective module manuals.
Fill each solvent bottle with 100 mL isopropanol
Switch the system on

WARNING

When opening capillary or tube fittings solvents may leak out.

The handling of toxic and hazardous solvents and reagents can bear health risks.

→ Please observe appropriate safety procedures (for example, goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the solvent vendor, especially when toxic or hazardous solvents are used.

NOTE

The purge tool of the LabAdvisor or Instrument Utilities can be used for automatically purging the pump.

NOTE

If the pump is not able to aspirate the solvent from the bottles, a syringe can be used to draw the solvent manually through tubing and degasser.

3 Installing the Pump

Priming the System

NOTE

When priming the vacuum degasser with a syringe, the solvent is drawn through the degasser tubes very quickly. The solvent at the degasser outlet will therefore not be fully degassed. Pump for approximately 10 minutes at your desired flow rate before starting an analysis. This will allow the vacuum degasser to properly degas the solvent in the degasser tubes.

- 1 Open the purge valve of the pump
- 2 Set the flow rate to 5 mL/min.
- 3 Select channel A1
- 4 Turn the flow on
- 5 Observe if the solvent in the tubing of channel A1 is advancing towards the pump. If it isn't, disconnect the solvent tubing from the solvent selection valve, attach a syringe with a syringe adapter and pull the liquid through the degasser. Reattach the tubing to the solvent selection valve.
- 6 Pump 30 mL isopropanol to remove residual air bubbles.
- 7 Switch to the next solvent channel and repeat steps 5 and 6 until all channels have been purged.
- 8 Turn the flow off and close the purge valve.

Regular Priming

When When the pumping system has been turned off for a certain time (for example, overnight) air will rediffuse into the solvent channel between the vacuum degasser and the pump. Solvents containing volatile ingredients will slightly lose these if left in the degasser without flow for a prolonged period of time.

Preparations Switch the system on

NOTE

The purge tool of the LabAdvisor or Instrument Utilities can be used for automatically purging the pump.

- 1 Open the purge valve of your pump by turning it counterclockwise and set the flow rate to 5 mL/min.
- 2 Flush the vacuum degasser and all tubes with at least 10 mL of solvent.
- 3 Repeat step 1 and 2 for the other channel(s) of the pump.
- 4 Set the required composition and flow rate for your application and close the purge valve.
- 5 Pump for approximately 10 minutes before starting your application.

Changing Solvents

When When the solvent of a channel is to be replaced by another solvent that is not compatible (solvents are immiscible or one solvent contains a buffer) it is necessary to follow the procedure below to prevent clogging of the pump by salt precipitation or residual liquid droplets in parts of the system.

Parts required	#	p/n	Description
	1		Purging solvent(s), see Table 4 on page 45
	1	5067-4741	ZDV union (Bio-inert)

Preparations Remove the column and replace it by a ZDV fitting.
Prepare bottles with appropriate intermediate solvents (see [Table 4](#) on page 45)

- 1 If the channel is not filled with buffer, proceed to step 4.
- 2 Place the solvent intake filter into a bottle of water.
- 3 Flush the channel at a flow rate suitable for the installed tubing (typically 3-5 mL/min) for 10 min.
- 4 Modify the flow path of your system as required for your application. For delay volume optimization see the Rapid Resolution System manual.

CAUTION

Buffer salt of aqueous buffers may precipitate in residual isopropanol.

Capillaries and filter may be clogged by precipitating salt.

→ Don't perform steps 5 to 7 for channels run with aqueous buffer as solvent.

- 5 Replace the solvent bottle by a bottle of isopropanol.
- 6 Flush the channel at a flow rate suitable for the installed tubing (typically 3-5 mL/min) for 5 min.
- 7 Swap the bottle of isopropanol with a bottle of solvent for your application.
- 8 Repeat steps 1 to 7 for the other channel(s) of the pump.
- 9 Install the desired column, set the required composition and flow rate for your application and equilibrate the system for approx. 10 minutes prior to starting a run.

Table 4 Choice of Priming Solvents for Different Purposes

Activity	Solvent	Comments
After an installation	Isopropanol	Best solvent to flush air out of the system
When switching between reverse phase and normal phase (both times)	Isopropanol	Miscible with almost all solvents
After an installation	Ethanol or methanol	Alternative to isopropanol (second choice) if no isopropanol is available
To clean the system when using buffers	HPLC grade water	Best solvent to re-dissolve buffer crystals
After changing aqueous solvents	HPLC grade water	Best solvent to re-dissolve buffer crystals
After the installation of normal phase seals (P/N 0905-1420)	Hexane + 5% isopropanol	Good wetting properties

3 **Installing the Pump** Priming the System



4 Using the Pump

Hints for Successful Use of the Quaternary Pump 48

Prevent Blocking of Solvent Filters 49

Algae Growth in HPLC Systems 50

How to Prevent and/or Reduce the Algae Problem 50

Solvent Information 51

This chapter explains the operational parameters of the module.



Hints for Successful Use of the Quaternary Pump

- Always place the solvent cabinet with the solvent bottles on top of the quaternary pump (or at a higher level).
- When using salt solutions and organic solvents in the quaternary pump it is recommended to connect the salt solution to one of the bottom gradient valve ports and the organic solvent to one of the upper gradient valve ports. It is best to have the organic channel directly above the salt solution channel. Regular flushing of all MCGV channels with water is recommended to remove all possible salt deposits in the valve ports.
- Before operating the quaternary pump, flush the pump and vacuum degasser, see “[Regular Priming](#)” on page 43). This is especially recommended if it has been turned off for some time (for example, overnight) and volatile solvent mixtures are used in the channels.
- Prevent blocking of solvent inlet filters. Never use the pump without solvent inlet filter. Prevent the growth of algae, see “[Prevent Blocking of Solvent Filters](#)” on page 49).
- Regularly check the purge valve frit and column frit. A blocked purge valve frit can be identified by a black or yellow surface, deposits or by a pressure greater than 10 bar, when pumping distilled water at a rate of 5 mL/min with an open purge valve.
- When using the quaternary pump at low flow rates (for example, 0.2 mL/min) check all 1/16-inch fittings for any signs of leaks.
- Whenever exchanging the pump seals the purge valve frit should be exchanged, too.
- When using buffers or other salt solutions, flush the system with water before switching it off. The seal wash option should be used when salt concentrations of 0.1 M or higher will be used for long time periods.
- Check the pump pistons for scratches when changing the piston seals. Scratched pistons will cause micro leaks and will decrease the lifetime of the seal.
- Pressurize the system according to the wear in procedure after changing the piston seals (see “[Maintenance of the Pump Head](#)” on page 126).
- Consider recommendations given in the solvent information section, see “[Solvent Information](#)” on page 51.

Prevent Blocking of Solvent Filters

Contaminated solvents or algae growth in the solvent bottle will reduce the lifetime of the solvent filter and will influence the performance of the module. This is especially true for aqueous solvents or phosphate buffers (pH 4 to 7). The following suggestions will prolong lifetime of the solvent filter and will maintain the performance of the module.

- Use a sterile, if possible amber, solvent bottle to slow down algae growth.
- Filter solvents through filters or membranes that remove algae.
- Exchange solvents every two days or refilter.
- If the application permits add 0.0001-0.001M sodium azide to the solvent.
- Place a layer of argon on top of your solvent.
- Avoid exposure of the solvent bottle to direct sunlight.

NOTE

Never use the system without solvent filter installed.

Algae Growth in HPLC Systems

The presence of algae in HPLC systems can cause a variety of problems that may be incorrectly diagnosed as instrument or application problems. Algae grow in aqueous media, preferably in a pH range of 4-8. Their growth is accelerated by buffers, for example phosphate or acetate. Since algae grow through photosynthesis, light will also stimulate their growth. Even in distilled water small-sized algae grow after some time.

Instrumental Problems Associated With Algae

Algae deposit and grow everywhere within the HPLC system causing:

- Blocked solvent filters or deposits on inlet or outlet valves resulting in unstable flow, composition or gradient problems or a complete failure of the pump.
- Small pore high pressure solvent filters, usually placed before the injector to plug resulting in high system pressure.
- PTFE frits blockage leading to increased system pressure.
- Column filters to plug giving high system pressure.
- Flow cell windows of detectors to become dirty resulting in higher noise levels (since the detector is the last module in the flow path, this problem is less common).

How to Prevent and/or Reduce the Algae Problem

- Always use freshly prepared solvents, especially use demineralized water which was filtered through about 0.2 μm filters.
- Never leave mobile phase in the instrument for several days without flow.
- Always discard old mobile phase.
- Use the amber solvent bottle (Solvent bottle, amber (p/n 9301-1450)) supplied with the instrument for your aqueous mobile phase.
- If possible add a few mg/l sodium azide or a few percent organic solvent to the aqueous mobile phase.

Solvent Information

Solvent Information

Observe the following recommendations on the use of solvents.

- Follow recommendations for avoiding the growth of algae, see [“Algae Growth in HPLC Systems”](#) on page 50
- Small particles can permanently block capillaries and valves. Therefore always filter solvents through 0.4 µm filters.
- Avoid or minimize the use of solvents, which may corrode parts in the flow path. Consider specifications for the pH range given for different materials like flow cells, valve materials etc. and recommendations in subsequent sections.

Solvent information for parts of the 1260 Infinity Bio-inert LC system

For the Agilent 1260 Infinity Bio-inert LC system, Agilent Technologies uses highest quality materials (see [“Bio-inert Materials”](#) on page 17) in the flow path (also referred to as wetted parts), which are widely accepted by life scientists, as they are known for optimum inertness towards biological samples and ensure best compatibility to common samples and solvents over a wide pH range. Explicitly, the complete flow path is free of stainless steel and free of other alloys containing metals like iron, nickel, cobalt, chromium, molybdenum or copper, which can interfere with biological samples. The flow downstream of the sample introduction does not contain metals at all.

However there are no materials, which combine suitability for versatile HPLC instrumentation (valves, capillaries, springs, pump heads, flow cells etc.) with complete compatibility to all possible chemicals and application conditions. This section gives recommendations, which solvents should be preferred. Chemicals which are known to cause issues should be avoided or exposure should be minimized, e.g. for short term cleaning procedures. After using potentially aggressive chemicals, the system should be flushed with compatible standard HPLC solvents.

PEEK

PEEK (Polyether-Ether Ketones) combines excellent properties regarding biocompatibility, chemical resistance, mechanical and thermal stability and is therefore the material of choice for biochemical instrumentation. It is stable in the specified pH range and inert to many common solvents. There is still a number of known incompatibilities to chemicals like chloroform, methylene chloride, THF, DMSO, strong acids (nitric acid > 10 %, sulphuric acid > 10 %, sulfonic acids, trichloroacetic acid), halogenes or aqueous halogene solutions, phenol and derivatives (cresols, salicylic acid etc.).

When used above room temperature, it is sensitive to bases and various organic solvents, which can cause PEEK to swell. As normal PEEK capillaries are very sensitive to high pressure especially under such conditions, Agilent uses stainless-steel clad PEEK capillaries in order to keep the flow path free of steel and to ensure pressure stability to at least 600 bar. In case of doubt, please consult common literature about the chemical compatibility of PEEK.

Titanium

Titanium is highly resistant to oxidizing acids (e.g. nitric, perchloric and hypochlorous acid) over a wide range of concentrations and temperatures, which is due to a thin oxide layer on the surface, which is stabilized by oxidizing compounds. Reducing acids (e.g. hydrochloric, sulfuric and phosphoric acid) can slightly corrode titanium, which increases with acid concentration and temperature. For example the corrosion rate of 3 % HCl (about pH 0.1) at room temperature is about 13 $\mu\text{m}/\text{year}$. At room temperature, titanium is resistant to concentrations of about 5 % sulfuric acid (about pH 0.3). The addition of nitric acid to hydrochloric or sulfuric acids significantly reduces corrosion rates. Titanium is subject to corrosion in anhydrous methanol, which can be avoided by adding a small amount of water (about 3 %). Slight corrosion is possible with ammonia > 10 %.

Fused silica

Fused silica is inert against all common solvents and acids except hydrofluoric acid. It is corroded by strong bases and should not be used above pH 12 at room temperature. The corrosion of flow cell windows can negatively affect measurement results. For a pH larger than 12, the use of flow cells with sapphire windows is recommended.

Gold

Gold is inert against all common HPLC solvents, acids and bases within the specified pH range. It can be corroded by complexing cyanides and concentrated acids like aqua regia (a mixture of concentrated hydrochloric and nitric acid).

Zirconium Oxide

Zirconium Oxide (ZrO_2) is inert against almost all common acids, bases and solvents. There are no documented incompatibilities for HPLC applications.

Platinum/Iridium

Platinum/Iridium is inert against almost all common acids, bases and solvents. There are no documented incompatibilities for HPLC applications.

PTFE

PTFE (polytetrafluorethen) is inert against almost all common acids, bases and solvents. There are no documented incompatibilities for HPLC applications.

Sapphire, Ruby and Al_2O_3 -based ceramics

Sapphire, Ruby and ceramics based on Al_2O_3 are inert against almost all common acids, bases and solvents. There are no documented incompatibilities for HPLC applications.

Data above were collected from external resources and are meant as a reference. Agilent cannot guarantee for the completeness and correctness of such information. Information can also not be generalized due to catalytic effects of impurities like metal ions, complexing agents, oxygen etc. Most data available refers to room temperature (typically 20 – 25 °C, 68 – 77 °F). If corrosion is possible it usually increases at higher temperatures. In case of doubt, additional resources should be consulted.

4 Using the Pump

Solvent Information



5 Optimizing Performance

When to Use a Degasser 56

Operational Hints for the Multi Channel Gradient Valve (MCGV) 57

When to use the Seal Wash Option 58

Choosing the Right Pump Seals 59

Optimize the Compressibility Compensation Setting 60

This chapter gives hints on how to optimize the performance or use additional devices.



When to Use a Degasser

The bio-inert quaternary pump has a built-in degasser, which should always be included to the flow path.

Operational Hints for the Internal Degasser

If you are using the degasser for the first time, if the degasser was switched off for some time (for example, overnight), or if the degasser lines are empty, you should prime the degasser before running an analysis.

The degasser can be primed either by drawing solvent through the degasser with a syringe or by pumping with the pump.

Priming the degasser with a syringe is recommended, when:

- degasser is used for the first time, or solvent tubes are empty, or
- changing to solvents that are immiscible with the solvent currently in the vacuum tubes.

Priming the degasser by using the pump at high flow rate (3 – 5 mL/min) is recommended, when:

- pump was turned off for some time (for example, during night) and volatile solvent mixtures are used, or
- solvents have been changed.

Operational Hints for the Multi Channel Gradient Valve (MCGV)

In a mixture of salt solutions and organic solvent the salt solution might be well dissolved in the organic solvent without showing precipitations. However in the mixing point of the gradient valve, at the boundary between the two solvents, micro precipitation is possible. Gravity forces the salt particles to fall down. Normally the A channel of the valve is used for the aqueous/salt solution and the B channel of the pump is used for the organic solvent. If used in this configuration the salt will fall back into the aqueous solution and will be dissolved. When using the pump in a different configuration (e.g., D - salt solution, A - organic solvent) the salt can fall into the port of the organic solvent and may lead to performance problems.

NOTE

When using salt solutions and organic solvents it is recommended to connect the salt solution to one of the bottom ports of the MCGV and the organic solvent to one of the upper gradient valve ports. It is best to have the organic channel directly above the salt solution channel. Regular flushing with water of all MCGV channels is recommended to remove all possible salt deposits in the valve ports.

When to use the Seal Wash Option

Highly concentrated buffer solutions will reduce the lifetime of the seals and pistons in your pump. The seal wash option allows to maintain the seal lifetime by flushing the back side of the seal with a wash solvent.

The seal wash option is strongly recommended when buffer concentrations of 0.1 M or higher will be used for long time periods in the pump.

The seal wash option comprises a support ring, secondary seal, gasket and seal holder for both piston sides. Place a wash bottle filled with 90 % water / 10 % isopropanol above the pump in the solvent cabinet. The peristaltic pump moves a flow through the pump head removing all possible buffer crystals from the back of the pump seal. This mixture prevents growth of algae or bacteria in the wash bottle and reduces the surface tension of the water.

Choosing the Right Pump Seals

The standard seal for the pump can be used for most applications. However applications that use normal phase solvents (for example, hexane) are not suited for the standard seal and require a different seal when used for a longer time in the pump.

For applications that use normal phase solvents (for example, hexane) we recommend using bio-inert PE piston seals and bio-inert PE wash seals. For normal phase applications, these seals have less abrasion compared to the standard seals.

NOTE

Polyethylene seals have a limited pressure range of 0 – 200 bar. When used above 200 bar their lifetime is reduced significantly. *DO NOT* apply the seal wear-in procedure performed with new standard seals at 600 bar.

Optimize the Compressibility Compensation Setting

The compressibility compensation default setting is 100×10^{-6} /bar for the pump. This setting represents an average value. Under normal conditions the default setting typically reduces the pressure pulsation to values below 1 % of system pressure that are sufficient for most applications and for all gradient analyses. The compressibility settings can be optimized by using the values for the various solvents described in [Table 5](#) on page 61. If the solvent in use is not listed in the compressibility tables, when using isocratic mixtures of solvents and if the default settings are not sufficient for your application the following procedure can be used to optimize the compressibility settings.

NOTE

When using mixtures of solvents it is not possible to calculate the compressibility of the mixture by interpolating the compressibility values of the pure solvents used in that mixture or by applying any other calculation. In these cases the following empirical procedure has to be applied to optimize your compressibility setting.

Inappropriate settings would mainly affect retention times of peaks eluted at the beginning of a gradient. Therefore optimize settings for the solvent at the beginning of the gradient. For mixtures containing up to 50 % water, use compressibility settings of water.

- 1 Start the pump with the required flow rate.
- 2 Before starting the optimization procedure, the flow must be stable. Check the tightness of the system with the pressure test.
- 3 Your pump must be connected to a data system or Instant Pilot with which the pressure and %-ripple can be monitored, or connect an external measurement device to the analog pressure output (see [“Electrical Connections”](#) on page 168)
- 4 Starting with a compressibility setting of 10×10^{-6} /bar increase the value in steps of 10. Re-zero the signal display as required. The compressibility compensation setting that generates the smallest pressure ripple is the optimum value for your solvent composition. If ChemStation (classic view) is used and the ripple shown is positive, then the compressibility setting should be decreased. If it is negative, it should be increased.

Table 5 Solvent Compressibility

Solvent (pure)	Compressibility (10-6/bar)
Acetone	126
Acetonitrile	115
Benzene	95
Carbon tetrachloride	110
Chloroform	100
Cyclohexane	118
Ethanol	114
Ethyl acetate	104
Heptane	120
Hexane	150
Isobutanol	100
Isopropanol	100
Methanol	120
1-Propanol	100
Toluene	87
Water	46

5 Optimizing Performance

Optimize the Compressibility Compensation Setting



6 Troubleshooting and Diagnostics

Overview of the Module's Indicators and Test Functions	64
Status Indicators	66
Power Supply Indicator	66
Module Status Indicator	67
User Interfaces	68
Agilent Lab Advisor Software	69

This chapter gives an overview about the troubleshooting and diagnostic features and the different user interfaces.



Overview of the Module's Indicators and Test Functions

Status Indicators

The module is provided with two status indicators which indicate the operational state (prerun, run, and error states) of the module. The status indicators provide a quick visual check of the operation of the module.

Error Messages

In the event of an electronic, mechanical or hydraulic failure, the module generates an error message in the user interface. For each message, a short description of the failure, a list of probable causes of the problem, and a list of suggested actions to fix the problem are provided (see chapter Error Information).

Test Functions

A series of test functions are available for troubleshooting and operational verification after exchanging internal components (see Tests and Calibrations).

System Pressure Test

The **System Pressure Test** is a quick test designed to determine the pressure tightness of the system (i.e. the high pressure flow path between pump and column). After exchanging flow path components (e.g. pump seals or injection seal), use this test to verify the system is pressure tight, see “[System Pressure Test](#)” on page 97.

Leak Rate Test

The **Leak Rate Test** is a diagnostic test designed to determine the pressure tightness of the pump components. When a problem with the pump is suspected, use this test to help troubleshoot the pump and its pumping performance, see “[Leak Rate Test](#)” on page 102.

Status Indicators

Two status indicators are located on the front of the module. The lower left indicates the power supply status, the upper right indicates the module status.

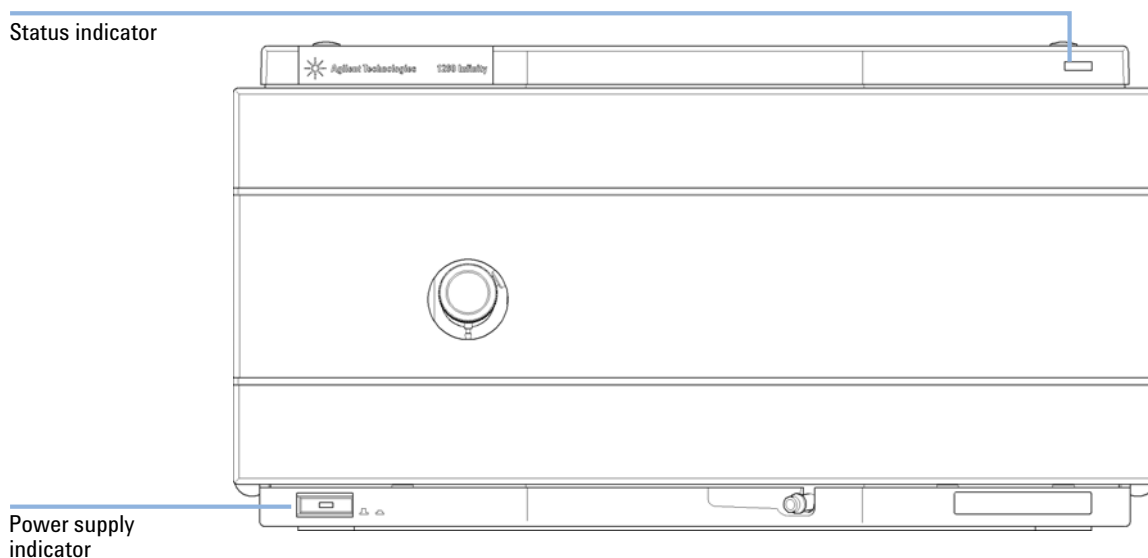


Figure 9 Location of Status Indicators

Power Supply Indicator

The power supply indicator is integrated into the main power switch. When the indicator is illuminated (*green*) the power is *ON*.

Module Status Indicator

The module status indicator indicates one of six possible module conditions:

- When the status indicator is *OFF* (and power switch light is on), the module is in a *prerun* condition, and is ready to begin an analysis.
- A *green* status indicator, indicates the module is performing an analysis (*run mode*).
- A *yellow* indicator indicates a *not-ready* condition. The module is in a not-ready state when it is waiting for a specific condition to be reached or completed (for example, immediately after changing a set point), or while a self-test procedure is running.
- An *error* condition is indicated when the status indicator is *red*. An error condition indicates the module has detected an internal problem which affects correct operation of the module. Usually, an error condition requires attention (e.g. leak, defective internal components). An error condition always interrupts the analysis.
- A *red-blinking* (modules with on-board LAN) or *yellow-blinking* (modules without on-board LAN) indicator indicates that the module is in resident mode (e.g. during update of main firmware).
- A *fast red-blinking* (modules with on-board LAN) or *fast yellow-blinking* (modules without on-board LAN) indicator indicates that the module is in boot loader mode (e.g. during update of main firmware). In such a case try to re-boot the module or try a cold-start.

User Interfaces

Depending on the user interface, the available tests vary. Some descriptions are only available in the service manual.

Table 6 Test functions available vs. user interface

Test	Instant Pilot G4208A	Agilent Lab Advisor
System Pressure Test	Yes (B.02.11)	Yes (B.01.04)
Leak Rate Test	No	Yes (B.01.04. SP1)

Agilent Lab Advisor Software

The Agilent Lab Advisor software is a standalone product that can be used with or without data system. Agilent Lab Advisor software helps to manage the lab for high quality chromatographic results and can monitor in real time a single Agilent LC or all the Agilent GCs and LCs configured on the lab intranet.

Agilent Lab Advisor software provides diagnostic capabilities for all Agilent 1200 Infinity Series modules. This includes diagnostic capabilities, calibration procedures and maintenance routines for all the maintenance routines.

The Agilent Lab Advisor software also allows users to monitor the status of their LC instruments. The Early Maintenance Feedback (EMF) feature helps to carry out preventive maintenance. In addition, users can generate a status report for each individual LC instrument. The tests and diagnostic features as provided by the Agilent Lab Advisor software may differ from the descriptions in this manual. For details refer to the Agilent Lab Advisor software help files.

This manual provides lists with the names of Error Messages, Not Ready messages, and other common issues.

The Instrument Utilities is a basic version of the LabAdvisor with limited functionality required for installation, use and maintenance. No advanced repair, troubleshooting and monitoring functionality is included.

6 Troubleshooting and Diagnostics
Agilent Lab Advisor Software



7 Error Information

What Are Error Messages	73
General Error Messages	74
Timeout	74
Shut-Down	74
Remote Timeout	75
Synchronization Lost	76
Leak	76
Leak Sensor Open	77
Leak Sensor Short	77
Compensation Sensor Open	78
Compensation Sensor Short	78
Fan Failed	79
Open Cover	79
Module Error Messages	80
Restart Without Cover	80
Solvent Zero Counter	80
Pressure Above Upper Limit	81
Pressure Below Lower Limit	82
Pressure Signal Missing	82
Missing Pressure Reading	83
Pump Configuration	83
Valve Fuse	84
Inlet-Valve Fuse	84
Valve Failed (MCGV/SSV)	85
Motor-Drive Power	86
Encoder Missing	87
Inlet-Valve Missing	87
Temperature Out of Range	88



7 Error Information

Agilent Lab Advisor Software

Temperature Limit Exceeded	88
Servo Restart Failed	89
Pump Head Missing	89
Index Limit	90
Index Adjustment	90
Index Missing	91
Stroke Length	91
Initialization Failed	92
Wait Timeout	93
Degasser: cannot read signal	94
Degasser: limit not reached	94

This chapter describes the meaning of error messages, and provides information on probable causes and suggested actions how to recover from error conditions.

What Are Error Messages

Error messages are displayed in the user interface when an electronic, mechanical, or hydraulic (flow path) failure occurs which requires attention before the analysis can be continued (for example, repair, or exchange of consumables is necessary). In the event of such a failure, the red status indicator at the front of the module is switched on, and an entry is written into the module logbook.

General Error Messages

General error messages are generic to all Agilent series HPLC modules and may show up on other modules as well.

Timeout

The timeout threshold was exceeded.

Probable cause

- 1 The analysis was completed successfully, and the timeout function switched off the module as requested.
- 2 A not-ready condition was present during a sequence or multiple-injection run for a period longer than the timeout threshold.

Suggested actions

- Check the logbook for the occurrence and source of a not-ready condition. Restart the analysis where required.
- Check the logbook for the occurrence and source of a not-ready condition. Restart the analysis where required.

Shut-Down

An external instrument has generated a shut-down signal on the remote line.

The module continually monitors the remote input connectors for status signals. A LOW signal input on pin 4 of the remote connector generates the error message.

Probable cause

- 1 Leak detected in another module with a CAN connection to the system.
- 2 Leak detected in an external instrument with a remote connection to the system.

Suggested actions

- Fix the leak in the external instrument before restarting the module.
- Fix the leak in the external instrument before restarting the module.

Probable cause

- 3 Shut-down in an external instrument with a remote connection to the system.
- 4 The degasser failed to generate sufficient vacuum for solvent degassing.

Suggested actions

- Check external instruments for a shut-down condition.
- Check the vacuum degasser for an error condition. Refer to the *Service Manual* for the degasser or the 1260 pump that has the degasser built-in.

Remote Timeout

A not-ready condition is still present on the remote input. When an analysis is started, the system expects all not-ready conditions (for example, a not-ready condition during detector balance) to switch to run conditions within one minute of starting the analysis. If a not-ready condition is still present on the remote line after one minute the error message is generated.

Probable cause

- 1 Not-ready condition in one of the instruments connected to the remote line.
- 2 Defective remote cable.
- 3 Defective components in the instrument showing the not-ready condition.

Suggested actions

- Ensure the instrument showing the not-ready condition is installed correctly, and is set up correctly for analysis.
- Exchange the remote cable.
- Check the instrument for defects (refer to the instrument's documentation).

Synchronization Lost

During an analysis, the internal synchronization or communication between one or more of the modules in the system has failed.

The system processors continually monitor the system configuration. If one or more of the modules is no longer recognized as being connected to the system, the error message is generated.

Probable cause

- 1 CAN cable disconnected.
- 2 Defective CAN cable.
- 3 Defective main board in another module.

Suggested actions

- Ensure all the CAN cables are connected correctly.
 - Ensure all CAN cables are installed correctly.
- Exchange the CAN cable.
- Switch off the system. Restart the system, and determine which module or modules are not recognized by the system.

Leak

A leak was detected in the module.

The signals from the two temperature sensors (leak sensor and board-mounted temperature-compensation sensor) are used by the leak algorithm to determine whether a leak is present. When a leak occurs, the leak sensor is cooled by the solvent. This changes the resistance of the leak sensor which is sensed by the leak-sensor circuit on the main board.

Probable cause

- 1 Loose fittings.
- 2 Broken capillary.
- 3 Loose or leaking purge valve, inlet valve, or outlet valve.
- 4 Defective pump seals.

Suggested actions

- Ensure all fittings are tight.
- Exchange defective capillaries.
- Ensure pump components are seated correctly. If there are still signs of a leak, exchange the appropriate seal (purge valve, inlet valve, outlet valve).
- Exchange the pump seals.

Leak Sensor Open

The leak sensor in the module has failed (open circuit).

The current through the leak sensor is dependent on temperature. A leak is detected when solvent cools the leak sensor, causing the leak-sensor current to change within defined limits. If the current falls outside the lower limit, the error message is generated.

Probable cause

- 1 Leak sensor not connected to the main board.
- 2 Defective leak sensor.
- 3 Leak sensor incorrectly routed, being pinched by a metal component.

Suggested actions

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Leak Sensor Short

The leak sensor in the module has failed (short circuit).

The current through the leak sensor is dependent on temperature. A leak is detected when solvent cools the leak sensor, causing the leak-sensor current to change within defined limits. If the current increases above the upper limit, the error message is generated.

Probable cause

- 1 Defective flow sensor.
- 2 Leak sensor incorrectly routed, being pinched by a metal component.

Suggested actions

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Compensation Sensor Open

The ambient-compensation sensor (NTC) on the main board in the module has failed (open circuit).

The resistance across the temperature compensation sensor (NTC) on the main board is dependent on ambient temperature. The change in resistance is used by the leak circuit to compensate for ambient temperature changes. If the resistance across the sensor increases above the upper limit, the error message is generated.

Probable cause

- 1 Defective main board.

Suggested actions

Please contact your Agilent service representative.

Compensation Sensor Short

The ambient-compensation sensor (NTC) on the main board in the module has failed (short circuit).

The resistance across the temperature compensation sensor (NTC) on the main board is dependent on ambient temperature. The change in resistance is used by the leak circuit to compensate for ambient temperature changes. If the resistance across the sensor falls below the lower limit, the error message is generated.

Probable cause

- 1 Defective main board.

Suggested actions

Please contact your Agilent service representative.

Fan Failed

The cooling fan in the module has failed.

The hall sensor on the fan shaft is used by the main board to monitor the fan speed. If the fan speed falls below a certain limit for a certain length of time, the error message is generated.

This limit is given by 2 revolutions/second for longer than 5 seconds.

Probable cause

- 1 Fan cable disconnected.
- 2 Defective fan.
- 3 Defective main board.
- 4 Improperly positioned cables or wires obstructing fan blades.

Suggested actions

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Ensure the fan is not mechanically blocked.

Open Cover

The top foam has been removed.

The sensor on the main board detects when the top foam is in place. If the foam is removed, the fan is switched off, and the error message is generated.

Probable cause

- 1 The top foam was removed during operation.
- 2 Foam not activating the sensor.
- 3 Dirty or defective sensor.
- 4 Rear of the module is exposed to strong direct sunlight.

Suggested actions

- Reinstall the top foam.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Ensure that the rear of module is not directly exposed to strong sunlight.

Module Error Messages

These errors are pump specific.

Restart Without Cover

The module was restarted with the top cover and foam open.

The sensor on the main board detects when the top foam is in place. If the module is restarted with the foam removed, the module switches off within 30 s, and the error message is generated.

Probable cause

- 1 The module started with the top cover and foam removed.
- 2 Rear of the module is exposed to strong direct sunlight.

Suggested actions

- Please contact your Agilent service representative.
- Ensure that the rear of module is not directly exposed to strong sunlight.

Solvent Zero Counter

Pump firmware version A.02.32 and higher allow to set solvent bottle fillings in the data system. If the volume level in the bottle falls below the specified value the error message appears when the feature is configured accordingly.

Probable cause

- 1 Volume in bottle below specified volume.
- 2 Incorrect setting.

Suggested actions

- Refill bottles and reset solvent counters.
- Make sure the limits are set correctly.

Pressure Above Upper Limit

The system pressure has exceeded the upper pressure limit.

Probable cause

- 1 Upper pressure limit set too low.
- 2 Blockage in the flowpath (after the damper).
- 3 Defective damper.
- 4 Defective main board.

Suggested actions

- Ensure the upper pressure limit is set to a value suitable for the analysis.
- Check for blockage in the flowpath. The following components are particularly subject to blockage: inline filter frit, needle (autosampler), seat capillary (autosampler), sample loop (autosampler), column frits and capillaries with small internal diameters (e.g. 50 µm ID).
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Pressure Below Lower Limit

The system pressure has fallen below the lower pressure limit.

Probable cause

- 1 Lower pressure limit set too high.
- 2 Air bubbles in the mobile phase.
- 3 Leak.
- 4 Defective damper.
- 5 Defective main board.
- 6 Defective active inlet valve (AIV)

Suggested actions

- Ensure the lower pressure limit is set to a value suitable for the analysis.
- Ensure solvents are degassed. Purge the module.
 - Ensure solvent inlet filters are not blocked.
 - Inspect the pump head, capillaries and fittings for signs of a leak.
 - Purge the module. Run a pressure test to determine whether the seals or other module components are defective.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Replace AIV cartridge.

Pressure Signal Missing

The pressure signal of the damper is missing.

The pressure signal of the damper must be within a specific voltage range. If the pressure signal is missing, the processor detects a voltage of approximately -120mV across the damper connector.

Probable cause

- 1 Damper disconnected.
- 2 Defective damper.

Suggested actions

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Missing Pressure Reading

The pressure readings read by the pump ADC (analog-digital converter) are missing.

The ADC reads the pressure signal of from the damper every 1ms. If the readings are missing for longer than 10 seconds, the error message is generated.

Probable cause

- 1 Damper disconnected.
- 2 Defective damper.
- 3 Defective main board.

Suggested actions

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Pump Configuration

At switch-on, the quaternary pump has recognized a new pump configuration.

The quaternary pump is assigned its configuration at the factory. If the gradient valve is disconnected, and the quaternary pump is rebooted, the error message is generated. However, the pump will function as an isocratic pump in this configuration. The error message reappears after each switch-on.

Probable cause

- 1 Gradient valve disconnected.

Suggested actions

- Reconnect the gradient valve.

Valve Fuse

Valve Fuse 0: Channels A and B

Valve Fuse 1: Channels C and D

The gradient valve in the quaternary pump has drawn excessive current causing the electronic fuse to open.

Probable cause

- 1 Defective gradient valve.
- 2 Defective connection cable (front panel to main board).
- 3 Defective main board.

Suggested actions

- Restart the quaternary pump. If the error message appears again, exchange the gradient valve.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Inlet-Valve Fuse

The active-inlet valve in the module has drawn excessive current causing the inlet-valve electronic fuse to open.

Probable cause

- 1 Defective active inlet valve.
- 2 Defective connection cable (front panel to main board).
- 3 Defective main board.

Suggested actions

- Restart the module. If the error message appears again, exchange the active inlet valve.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Valve Failed (MCGV/SSV)

Valve 0 Failed: valve A

Valve 1 Failed: valve B

Valve 2 Failed: valve C

Valve 3 Failed: valve D

One of the valves of the multi-channel gradient valve has failed to switch correctly.

The processor monitors the valve voltage before and after each switching cycle. If the voltages are outside expected limits, the error message is generated.

Probable cause

- 1 Gradient valve disconnected.
- 2 Connection cable (inside instrument) not connected.
- 3 Connection cable (inside instrument) defective.
- 4 Gradient valve defective.

Suggested actions

- Ensure the gradient valve is connected correctly.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Exchange the gradient valve.

Motor-Drive Power

The current drawn by the pump motor exceeded the maximum limit.

Blockages in the flow path are usually detected by the pressure sensor in the damper, which result in the pump switching off when the upper pressure limit is exceeded. If a blockage occurs before the damper, the pressure increase cannot be detected by the pressure sensor and the module will continue to pump. As pressure increases, the pump drive draws more current. When the current reaches the maximum limit, the module is switched off, and the error message is generated.

Probable cause

- 1** Flow path blockage in front of the damper.
- 2** Blocked active inlet valve.
- 3** Blocked outlet valve.
- 4** High friction (partial mechanical blockage) in the pump drive assembly.
- 5** Defective pump drive assembly.
- 6** Defective main board.

Suggested actions

- Ensure the capillaries and frits between the pump head and damper inlet are free from blockage.
- Exchange the active inlet valve.
- Exchange the outlet valve.
- Remove the pump-head assembly. Ensure there is no mechanical blockage of the pump-head assembly or pump drive assembly.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Encoder Missing

The optical encoder on the pump motor in the module is missing or defective.

The processor checks the presence of the pump encoder connector every 2 seconds. If the connector is not detected by the processor, the error message is generated.

Probable cause

- 1 Defective or disconnected pump encoder connector.
- 2 Defective pump drive assembly.

Suggested actions

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Inlet-Valve Missing

The active-inlet valve in the module is missing or defective.

The processor checks the presence of the active-inlet valve connector every 2 seconds. If the connector is not detected by the processor, the error message is generated.

Probable cause

- 1 Disconnected or defective cable.
- 2 Disconnected or defective connection cable (front panel to main board).
- 3 Defective active inlet valve.

Suggested actions

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.
- Exchange the active inlet valve.

Temperature Out of Range

The temperature sensor readings in the motor-drive circuit are out of range.

The values supplied to the ADC by the hybrid sensors must be between 0.5 V and 4.3 V. If the values are outside this range, the error message is generated.

Probable cause

- 1 Defective main board.

Suggested actions

Please contact your Agilent service representative.

Temperature Limit Exceeded

The temperature of one of the motor-drive circuits is too high.

The processor continually monitors the temperature of the drive circuits on the main board. If excessive current is being drawn for long periods, the temperature of the circuits increases. If the temperature exceeds the upper limit, the error message is generated.

Probable cause

- 1 High friction (partial mechanical blockage) in the pump drive assembly.
- 2 Partial blockage of the flowpath in front of the damper.
- 3 Defective pump drive assembly.
- 4 Defective main board.

Suggested actions

Remove the pump-head assembly. Ensure there is no mechanical blockage of the pump-head assembly or pump drive assembly.

Ensure the outlet valve is not blocked.

Please contact your Agilent service representative.

Please contact your Agilent service representative.

Servo Restart Failed

The pump motor in the module was unable to move into the correct position for restarting.

When the module is switched on, the first step is to switch on the C phase of the variable reluctance motor. The rotor should move to one of the C positions. The C position is required for the servo to be able to take control of the phase sequencing with the commutator. If the rotor is unable to move, or if the C position cannot be reached, the error message is generated.

Probable cause	Suggested actions
1 Disconnected or defective cable.	Please contact your Agilent service representative.
2 Blocked active inlet valve.	Exchange the active inlet valve.
3 Mechanical blockage of the module.	Remove the pump-head assembly. Ensure there is no mechanical blockage of the pump-head assembly or pump drive assembly.
4 Defective pump drive assembly.	Please contact your Agilent service representative.
5 Defective main board.	Please contact your Agilent service representative.

Pump Head Missing

The pump-head end stop in the pump was not found.

When the pump restarts, the metering drive moves forward to the mechanical end stop. Normally, the end stop is reached within 20 seconds, indicated by an increase in motor current. If the end point is not found within 20 seconds, the error message is generated.

Probable cause	Suggested actions
1 Pump head not installed correctly (screws not secured, or pump head not seated correctly).	Install the pump head correctly. Ensure nothing (e.g. capillary) is trapped between the pump head and body.
2 Broken piston.	Exchange the piston.

Index Limit

The time required by the piston to reach the encoder index position was too short (pump).

During initialization, the first piston is moved to the mechanical stop. After reaching the mechanical stop, the piston reverses direction until the encoder index position is reached. If the index position is reached too fast, the error message is generated.

Probable cause

- 1 Irregular or sticking drive movement.
- 2 Defective pump drive assembly.

Suggested actions

- Remove the pump head, and examine the seals, pistons, and internal components for signs of wear, contamination or damage. Exchange components as required.
- Please contact your Agilent service representative.

Index Adjustment

The encoder index position in the module is out of adjustment.

During initialization, the first piston is moved to the mechanical stop. After reaching the mechanical stop, the piston reverses direction until the encoder index position is reached. If the time to reach the index position is too long, the error message is generated.

Probable cause

- 1 Irregular or sticking drive movement.
- 2 Defective pump drive assembly.

Suggested actions

- Remove the pump head, and examine the seals, pistons, and internal components for signs of wear, contamination or damage. Exchange components as required.
- Please contact your Agilent service representative.

Index Missing

The encoder index position in the module was not found during initialization.

During initialization, the first piston is moved to the mechanical stop. After reaching the mechanical stop, the piston reverses direction until the encoder index position is reached. If the index position is not recognized within a defined time, the error message is generated.

Probable cause

- 1 Disconnected or defective encoder cable.
- 2 Defective pump drive assembly.

Suggested actions

- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Stroke Length

The distance between the lower piston position and the upper mechanical stop is out of limits (pump).

During initialization, the module monitors the drive current. If the piston reaches the upper mechanical stop position before expected, the motor current increases as the module attempts to drive the piston beyond the mechanical stop. This current increase causes the error message to be generated.

Probable cause

- 1 Defective pump drive assembly.

Suggested actions

- Please contact your Agilent service representative.

Initialization Failed

The module failed to initialize successfully within the maximum time window.

A maximum time is assigned for the complete pump-initialization cycle. If the time is exceeded before initialization is complete, the error message is generated.

Probable cause

- 1 Blocked active inlet valve.
- 2 Defective pump drive assembly.
- 3 Defective main board.

Suggested actions

- Exchange the active inlet valve.
- Please contact your Agilent service representative.
- Please contact your Agilent service representative.

Wait Timeout

When running certain tests in the diagnostics mode or other special applications, the pump must wait for the pistons to reach a specific position, or must wait for a certain pressure or flow to be reached. Each action or state must be completed within the timeout period, otherwise the error message is generated.

Possible Reasons for a Wait Timeout:

- Pressure not reached.
- Pump channel A did not reach the delivery phase.
- Pump channel B did not reach the delivery phase.
- Pump channel A did not reach the take-in phase.
- Pump channel B did not reach the take-in phase.
- Solvent volume not delivered within the specified time.

Probable cause

- 1** Purge valve open.
- 2** Leak at fittings, purge valve, active inlet valve, outlet valve or piston seals.
- 3** Flow changed after starting test.
- 4** Defective pump drive assembly.

Suggested actions

- Ensure that purge valve is closed.
- Ensure pump components are seated correctly. If there are still signs of a leak, exchange the appropriate seal (purge valve, active inlet valve, outlet valve, piston seal).
- Ensure correct operating condition for the special application in use.
- Please contact your Agilent service representative.

Degasser: cannot read signal

The pump board gets no or wrong pressure signals from the built-in degasser.

Probable cause	Suggested actions
1 Degasser board defect, missing or not connected to the pump main board.	Please contact your Agilent service representative.
2 Degasser sensor defect or not connected to degasser board	Please contact your Agilent service representative.

Degasser: limit not reached

This error is thrown, if the degasser does not become ready after 8 min, i.e. is higher than 180 mbar.

Probable cause	Suggested actions
1 Liquid in degasser tubing.	Please contact your Agilent service representative.
2 Leak in degasser tubing or chamber.	Please contact your Agilent service representative.
3 Degasser vacuum pump defect.	Please contact your Agilent service representative.



8 Test Functions and Calibration

Introduction	96
System Pressure Test	97
Running the Test	99
Evaluating the Results	99
System Pressure Test failed	101
Leak Rate Test	102
Running the Test	104
Evaluating the Results	104
Potential Causes of Leak Rate Test Failure	105

This chapter describes the tests for the module.



Introduction

All tests are described based on the Agilent Lab Advisor Software B.01.04. Other user interfaces may not provide all tests listed here.

For details on the use of the interface refer to the interface documentation.

Table 7 Interfaces and available test functions

Interface	Comment	Available Function
Agilent Instrument Utilities		System Pressure Test
Agilent Lab Advisor	All tests are available	<ul style="list-style-type: none">• System Pressure Test• Leak Rate Test
Agilent ChemStation	No tests available	
Agilent Instant Pilot		System Pressure Test

System Pressure Test

Introduction

The **System Pressure Test** is used for checking the tightness of the LC system and identifying leaks between the pump and a position in the flow path following the pump blocked by a blank nut.


Test Principle

A solvent can be chosen from available solvent channels and a maximum pressure can be defined at which the test will be run. In contrast to older revisions of this test, any solvent can be used.

Before the test, the pump and system are flushed with solvent in order to remove air bubbles, as air bubbles are compressed during the test and therefore would appear as leaks. Using a degasser is highly recommended. Then the flow path is blocked by a blank nut at any position between the purge valve and the TCC outlet.

8 Test Functions and Calibration

System Pressure Test

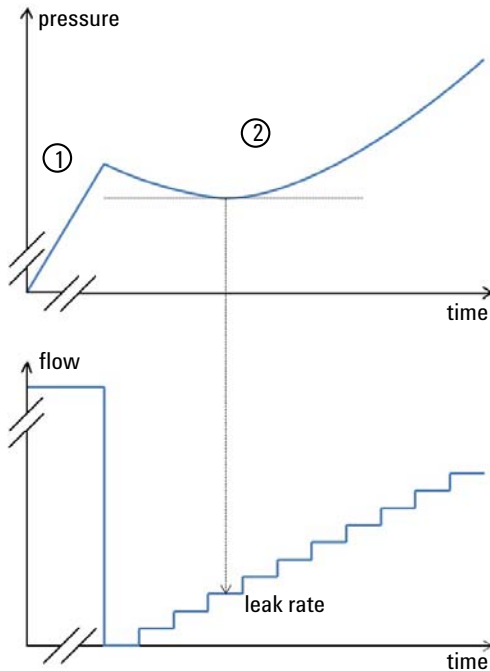

In the first phase of the test, the pump delivers flow at a rate of 200 μL until a pressure of 50 bar below the defined maximum pressure is reached. In the second phase, the pump delivers a small flow which is increased stepwise. If there is a leak in the system, the pressure will drop initially, as the low flow cannot compensate the leak flow. As soon as the pump flow rate exceeds the leak flow rate, the pressure will increase again and the test is stopped at about 20 bar below the maximum pressure. The point in phase 2, where the lowest pressure is reached and stays constant for a short time corresponds to the leak rate, that is provided as a test result. A leak rate smaller than 3 $\mu\text{L}/\text{min}$ is good enough for operating the pump reliably.

Running the Test

Parts required	p/n	Description
	01080-83202	Blank nut

CAUTION

Damage to pressure sensitive parts

Even columns that are suitable for high pressures are sensitive to pressure drops that occur during this test.

- Do not include any pressure sensitive parts to the flow path and choose a maximum pressure that is compatible to your system. For example, do not include columns, a standard pressure flow cell (up to 20 bar) or a 400 bar autosampler to a 600 bar pressure test.

Running the test from the Agilent Lab Advisor

- 1 Select the **System Pressure Test** from the **Test Selection** menu.
- 2 Start the test and follow the instructions.

NOTE

Make sure to release the pressure by slowly opening the purge valve when the test has finished.

“[Evaluating the Results](#)” on page 99 describes the evaluation and interpretation of the **System Pressure Test** results. For detailed instructions refer to the Agilent Lab Advisor software.

Evaluating the Results

The test fails, if the leak rate between pump and blank nut is higher than the limit of 3 $\mu\text{L}/\text{min}$.

If the **System Pressure Test** fails:

- Ensure that all fittings between the pump and the blank nut are tight.
- Repeat the test.

8 Test Functions and Calibration

System Pressure Test

NOTE

Often it is only a damaged blank nut itself (poorly shaped from overtightening) that causes a failure of the test. Before investigating on any other possible sources of failure make sure that the blank nut you are using is in good condition and properly tightened!

If the test fails again, insert the blank nut at the outlet of the previous module in the stack (e.g. autosampler outlet if TCC has been tested before), and repeat the test. Exclude each module one by one to determine which module is leaking.

If the pump is determined to be the source of the leak, run the **Pump Leak Rate Test**.

System Pressure Test failed

The test will fail, if the sum of all leaks in the system (pump, autosampler or column compartment and connections) exceeds the test limit. After isolating and fixing the cause of the leak, repeat the **System Pressure Test** to confirm the system is pressure tight.

Probable cause	Suggested actions
1 Purge valve open.	Close the purge valve.
2 Loose or leaky fittings.	Tighten the fitting or exchange the capillary.
3 Pump: Damaged pump seals or pistons.	Run the Leak Rate Test to confirm the leak.
4 Loose purge valve.	Tighten the purge valve nut (14 mm wrench).
5 Autosampler: Loose or leaky fitting.	Tighten or exchange the fitting or capillary.
6 Autosampler: Rotor seal (injection valve).	Exchange the rotor seal.
7 Autosampler: Damaged metering seal or piston.	Exchange the metering seal. Check the piston for scratches. Exchange the piston if required.
8 Autosampler: Needle seat.	Exchange the needle seat.
9 Column compartment: Loose or leaky fitting.	Tighten or exchange the fitting or capillary.
10 Column compartment: Rotor seal in optional valve.	Exchange the rotor seal.

Leak Rate Test

Introduction

The **Leak Rate Test** is used for verifying the internal tightness of the pump and helps identifying parts which may have caused a leak.

System requirements

Minimum software revisions:

- Lab Advisor B.01.04. SP1 (G1310B 1260 Isocratic Pump, G1311B 1260 Quaternary Pump, G4280B 1220 Isocratic Pump, G4281B 1220 Gradient Pump)
- Lab Advisor B.01.04. SP2 (G1311C 1260 Quaternary Pump VL, G4220A 1290 Binary Pump, G4220B 1290 Binary Pump VL, G5611A 1260 Bio-inert Quaternary Pump)

Minimum firmware revision: A.06.34 for G5611A and A.06.33 for all other pumps.


NOTE

This test does not work in emulation mode. In case of an emulated module, convert to the original type first.

Test Principle

A solvent can be chosen from available solvent channels and a maximum target pressure can be defined at which the test will be run. Typically, this is the maximum pressure specified for the pump. The test can be run with any solvent compatible to the pump.

Before the test, the pump is flushed with solvent in order to remove air bubbles, as air bubbles are compressed during the test and therefore would appear as leaks. Using a degasser is highly recommended.

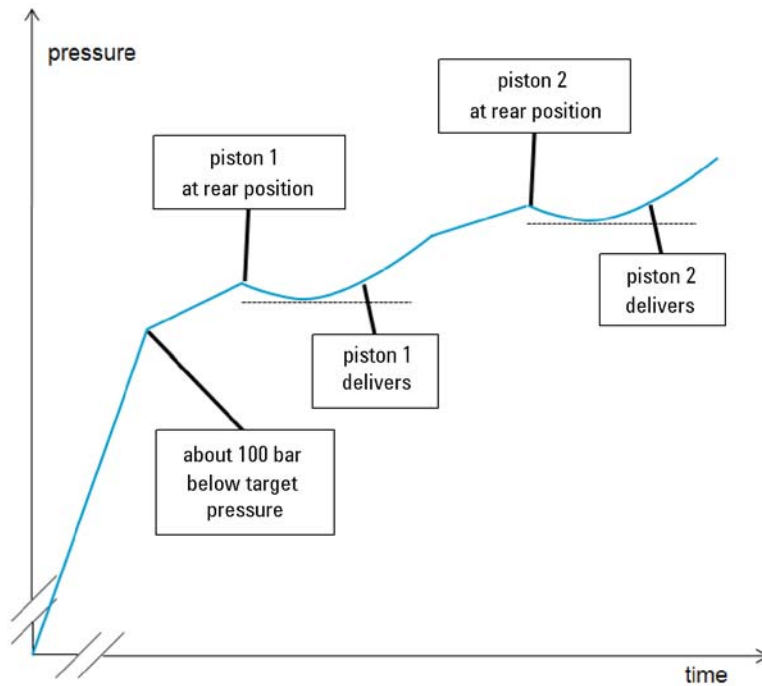

Initially, the pressure is increased to about 100 bar below the target pressure, which has been set for the test.

Then piston 1 is brought to its rear position. An increasing flow is delivered by piston 1. In case of a leak, the pressure will drop initially as long as the flow rate delivered by the piston is lower than the leak rate. As soon as the flow rate of the piston exceeds the leak rate, the measured pressure will increase again. Therefore the minimum pressure of that curve segment corresponds to the flow and leak rate at that time and the leak rate is measured. Compare to the description of the system pressure test “[System Pressure Test](#)” on page 97.

Subsequently, piston 2 is moved to its rear position, then piston 2 delivers and the measurement is done as described for piston 1.

Running the Test

Parts required	p/n	Description
	01080-83202	Blank nut

Running the test from the Agilent Lab Advisor


- 1 Select the **Leak Rate Test** from the **Test Selection** menu.
- 2 Start the test and follow the instructions.

NOTE

Make sure to release the pressure by slowly opening the purge valve when the test has finished.

Evaluating the Results

Results of the leak rate test are the leak rates measured for pistons 1 and 2 as described for the test principle. If any of the leak rates exceeds $3 \mu\text{L}/\text{min}$, the test will fail.

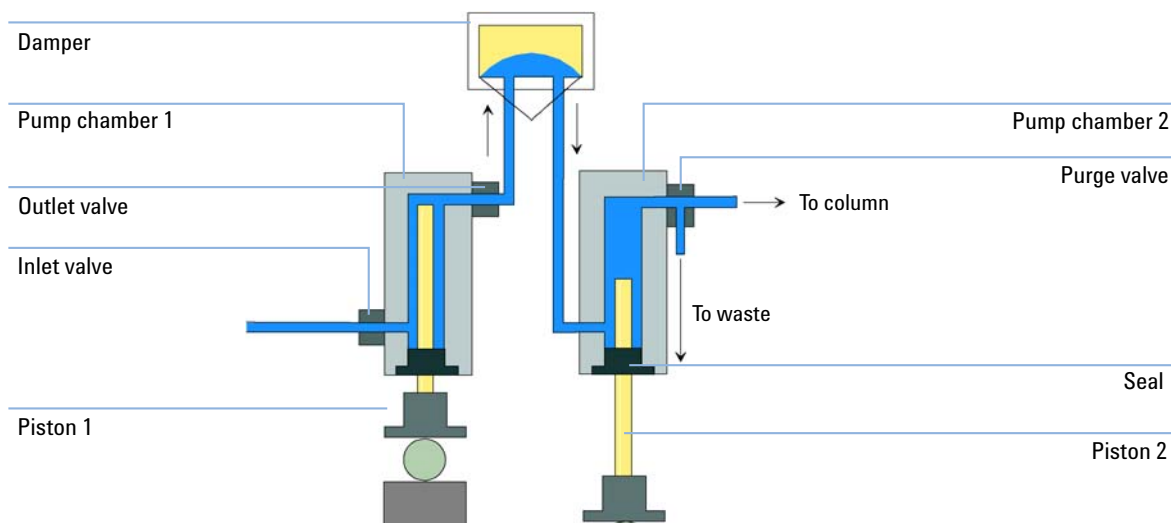

Potential Causes of Leak Rate Test Failure

Secondary Leak

If a leak is found for movement of piston 2 (secondary leak), the following reasons are possible:

Probable cause	Suggested actions
1 System not flushed properly	Flush system for several minutes
2 Degassing efficiency is low	Check degasser performance
3 Purge valve not closed or defect	Check purge valve
4 Blank nut not installed tightly	Tighten or replace blank nut
5 Outlet valve leaking (read below)	Replace outlet valve
6 Leak at piston 2 or seal in chamber 2	Inspect piston, replace piston and/or seal

Primary Leak

If a leak is found for movement of piston 1 (primary leak), any leak described for piston movement 2 will cause a failure for piston 1 as well, as the liquid can move through the outlet valve to chamber 2. Such cases need to be identified as described before. Additionally, following causes are possible:

Probable cause	Suggested actions
1 Leak at piston 1 or seal in chamber 1	Inspect piston, replace piston and/or seal
2 Leak at inlet valve	Replace inlet valve or inlet valve cartridge (AIV only)


Internal Outlet Valve Leak

A leak of the outlet valve will be identified separately (internal outlet valve leak) by calculating the difference between leak rate 1 and leak rate 2. If the second leak rate is higher than the first one, this is due to a flow back through the outlet valve.

Probable cause	Suggested actions
1 Leak at outlet valve	Replace the part which has failed and re-run the test.

8 Test Functions and Calibration

Leak Rate Test

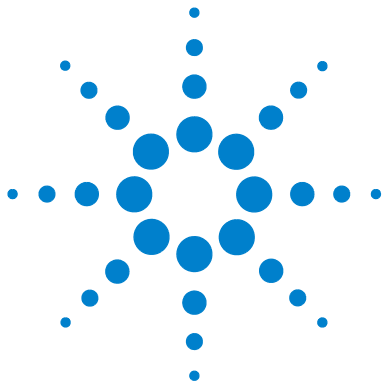

9 Maintenance

Introduction to Maintenance	108
Warnings and Cautions	109
Overview of Maintenance	111
Cleaning the Module	112
Checking and Replacing the Solvent Filter	113
Exchanging the Active Inlet Valve (AIV) or its Cartridge	114
Exchanging the Outlet Valve	117
Exchanging the Purge Valve Frit or the Purge Valve	119
Removing the Pump Head Assembly	122
Seal Wear-in Procedure	125
Maintenance of the Pump Head	126
Reinstalling the Pump Head Assembly	129
Exchanging the Multi-Channel Gradient Valve (MCGV)	131
Exchanging the Optional Interface Board	134
Replacing the Module's Firmware	136

This chapter describes the maintenance of the module.



Introduction to Maintenance

The module is designed for easy repair. The most frequent repairs such as piston seal change and purge valve frit change can be done from the front of the module with the module in place in the system stack.

Warnings and Cautions

WARNING

Toxic, flammable and hazardous solvents, samples and reagents

The handling of solvents, samples and reagents can hold health and safety risks.

- When working with these substances observe appropriate safety procedures (for example by wearing goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet supplied by the vendor, and follow good laboratory practice.
 - The volume of substances should be reduced to the minimum required for the analysis.
 - Do not operate the instrument in an explosive atmosphere.
-

WARNING

Electrical shock

Repair work at the module can lead to personal injuries, e.g. shock hazard, when the cover is opened.

- Do not remove the metal top cover of the module. No serviceable parts inside.
 - Only certified persons are authorized to carry out repairs inside the module.
-

WARNING

Personal injury or damage to the product

Agilent is not responsible for any damages caused, in whole or in part, by improper use of the products, unauthorized alterations, adjustments or modifications to the products, failure to comply with procedures in Agilent product user guides, or use of the products in violation of applicable laws, rules or regulations.

- Use your Agilent products only in the manner described in the Agilent product user guides.
-

CAUTION

Safety standards for external equipment

- If you connect external equipment to the instrument, make sure that you only use accessory units tested and approved according to the safety standards appropriate for the type of external equipment.
-

CAUTION

Sample degradation and contamination of the instrument

Metal parts in the flow path can interact with the bio-molecules in the sample leading to sample degradation and contamination.

- For bio-inert applications, always use dedicated bio-inert parts, which can be identified by the bio-inert symbol or other markers described in this manual.
 - Do not mix bio-inert and non-inert modules or parts in a bio-inert system.
-

Overview of Maintenance

The following pages describe maintenance (simple repairs) of the pump that can be carried out without opening the main cover.

Table 8 Simple Repair Procedures

Procedure	Typical Frequency	Notes
“Checking and Replacing the Solvent Filter” on page 113	If solvent filter is blocked	Gradient performance problems, intermittent pressure fluctuations
“Exchanging the Active Inlet Valve (AIV) or its Cartridge” on page 114	If internally leaking	Pressure ripple unstable, run Leak Rate Test for verification
“Exchanging the Outlet Valve” on page 117	If internally leaking	Pressure ripple unstable, run Leak Rate Test for verification
“Exchanging the Purge Valve Frit or the Purge Valve” on page 119	If internally leaking	Solvent dripping out of waste outlet when valve closed
“Exchanging the Purge Valve Frit or the Purge Valve” on page 119	If the frit shows indication of contamination or blockage	A pressure drop of > 10 bar across the frit (at a water flow of 5 mL/min with open purge valve) indicates blockage
“Maintenance of the Pump Head” on page 126	If pump performance indicates seal wear	Leaks at lower pump head side, unstable retention times, pressure ripple unstable — run Leak Rate Test for verification
Exchanging pistons, see “Maintenance of the Pump Head” on page 126	If scratched	Seal life time shorter than usual — check pistons while changing the seals
“Exchanging the Optional Interface Board” on page 134	If defective	Error condition, indicated by red status indicator

Cleaning the Module

The module case should be kept clean. Cleaning should be done with a soft cloth slightly dampened with water or a solution of water and mild detergent. Do not use an excessively damp cloth as liquid could drip into the module.

WARNING

Liquid dripping into the electronic compartment of your module.

Liquid in the module electronics can cause shock hazard and damage the module.

- Do not use an excessively damp cloth during cleaning.
 - Drain all solvent lines before opening any fittings.
-

Checking and Replacing the Solvent Filter

A functional solvent filter is essential for a good pump performance and for protecting the LC system.

When If solvent filter is blocked.

Parts required	p/n	Description
	5041-2168	Solvent inlet filter, 20 µm pore size

See “[Bottle Head Assembly](#)” on page 148 for related parts.

CAUTION

Small particles can permanently block the capillaries and valves of the module.
 Damage of the module.

- Always filter solvents.
- Never use the module without solvent inlet filter.

NOTE

If the filter is in good condition the solvent will freely drip out of the solvent tube (hydrostatic pressure). If the solvent filter is partly blocked only very little solvent will drip out of the solvent tube.

- 1 Remove the solvent filter from the inlet filter adapter and replace it by a new one.

9 Maintenance

Exchanging the Active Inlet Valve (AIV) or its Cartridge

Exchanging the Active Inlet Valve (AIV) or its Cartridge

When If internally leaking (backflow)

Tools required

Description
Wrench, 14 mm
Pair of tweezers

Parts required

p/n	Description
G5611-60025	Active Inlet Valve (AIV, Bio-inert), without cartridge
G5611-60020	Cartridge for AIV (Bio-inert)
G1311-67304	Connecting tube, MCGV to AIV


Preparations

- Switch off pump at the main power switch and unplug the power cable.
- Use a solvent shutoff valve or lift up solvent filters in solvent reservoirs for avoiding leakages.

- 1 Remove the front cover.
- 2 Unplug the active inlet valve cable from the connector.
- 3 Disconnect the solvent inlet tube from the inlet valve (be aware that solvent may leak out of the tube due to hydrostatic flow).
- 4 Unscrew the adapter from the active inlet valve.

Exchanging the Active Inlet Valve (AIV) or its Cartridge

- Using a 14 mm wrench loosen the active inlet valve and remove the valve from the pump head.

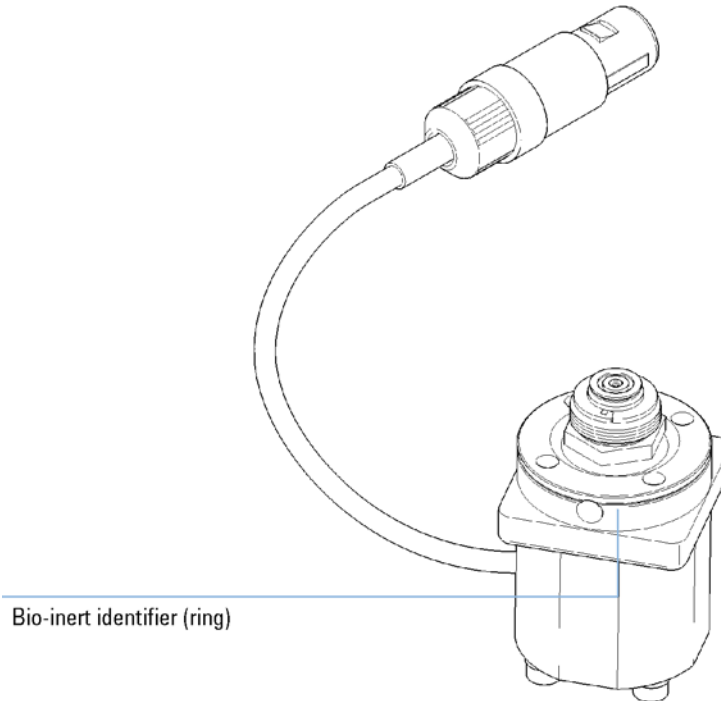

- Using a pair of tweezers remove the valve cartridge from the actuator assembly.
- Before inserting the new valve cartridge clean the area in the actuator assembly. Flush the cartridge area thoroughly with alcohol.
- Insert a new cartridge into the actuator assembly (make sure the valve cartridge is completely inserted into the actuator assembly).
- Insert the new valve into the pump head. Using the 14 mm wrench turn the nut until it is hand tight.
- Position the valve such that the solvent inlet tube connection points towards the front.
- Using the 14 mm wrench tighten the nut by turning the valve into its final position (not more than a quarter turn).
- Reconnect the adapter at the active inlet valve.

9 Maintenance

Exchanging the Active Inlet Valve (AIV) or its Cartridge

- 13 Reconnect the solvent inlet tube to the adapter. Reconnect the active inlet valve cable to the connector in the Z-panel.
- 14 Reinstall the front cover.
- 15 Purge the system with 30 mL of solvent in order to achieve a low pressure ripple, see “Regular Priming” on page 43.

Exchanging the Outlet Valve

When	If internally leaking	
Tools required	p/n	Description
	8710-0510	Wrench open 1/4 — 5/16 inch
	8710-1924	Wrench open 14 mm
Parts required	p/n	Description
	G5611-60067	Outlet Valve (Bio-inert)
Preparations	<ul style="list-style-type: none"> • Switch off pump at the main power switch • Remove the front cover 	

NOTE

Before exchanging the outlet valve you can try to clean it in an ultrasonicator for 5 – 10 min. Place the valve in an upright position in a small beaker with alcohol.


9 Maintenance

Exchanging the Outlet Valve

- 1 Using the 1/4 inch wrench disconnect the valve capillary from the outlet valve. Using the 14 mm wrench loosen the valve and remove it from the pump body.


Valve capillary

Outlet valve

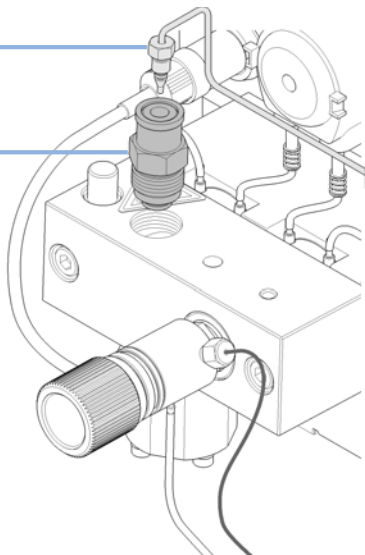

- 2 Do not disassemble the outlet valve, as this can damage the valve.

Bio-inert identifier (ring)

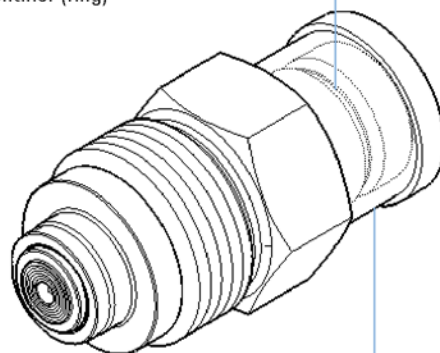


Plastic jacket

NOTE

The bio-inert outlet valve can be identified by a ring, which is normally covered by the plastic jacket.

- 3 Reinstall the outlet valve and tighten the valve. Reconnect the valve capillary.

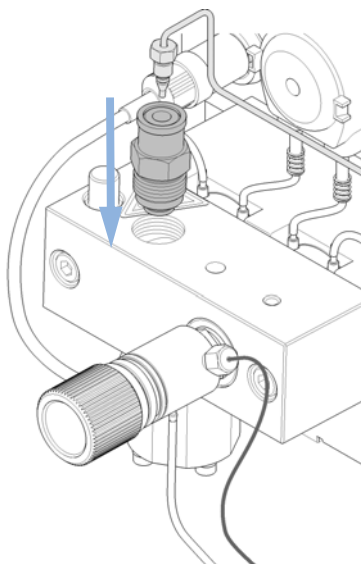

Exchanging the Purge Valve Frit or the Purge Valve

- When**
- Frit – when piston seals are exchanged or when contaminated or blocked (pressure drop of > 10 bar across the frit at a flow rate of 5 mL/min of water with purge valve opened)
 - Purge valve – if internally leaking

Tools required	p/n	Description
	8710-0510	Wrench open 1/4 — 5/16 inch
	8710-1924	Wrench open 14 mm
		Pair of tweezers
OR		Toothpick


Parts required	#	p/n	Description
	1	01018-22707	PTFE frits (pack of 5)
	1	G5611-60061	Purge Valve (Bio-inert)
	1	5067-4728	Seal cap (OPTIONAL)

- Preparations**
- Switch off pump at the main power switch
 - Remove the front cover
 - Use a solvent shutoff valve or lift up solvent filters in solvent reservoirs for avoiding leakages.


9 Maintenance

Exchanging the Purge Valve Frit or the Purge Valve

- 1** Using a 1/4 inch wrench disconnect the pump outlet capillary from the purge valve. Disconnect the waste tube. Beware of leaking solvents due to hydrostatic pressure.

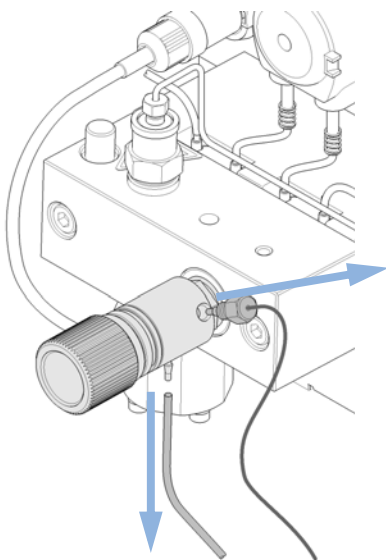


- 2** Using the 14 mm wrench unscrew the purge valve and remove it.

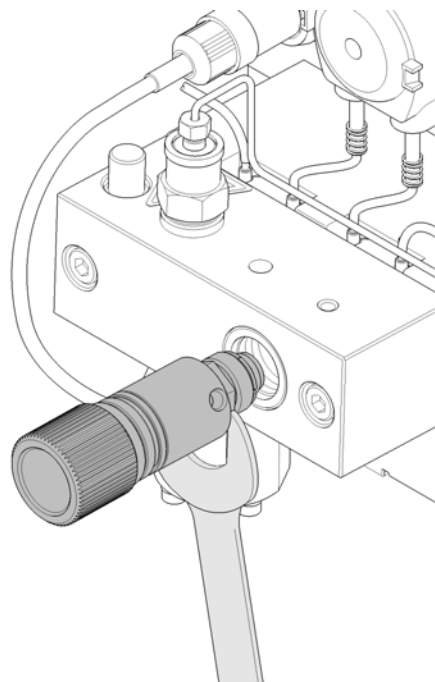


- 3** Remove the seal cap from the purge valve.

- 4** Using a pair of tweezers or a toothpick remove the frit.

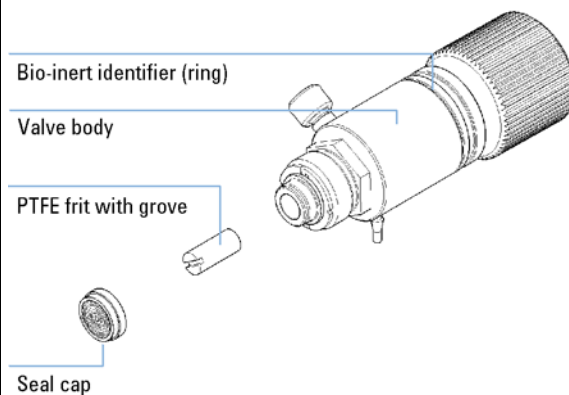


Exchanging the Purge Valve Frit or the Purge Valve

- 5** Place a new frit into the purge valve with the orientation of the frit as shown below (slit in frit points to the front). Reinstall the seal cap including the gold seal.

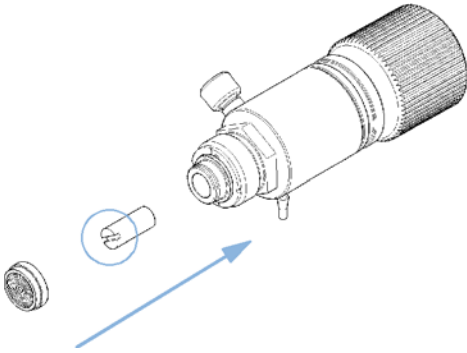**NOTE**

Before reinstallation always check the gold seal in the seal cap. A deformed seal cap should be exchanged.

- 6** Insert the purge valve into the pump head and locate the pump outlet capillary and the waste tube.

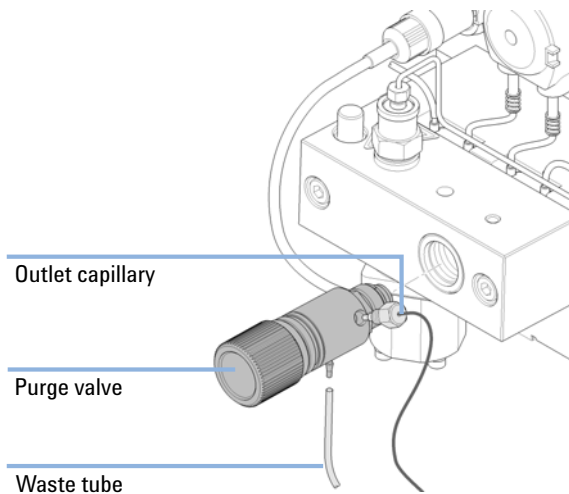

- 7** Tighten the purge valve and reconnect outlet capillary and waste tubing.

Removing the Pump Head Assembly

- When**
- Exchanging the seals
 - Exchanging the pistons
 - Exchanging seals of the seal wash option

Tools required	p/n	Description
		Wrench, 1/4 inch
	8710-2392	Hexagonal key, 4.0 mm, 15 cm long, T-handle

- Preparations**
- Switch off pump at the main power switch and unplug the power cable.
 - Use a solvent shutoff valve or lift up solvent filters in solvent reservoirs for avoiding leakages.

CAUTION


Damage of the pump drive

Starting the pump when the pump head is removed may damage the pump drive.


→ Never start the pump when the pump head is removed.

Removing the Pump Head Assembly


- 1** Remove the front cover by pressing the clip fasteners on both sides of the cover.

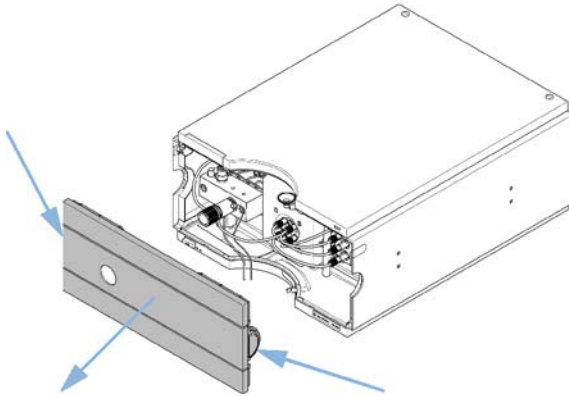

- 2** Disconnect the active inlet valve cable.

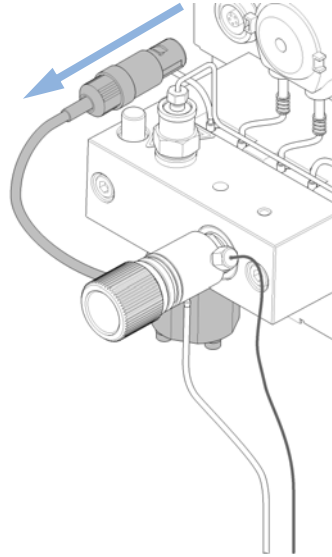

- 3** Using a 1/4 inch wrench remove the outlet capillary.

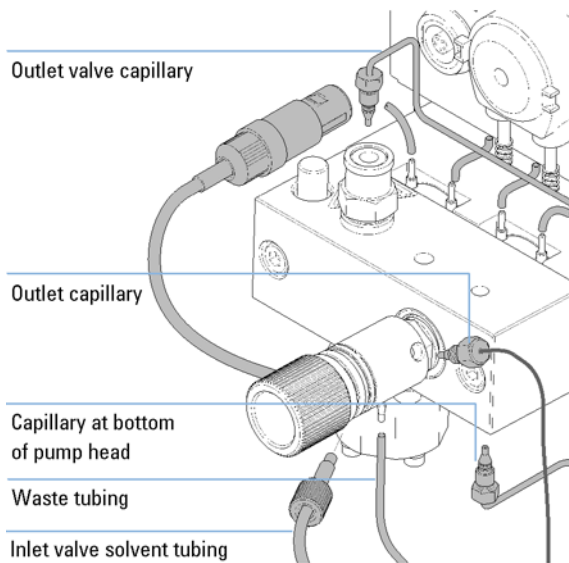


- 4** Remove the waste tubing and disconnect the solvent tubing from the inlet valve.

- 5** Remove the capillary at the bottom of the pump head.

9 Maintenance

Removing the Pump Head Assembly

- 6 Using a 4 mm hexagonal key, stepwise loosen the two pump head screws and remove the pump head from the pump drive.

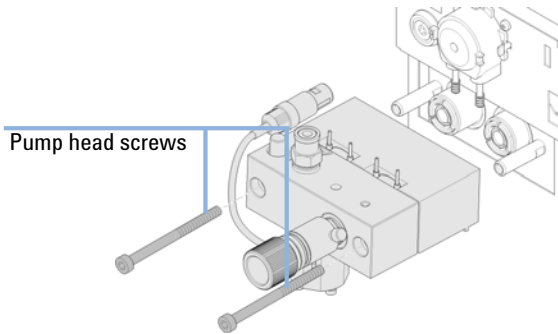

Seal Wear-in Procedure

Parts required	p/n	Description
	0100-1847	Adapter AIV to solvent inlet tubes
	5022-2159	Restriction capillary

CAUTION

Seal damage

→ This procedure is required for Standard seals (pack of 2) (p/n 5063-6589), but it will damage the PE seals.

NOTE

The stainless steel restriction capillary is acceptable for a temporary test within a bio-inert system. In case of concerns, a PEEK capillary (e.g. PEEK capillary, 150 cm length, 0.13 mm i.d. (p/n 0890-1915)) and PEEK fittings (e.g. PEEK Fittings 10/PK (p/n 5063-6591)) may be used instead.

- 1 Place a bottle with 100 mL of isopropanol in the solvent cabinet and put a tubing (including bottle head assembly) into the bottle.
- 2 If an active inlet valve is installed, screw the PEEK adapter 1/4-28 to 10-32 (p/n 0100-1847) to the AIV and connect the inlet tube from the bottle head directly to it.
- 3 Connect the Restriction capillary (p/n 5022-2159) to the purge valve. Connect its other end to a waste container.
- 4 Open the purge valve and purge the system for 5 minutes with isopropanol at a flow rate of 2 mL/min.
- 5 Close the purge valve and set the flow to a rate adequate to achieve a pressure of 350 bar. Pump 15 min at this pressure to wear in the seals. The pressure can be monitored using your instrument control software or tool.
- 6 Turn OFF the pump, slowly open the purge valve to release the pressure from the system, disconnect the restriction capillary and reinstall the bottle with the solvent for your application.
- 7 Rinse your system with the solvent used for your next application.
- 8 Replace the purge valve frit, see [“Exchanging the Purge Valve Frit or the Purge Valve”](#) on page 119.

Maintenance of the Pump Head

When When maintaining seal wash option

Tools required	p/n	Description
	8710-2392	Hexagonal key, 4.0 mm, 15 cm long, T-handle

Parts required	p/n	Description
	01018-23702	Insert tool
	0905-1731	Wash Seal PTFE for Bio-inert Pump 600 bar
	01018-07102	Gasket (Seal wash)
	G5611-21503	Piston Seal PTFE (Bio-inert)
	5067-4695	Sapphire piston

Preparations

- Switch off pump at the main power switch.
- Remove the front cover.
- Use a solvent shutoff valve or lift up solvent filters for avoiding leakages.
- Remove the pump head, see [“Removing the Pump Head Assembly”](#) on page 122.
- Remove the wash solvent tubings from the support ring inlet and outlet.


CAUTION

Seal life time


Wrongly inserting the backup ring will strongly reduce the life time of the seals and the pump head parts.

→ Note the correct installation of the backup ring inside the support ring.


1 Remove the seal holder and the seal wash support rings from the piston housing. Remove the seal holder from the support ring assembly.

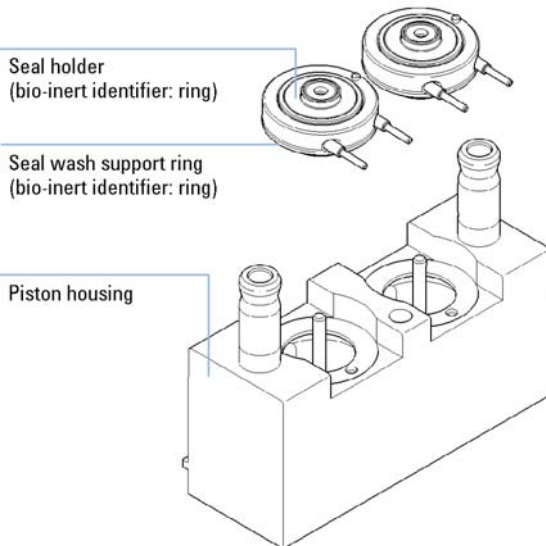


2 Using the steel side of the insert tool remove the seal wash gasket and the secondary seal from the support ring. The removed seal will be damaged and cannot be re-used!

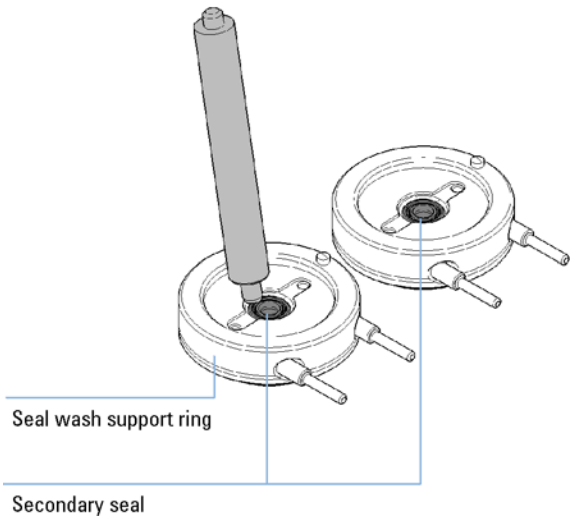

3 Using the plastic side of the insert tool press the wash seal (spring pointing upwards) into the recess of the support ring.

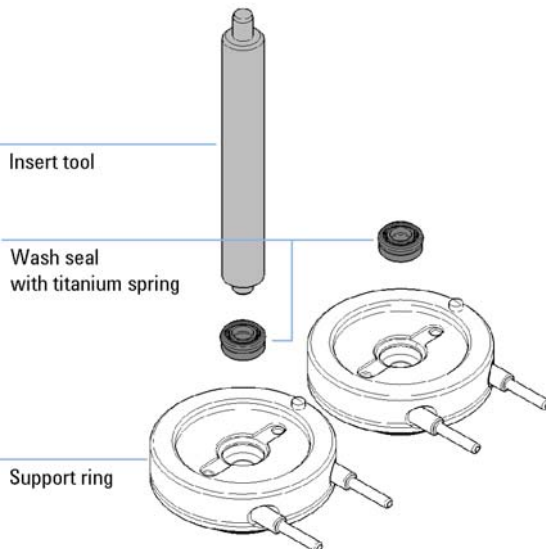


4 If the backup seal has been removed from the seal holder, insert it in the correct orientation as shown below. Please ensure that the cut end faces the seal holder.

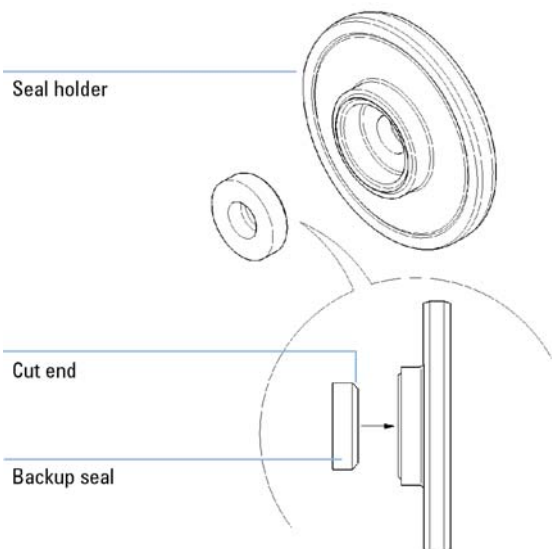


9 Maintenance

Maintenance of the Pump Head


- 5** Place a seal wash gasket in the recess of the support ring. Put the seal holder on top of the gasket.

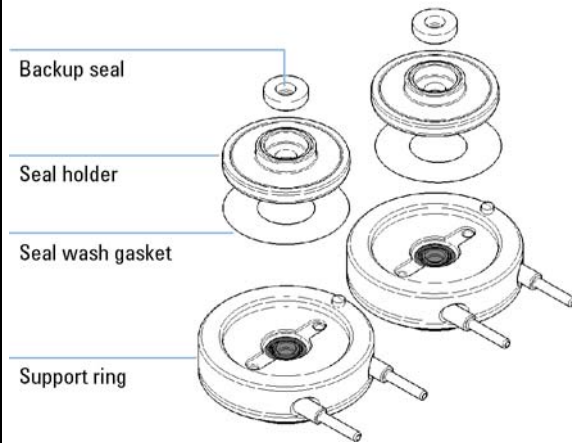


- 6** Place the support rings on the piston housing (pistons not installed) and snap the pump head and piston housing together. Note the correct position of the pin on the support ring.

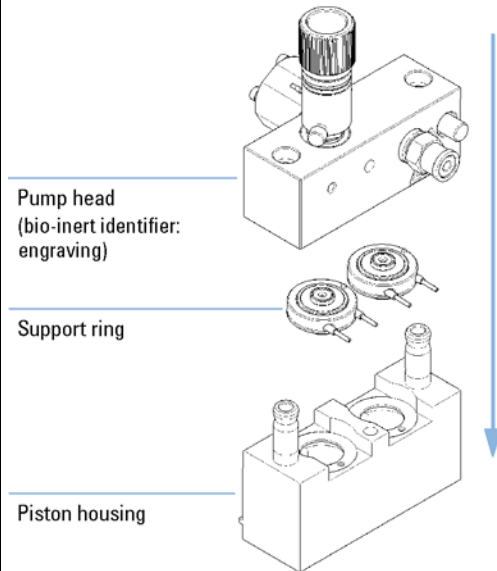

- 7** Insert the pistons and carefully press them into the seals.

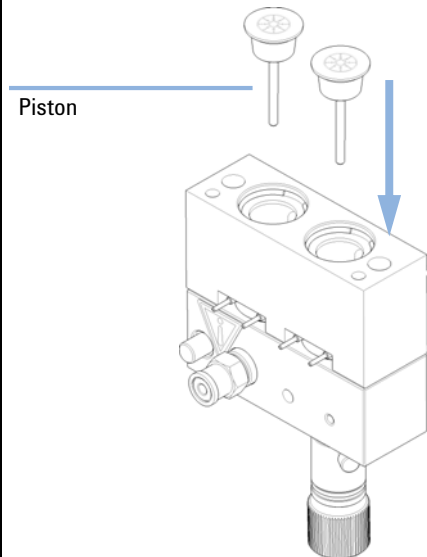

- 8** Tighten the lock screw.

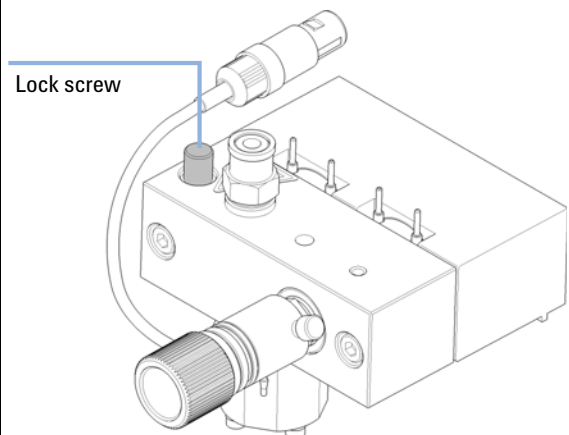

Reinstalling the Pump Head Assembly

When When reassembling the pump

Tools required


p/n	Description
8710-2392	Hexagonal key, 4.0 mm, 15 cm long, T-handle

Parts required

p/n	Description
G5611-60165	Pump Head with Seal Wash Option (Bio-inert)
79846-65501	Pump head grease

1 If needed, apply a small amount of grease on the back of the screws. Normally, the grease added during manufacturing is sufficient for a long time.


2 Slide the pump head assembly onto the pump drive and use a 4 mm hexagonal key to tighten the pump head screws stepwise with increasing torque.

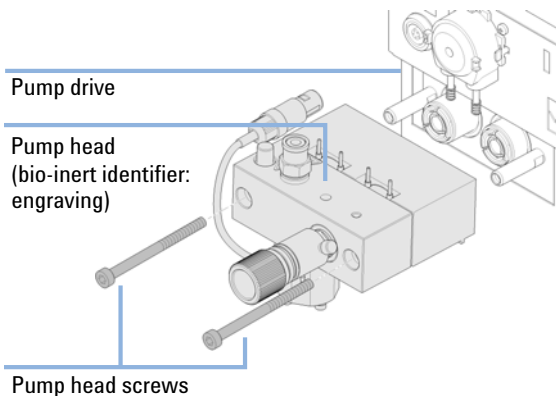


9 Maintenance

Reinstalling the Pump Head Assembly

- 3** Reconnect all capillaries, tubes and (if installed) the active inlet valve cable to its connector.

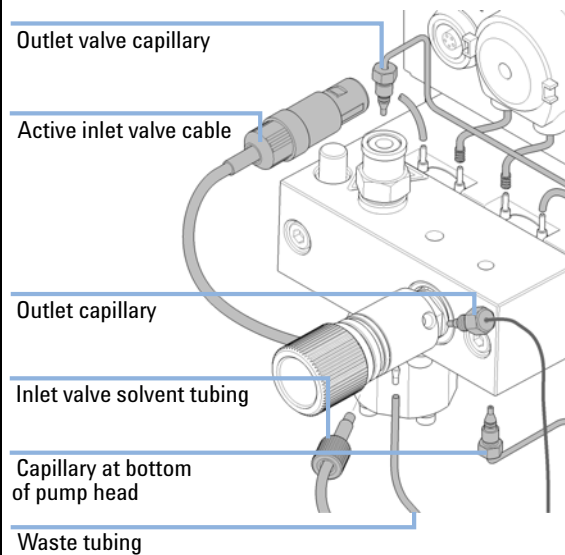

- 4** Reinstall the front cover.

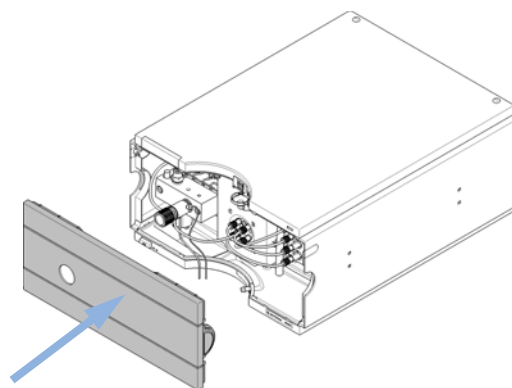

Exchanging the Multi-Channel Gradient Valve (MCGV)

Tools required	p/n	Description
	8710-0899	Screwdriver, Pozidriv #1
Parts required	p/n	Description
	G5611-67701	Multi-Channel Gradient Valve (MCGV) (Bio-inert)

- Preparations**
- Switch off pump at the main power switch
 - Remove the front cover
 - Use a solvent shutoff valve or lift up solvent filters in solvent reservoirs for avoiding leakages.


NOTE

The life time of the multi-channel gradient valve can be increased by regularly flushing the valve, especially when using buffers. If using buffers, flush all channels of the valve with water to prevent precipitation of the buffer, otherwise salt crystals could drop into an unused channel and form plugs that may cause leaks of that channel. Such leaks will interfere with the general performance of the valve. When using buffers in combination with organic solvents in the Agilent 1260 Infinity Bio-inert Quaternary Pump it is recommended to connect the aqueous solutions/buffers to one of the bottom ports (A and D) and the organic solvent to one of the upper gradient valve ports. It is best to have the organic channel directly above the buffer channel (e.g., A - buffer, B - organic solvent).


9 Maintenance

Exchanging the Multi-Channel Gradient Valve (MCGV)


- 1** Disconnect the connecting tube, waste tube and the solvent tubes from the MCGV.

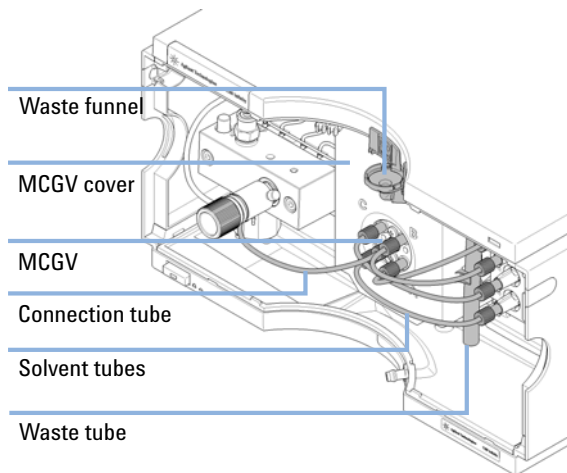

- 2** Press the lower sides of the cover to unclip it. Remove the cover.

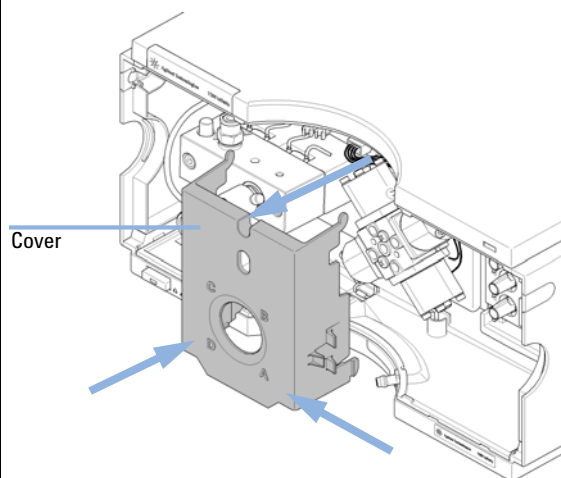

- 3** Disconnect the MCGV cable, unscrew the two screws and remove the valve.

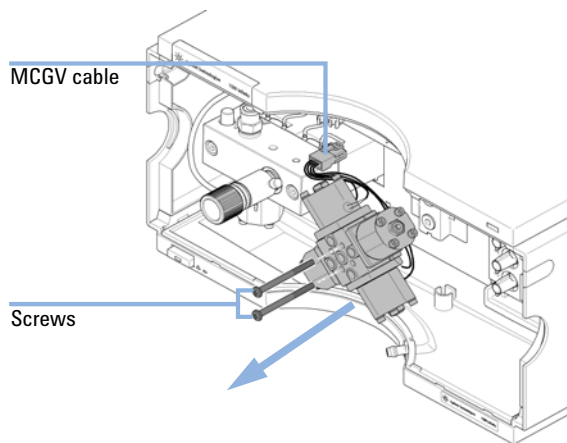


NOTE


The bio-inert version of the MCGV has no specific bio label and can be identified by its part number on actuator B.

Exchanging the Multi-Channel Gradient Valve (MCGV)


4 Place the new MCGV into position. Make sure that channel A of the MCGV is put at the bottom-right position. Tighten the two screws and connect the cable to its connector.

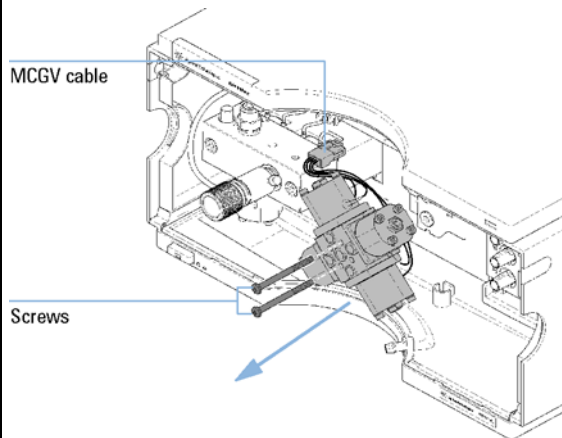


5 Install the MCGV cover.

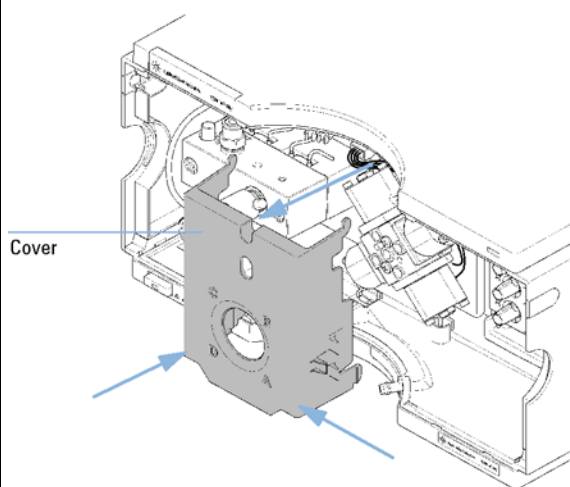

6 Reconnect the waste funnel with the waste tube holder in the top cover. Insert waste tube in the holder in the leak pan and clip tube to the MCGV cover.

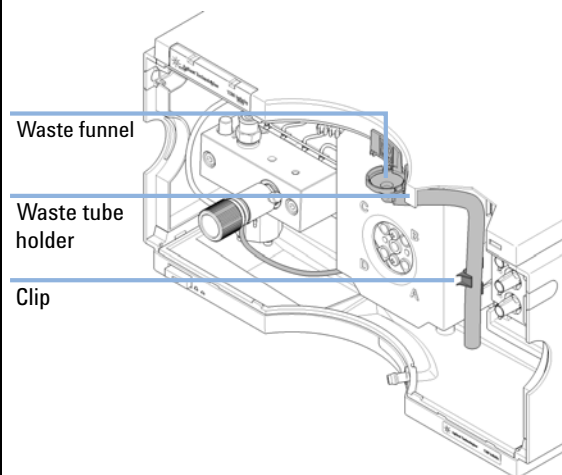

7 Reconnect the tube from the inlet valve to the middle position of the MCGV. Connect solvent tubes for channels A-D from the MCGV to the degasser outlets.

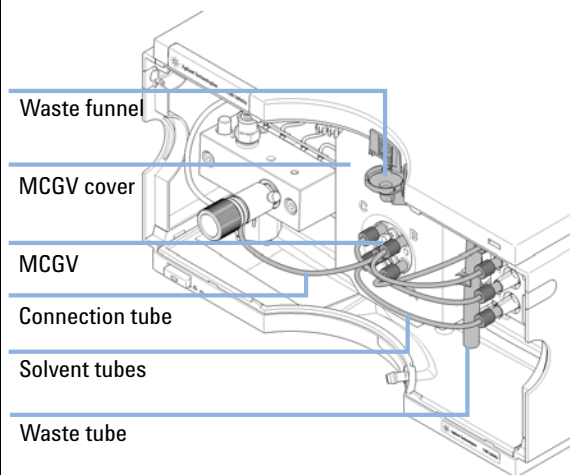

Exchanging the Optional Interface Board

When Board defective

Parts required	#	p/n	Description
	1	G1351-68701	Interface board (BCD) with external contacts and BCD outputs

CAUTION

Electronic boards are sensitive to electrostatic discharge (ESD) and should be handled with care so as not to damage them. Touching electronic boards and components can cause electrostatic discharge.

ESD can damage electronic boards and components.

- Be sure to hold the board by the edges and do not touch the electrical components. Always use an ESD protection (for example, an ESD wrist strap) when handling electronic boards and components.
-

Exchanging the Optional Interface Board

- 1 Switch off the pump at the main power switch, unplug the pump from line power.
- 2 Disconnect cables from the interface board connectors.
- 3 Loosen the screws. Slide out the interface board from the pump.

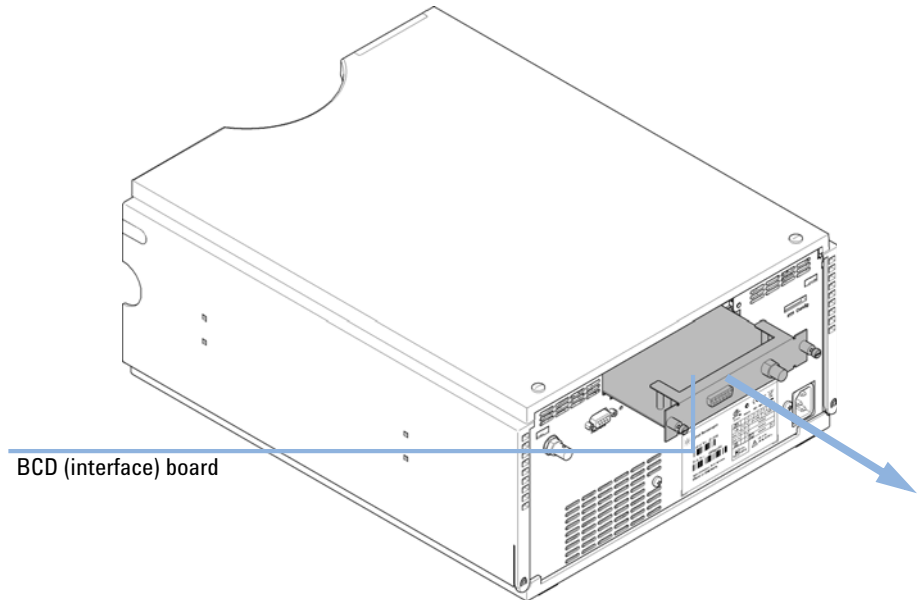


Figure 10 Exchanging the Interface Board

- 4 Install the new interface board. Secure screws.
- 5 Reconnect the cables to the board connector.
- 6 Reconnect the pump to line power.

Replacing the Module's Firmware

When	The installation of newer firmware might be necessary <ul style="list-style-type: none">• if a newer version solves problems of older versions or• to keep all systems on the same (validated) revision. The installation of older firmware might be necessary <ul style="list-style-type: none">• to keep all systems on the same (validated) revision or• if a new module with newer firmware is added to a system or• if third part control software requires a special version.
-------------	--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------

Tools required	Description
	LAN/RS-232 Firmware Update Tool
OR	Agilent Diagnostic Software
OR	Instant Pilot G4208A

Parts required	# Description
	1 Firmware, tools and documentation from Agilent web site

Preparations Read update documentation provided with the Firmware Update Tool.

To upgrade/downgrade the module's firmware carry out the following steps:

- 1 Download the required module firmware, the latest LAN/RS-232 FW Update Tool and the documentation from the Agilent web.
 - http://www.chem.agilent.com/scripts/cag_firmware.asp.
- 2 For loading the firmware into the module follow the instructions in the documentation.

Module Specific Information

There is no specific information for this module.



10 Parts for Maintenance

Pump Head Assembly with Seal Wash Option	138
Outlet Valve	140
Purge Valve Assembly	141
Active Inlet Valve	142
Accessory Kit, Bio-inert (G5611-68755)	144
Starter-Kit Bio-inert	145
Solvent Cabinet	146
Bottle Head Assembly	148
Hydraulic Path	149
System Tool Kit	151

This chapter provides information on parts for maintenance.



Pump Head Assembly with Seal Wash Option



For bio-inert modules use bio-inert parts only!

Item	p/n	Description
	G5611-60165	Pump Head with Seal Wash Option (Bio-inert)
1	5067-4695	Sapphire piston
2	G1312-60062	Piston housing (incl. spring)
3	G5611-63010	Support Ring with Seal Wash Option (Bio-inert)
4	0905-1731	Wash Seal PTFE for Bio-inert Pump 600 bar
	5065-9978	Silicone tubing, 1 mm i.d., 3 mm o.d., 5 m, re-order number
5	01018-07102	Gasket (Seal wash)
6	G5611-26210	Seal Holder (Bio-inert)
7	G4220-24013	Backup Seal for Support Ring
8	G5611-21503	Piston Seal PTFE (Bio-inert)
9	G5611-25260	Pump Chamber Housing (Bio-inert)
10	G5611-60025	Active Inlet Valve (AIV, Bio-inert), without cartridge
	G5611-60020	Cartridge for AIV (Bio-inert)
11	G5611-60067	Outlet Valve (Bio-inert)
12	5042-1303	Lock screw
13	G5611-60061	Purge Valve (Bio-inert)
14	0515-2118	Pump head screw (M5, 60 mm)
15	G1311-60161	1220/1260 Seal Wash Pump Assembly

Pump Head Assembly with Seal Wash Option

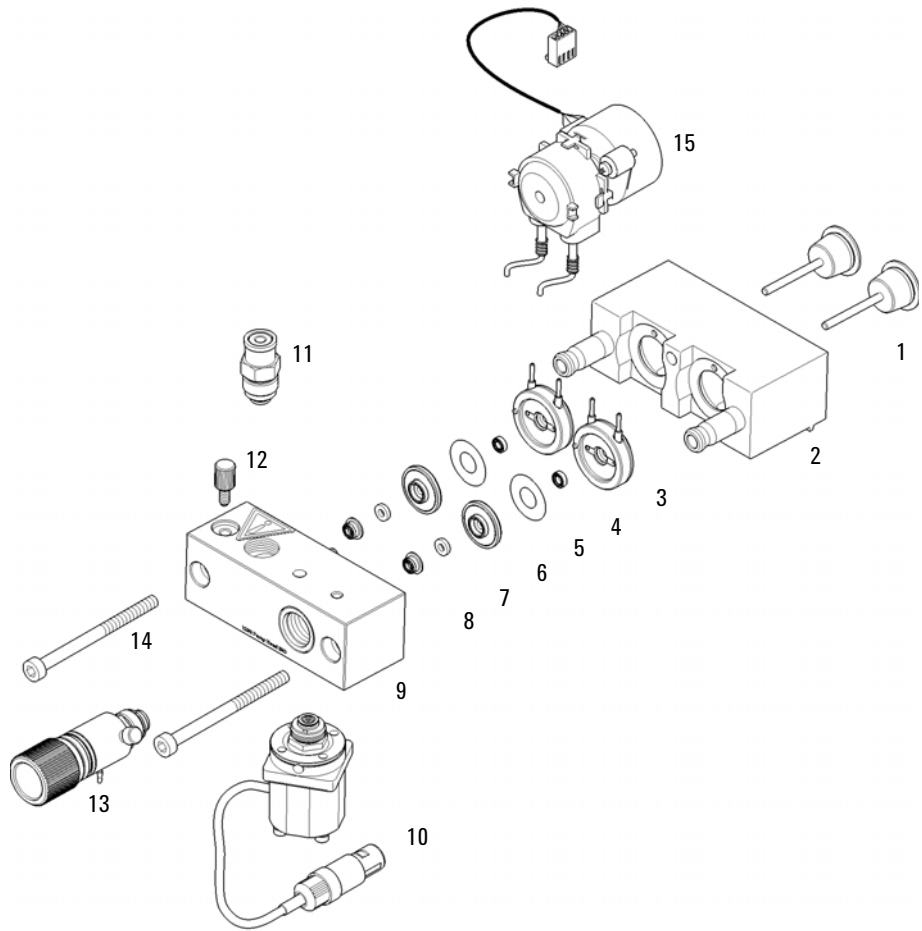


Figure 11 Pump Head with Seal Wash Option

Outlet Valve



For bio-inert modules use bio-inert parts only!

p/n	Description
G5611-60067	Outlet Valve (Bio-inert)

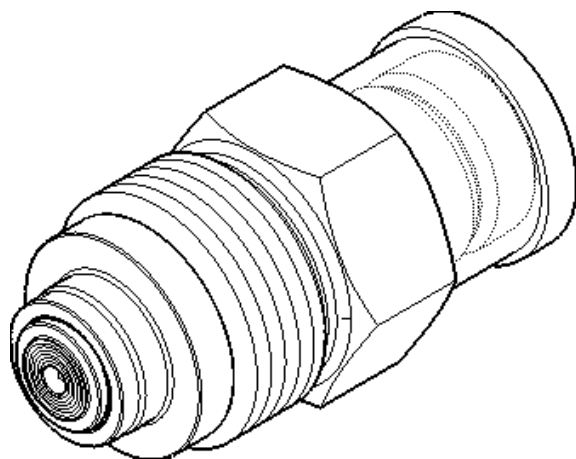


Figure 12 Outlet valve, bio-inert

Purge Valve Assembly



For bio-inert modules use bio-inert parts only!

Item	p/n	Description
1	G5611-60061	Purge Valve (Bio-inert)
2	01018-22707	PTFE frits (pack of 5)
3	5067-4728	Seal cap

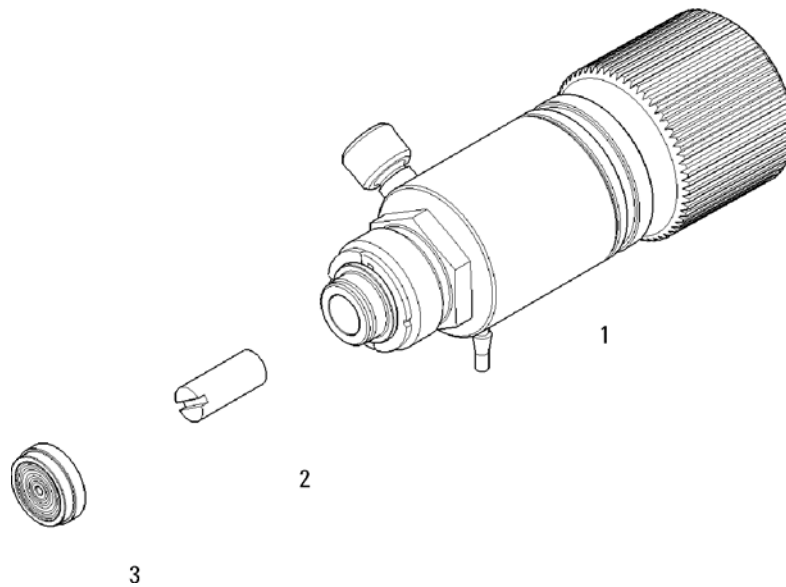


Figure 13 Purge valve assembly, bio-inert

Active Inlet Valve



For bio-inert modules use bio-inert parts only!

p/n	Description
G5611-60025	Active Inlet Valve (AIV, Bio-inert), without cartridge
G5611-60020	Cartridge for AIV (Bio-inert)
0100-1847	Adapter AIV to solvent inlet tubes

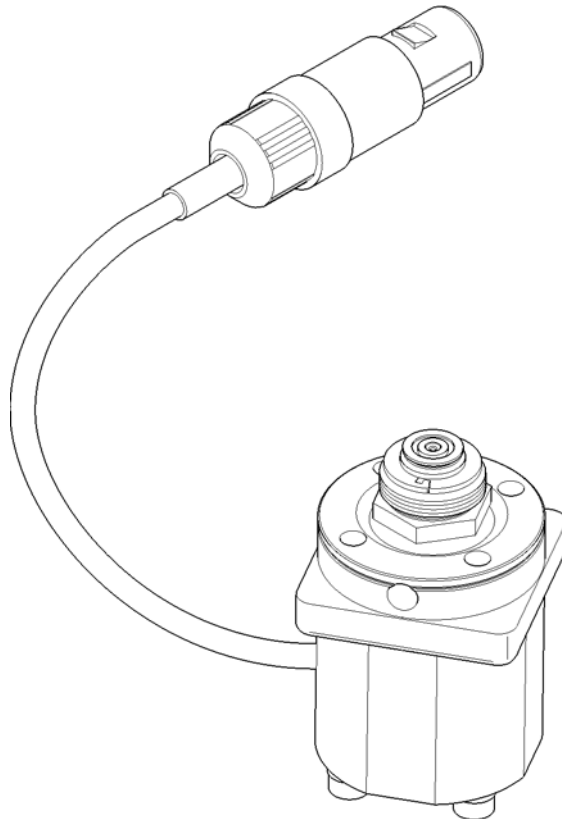


Figure 14 Active inlet valve

10 Parts for Maintenance

Accessory Kit, Bio-inert (G5611-68755)

Accessory Kit, Bio-inert (G5611-68755)

Accessory Kit (Bio-inert) (p/n G5611-68755)

p/n	Description
5062-2461	Waste tube, 5 m (reorder pack)
5063-6527	Tubing assembly, i.d. 6 mm, o.d. 9 mm, 1.2 m (to waste)
5181-1519	CAN cable, Agilent module to module, 1 m
G5611-60502	Capillary 900 x 0.17 mm, titanium (Bio-inert) pump to thermostatted autosampler
5042-9954	Tubing clip (2x), re-order 4/pk
G5611-60500	Capillary 400 x 0.17 mm, titanium (Bio-inert) pump to injector

Starter-Kit Bio-inert

Starter-Kit Bio-inert HPLC (p/n G5611-68707)

p/n	Description
9301-1420 (3x)	Solvent bottle, transparent
9301-1450	Solvent bottle, amber
01018-22707	PTFE frits (pack of 5)
5182-0716	Screw Cap Vial, 2 mL, amber glass, write-on spot, 100/pk
5182-0717	Blue screw caps 100/pk
5063-6507 (2x)	Chip, Column I.D. Assy
5041-2168 (2x)	Solvent inlet filter, 20 µm pore size
G5611-68710	Capillary/Fitting Starterkit ID. 17 mm (Bio-inert)

Solvent Cabinet

Item	p/n	Description
1	5065-9981	Solvent cabinet, including all plastic parts
2	5043-0207	Name plate 1260
3	5065-9954	Front panel, solvent cabinet
4	5042-8567	Leak pan
5	9301-1420	Solvent bottle, transparent
6	9301-1450	Solvent bottle, amber
7	G1311-60003	Bottle-head assembly

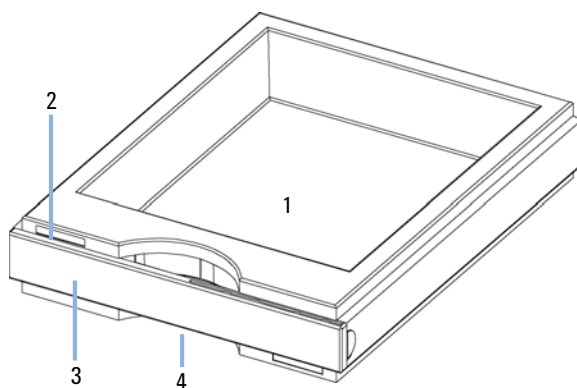




Figure 15 Solvent Cabinet Parts

Bottle Head Assembly

Item	p/n	Description
	G1311-60003	Bottle-head assembly
1	5063-6598	Ferrules with lock ring (10/Pk)
2	5063-6599	Tube screw (10/Pk)
3		Wire marker
4	5062-2483	Solvent tubing, 5 m
5	5062-8517	Inlet filter adapter (4/Pk)
6	5041-2168	Solvent inlet filter, 20 µm pore size

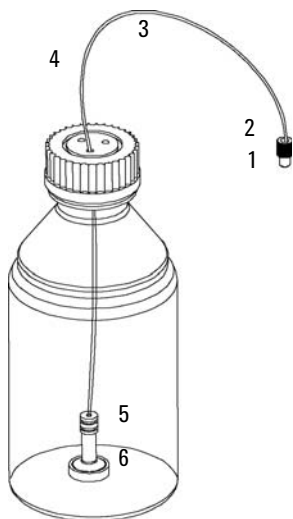


Figure 16 Bottle-Head Assembly Parts

Hydraulic Path



For bio-inert modules use bio-inert parts only!

Item	p/n	Description
1	G5611-60500	Capillary 400 x 0.17 mm, titanium (Bio-inert) pump to injector
1	G5611-60502	Capillary 900 x 0.17 mm, titanium (Bio-inert) pump to thermostatted autosampler
	G1311-60003	Bottle-head assembly
2	G1322-67300	Kit of 4 solvent tubes for connection degasser to MCGV including labels
3	G5611-67301	Capillary Chamber 1 to Damper (Bio-inert)
4	G5611-67300	Capillary Damper to Chamber 2 (Bio-inert)
5	G1311-67304	Connecting tube, MCGV to AIV
6	5062-2461	Waste tube, 5 m (reorder pack)
	0100-1847	Adapter AIV to solvent inlet tubes
	5067-4741	ZDV union (Bio-inert)

10 Parts for Maintenance

Hydraulic Path

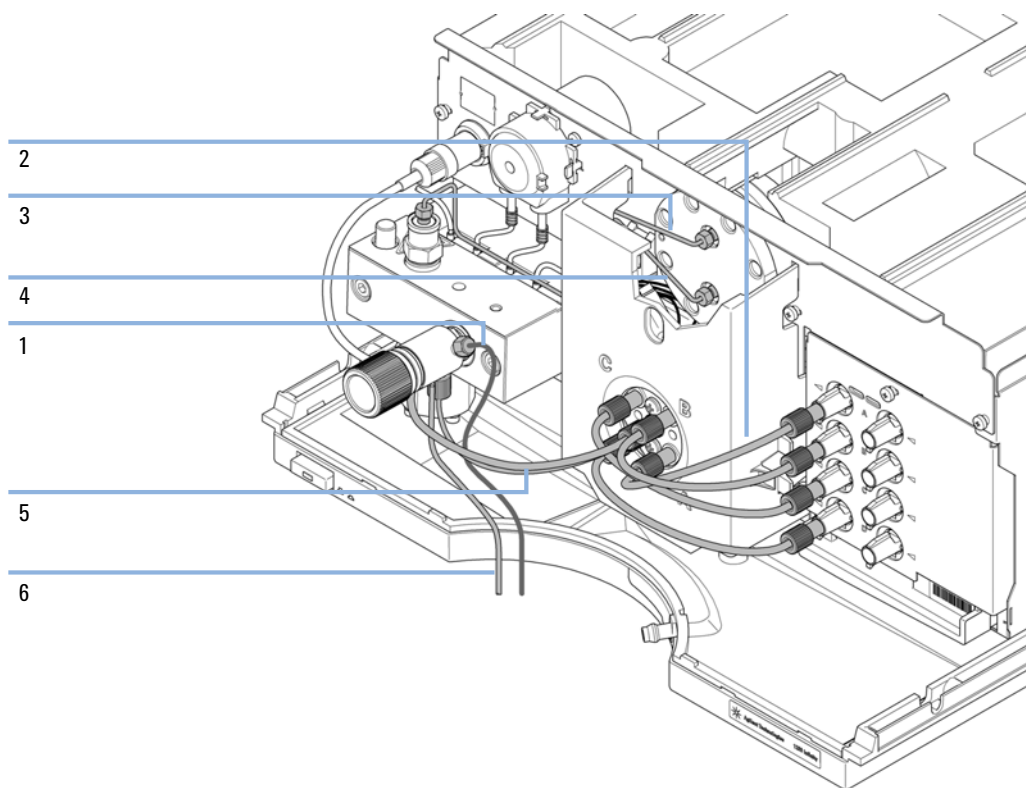


Figure 17 Hydraulic Flow Path

System Tool Kit

HPLC System Tool Kit (p/n G4203-68708)

p/n	Description
0100-1681	Adapter syringe/seal wash tube
0100-1710	Mounting Tool for Tubing Connections
01018-23702	Insert tool
5023-0240	Hex driver, ¼", slitted
8710-0060	Hex-key wrench, 9/64 inch
8710-0510 (2x)	Wrench open 1/4 — 5/16 inch
8710-0641	Hex key set 1 – 5 mm
8710-0899	Pozidriv screwdriver
8710-1534	Wrench, 4 mm both ends, open end
8710-1924	Wrench open 14 mm
8710-2392	Hex key 4 mm 15 cm long T-handle
8710-2393	Hex key 1.5 mm, straight handle 10 cm
8710-2394	Hex key 9/64 inch 15 cm long T-handle
8710-2409	Wrench open end, 5/16 – 3/8 inch
8710-2411	Hex key 3 mm 12 cm long
8710-2412	Hex key 2.5 mm, 15 cm long, straight handle
8710-2438	Hex key 2.0 mm
8710-2509	Screwdriver Torx TX8
8710-2594	Double open end wrench 4x5
9301-0411	Syringe, Plastic
9301-1337	Adapter syringe/solvent tube with fitting

Preventive Maintenance Kit



For bio-inert modules use bio-inert parts only!

p/n	Description
G5611-21503 (2x)	Piston Seal PTFE (Bio-inert)
01018-22707	PTFE frits (pack of 5)
5067-4728	Seal cap
0905-1731 (2x)	Wash Seal PTFE for Bio-inert Pump 600 bar
01018-07102 (2x)	Gasket (Seal wash)
5042-8507	Seal wash pump cartridge (silicone tubing)



11 Identifying Cables

Cable Overview	154
Analog Cables	156
Remote Cables	158
BCD Cables	161
CAN/LAN Cables	163
External Contact Cable	164
Agilent Module to PC	165
Agilent 1200 Module to Printer	166

This chapter provides information on cables used with the Agilent 1200 Infinity Series modules.



Cable Overview

NOTE

Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.

Analog cables

p/n	Description
35900-60750	Agilent module to 3394/6 integrators
35900-60750	Agilent 35900A A/D converter
01046-60105	Analog cable (BNC to general purpose, spade lugs)

Remote cables

p/n	Description
03394-60600	Agilent module to 3396A Series I integrators 3396 Series II / 3395A integrator, see details in section “Remote Cables” on page 158
03396-61010	Agilent module to 3396 Series III / 3395B integrators
5061-3378	Agilent module to Agilent 35900 A/D converters (or HP 1050/1046A/1049A)
01046-60201	Agilent module to general purpose

BCD cables

p/n	Description
03396-60560	Agilent module to 3396 integrators
G1351-81600	Agilent module to general purpose

CAN cables

p/n	Description
5181-1516	CAN cable, Agilent module to module, 0.5 m
5181-1519	CAN cable, Agilent module to module, 1 m

LAN cables

p/n	Description
5023-0203	Cross-over network cable, shielded, 3 m (for point to point connection)
5023-0202	Twisted pair network cable, shielded, 7 m (for point to point connection)


External Contact Cable

p/n	Description
G1103-61611	External contact cable - Agilent module interface board to general purposes

RS-232 cables

p/n	Description
G1530-60600	RS-232 cable, 2 m
RS232-61600	RS-232 cable, 2.5 m Instrument to PC, 9-to-9 pin (female). This cable has special pin-out, and is not compatible with connecting printers and plotters. It's also called "Null Modem Cable" with full handshaking where the wiring is made between pins 1-1, 2-3, 3-2, 4-6, 5-5, 6-4, 7-8, 8-7, 9-9.
5181-1561	RS-232 cable, 8 m

Analog Cables

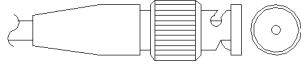

One end of these cables provides a BNC connector to be connected to Agilent modules. The other end depends on the instrument to which connection is being made.


Agilent Module to 3394/6 Integrators

p/n 35900-60750	Pin 3394/6	Pin Agilent module	Signal Name
	1		Not connected
	2	Shield	Analog -
	3	Center	Analog +


Agilent Module to BNC Connector

p/n 8120-1840	Pin BNC	Pin Agilent module	Signal Name
	Shield	Shield	Analog -
	Center	Center	Analog +

Agilent Module to General Purpose

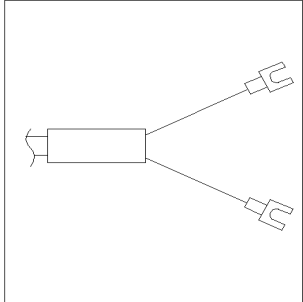p/n 01046-60105	Pin	Pin Agilent module	Signal Name
	1		Not connected
	2	Black	Analog -
	3	Red	Analog +

Remote Cables

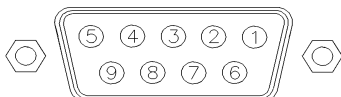

One end of these cables provides a Agilent Technologies APG (Analytical Products Group) remote connector to be connected to Agilent modules. The other end depends on the instrument to be connected to.

Agilent Module to 3396A Integrators

p/n 03394-60600	Pin 3396A	Pin Agilent module	Signal Name	Active (TTL)
	9	1 - White	Digital ground	
	NC	2 - Brown	Prepare run	Low
	3	3 - Gray	Start	Low
	NC	4 - Blue	Shut down	Low
	NC	5 - Pink	Not connected	
	NC	6 - Yellow	Power on	High
	5,14	7 - Red	Ready	High
	1	8 - Green	Stop	Low
	NC	9 - Black	Start request	Low
	13, 15		Not connected	

Agilent Module to 3396 Series II / 3395A Integrators

Use the cable Agilent module to 3396A Series I integrators (p/n 03394-60600) and cut pin #5 on the integrator side. Otherwise the integrator prints START; not ready.

Agilent Module to 3396 Series III / 3395B Integrators

p/n 03396-61010	Pin 33XX	Pin Agilent module	Signal Name	Active (TTL)
	9	1 - White	Digital ground	
	NC	2 - Brown	Prepare run	Low
	3	3 - Gray	Start	Low
	NC	4 - Blue	Shut down	Low
	NC	5 - Pink	Not connected	
	NC	6 - Yellow	Power on	High
	14	7 - Red	Ready	High
	4	8 - Green	Stop	Low
	NC	9 - Black	Start request	Low
	13, 15		Not connected	


Agilent Module to Agilent 35900 A/D Converters

p/n 5061-3378	Pin 35900 A/D	Pin Agilent module	Signal Name	Active (TTL)
	1 - White	1 - White	Digital ground	
	2 - Brown	2 - Brown	Prepare run	Low
	3 - Gray	3 - Gray	Start	Low
	4 - Blue	4 - Blue	Shut down	Low
	5 - Pink	5 - Pink	Not connected	
	6 - Yellow	6 - Yellow	Power on	High
	7 - Red	7 - Red	Ready	High
	8 - Green	8 - Green	Stop	Low
	9 - Black	9 - Black	Start request	Low

11 Identifying Cables

Remote Cables

Agilent Module to General Purpose

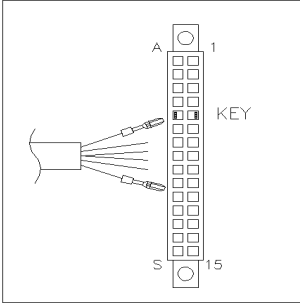p/n 01046-60201	Pin Universal	Pin Agilent module	Signal Name	Active (TTL)
		1 - White	Digital ground	
		2 - Brown	Prepare run	Low
		3 - Gray	Start	Low
		4 - Blue	Shut down	Low
		5 - Pink	Not connected	
		6 - Yellow	Power on	High
		7 - Red	Ready	High
		8 - Green	Stop	Low
		9 - Black	Start request	Low

BCD Cables



One end of these cables provides a 15-pin BCD connector to be connected to the Agilent modules. The other end depends on the instrument to be connected to


Agilent Module to General Purpose

p/n G1351-81600	Wire Color	Pin Agilent module	Signal Name	BCD Digit
	Green	1	BCD 5	20
	Violet	2	BCD 7	80
	Blue	3	BCD 6	40
	Yellow	4	BCD 4	10
	Black	5	BCD 0	1
	Orange	6	BCD 3	8
	Red	7	BCD 2	4
	Brown	8	BCD 1	2
	Gray	9	Digital ground	Gray
	Gray/pink	10	BCD 11	800
	Red/blue	11	BCD 10	400
	White/green	12	BCD 9	200
	Brown/green	13	BCD 8	100
	not connected	14		
	not connected	15	+ 5 V	Low


11 Identifying Cables

BCD Cables

Agilent Module to 3396 Integrators

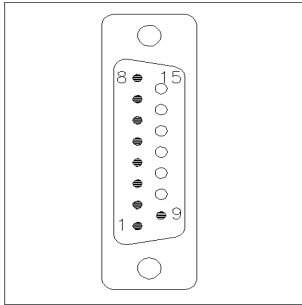p/n 03396-60560	Pin 3396	Pin Agilent module	Signal Name	BCD Digit
	1	1	BCD 5	20
	2	2	BCD 7	80
	3	3	BCD 6	40
	4	4	BCD 4	10
	5	5	BCD0	1
	6	6	BCD 3	8
	7	7	BCD 2	4
	8	8	BCD 1	2
	9	9	Digital ground	
	NC	15	+ 5 V	Low

CAN/LAN Cables

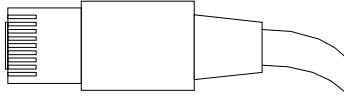

Both ends of this cable provide a modular plug to be connected to Agilent modules CAN or LAN connectors.


CAN Cables

p/n	Description
5181-1516	CAN cable, Agilent module to module, 0.5 m
5181-1519	CAN cable, Agilent module to module, 1 m

LAN Cables

p/n	Description
5023-0203	Cross-over network cable, shielded, 3 m (for point to point connection)
5023-0202	Twisted pair network cable, shielded, 7 m (for point to point connection)

External Contact Cable

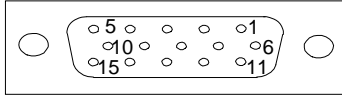

One end of this cable provides a 15-pin plug to be connected to Agilent modules interface board. The other end is for general purpose.

Agilent Module Interface Board to general purposes

p/n G1103-61611	Color	Pin Agilent module	Signal Name
	White	1	EXT 1
	Brown	2	EXT 1
	Green	3	EXT 2
	Yellow	4	EXT 2
	Grey	5	EXT 3
	Pink	6	EXT 3
	Blue	7	EXT 4
	Red	8	EXT 4
	Black	9	Not connected
	Violet	10	Not connected
	Grey/pink	11	Not connected
	Red/blue	12	Not connected
	White/green	13	Not connected
	Brown/green	14	Not connected
	White/yellow	15	Not connected

Agilent Module to PC

p/n	Description
G1530-60600	RS-232 cable, 2 m
RS232-61600	RS-232 cable, 2.5 m Instrument to PC, 9-to-9 pin (female). This cable has special pin-out, and is not compatible with connecting printers and plotters. It's also called "Null Modem Cable" with full handshaking where the wiring is made between pins 1-1, 2-3, 3-2, 4-6, 5-5, 6-4, 7-8, 8-7, 9-9.
5181-1561	RS-232 cable, 8 m

11 Identifying Cables

Agilent 1200 Module to Printer

Agilent 1200 Module to Printer

p/n	Description
5181-1529	Cable Printer Serial & Parallel, is a SUB-D 9 pin female vs. Centronics connector on the other end (NOT FOR FW UPDATE). For use with G1323 Control Module.



12 Hardware Information

Electrical Connections	168
Rear View of the Module	169
Interfaces	170
Overview Interfaces	172
Setting the 8-bit Configuration Switch (without On-board) LAN	176
Communication Settings for RS-232C	177
Special Settings	179

This chapter describes the pump in more detail on hardware and electronics.



Electrical Connections

- The CAN bus is a serial bus with high speed data transfer. The two connectors for the CAN bus are used for internal module data transfer and synchronization.
- One analog output provides signals for integrators or data handling systems.
- The REMOTE connector may be used in combination with other analytical instruments from Agilent Technologies if you want to use features such as start, stop, common shut down, prepare, and so on.
- With the appropriate software, the RS-232C connector may be used to control the module from a computer through a RS-232C connection. This connector is activated and can be configured with the configuration switch.
- The power input socket accepts a line voltage of 100 – 240 VAC \pm 10 % with a line frequency of 50 or 60 Hz. Maximum power consumption varies by module. There is no voltage selector on your module because the power supply has wide-ranging capability. There are no externally accessible fuses, because automatic electronic fuses are implemented in the power supply.

NOTE

Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations.

Rear View of the Module

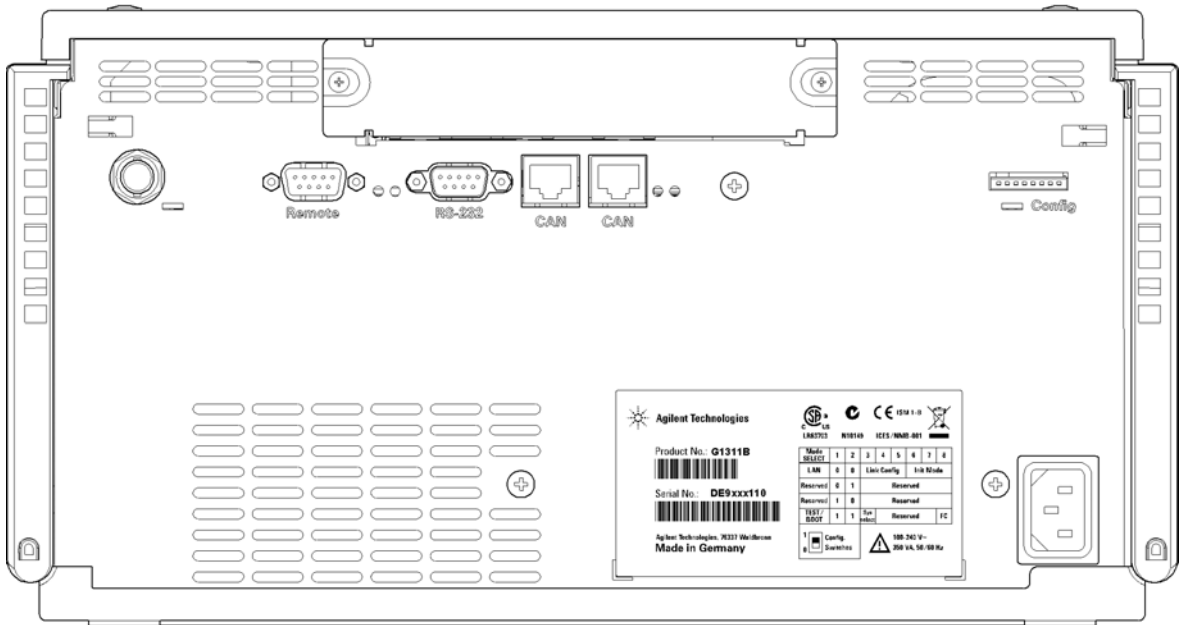


Figure 18 Rear view of the pump

Interfaces

The Agilent 1200 Infinity Series modules provide the following interfaces:

Table 9 Agilent 1200 Infinity Series Interfaces

Module	CAN	LAN/BCD (optional)	LAN (on-board)	RS-232	Analog	APG Remote	Special
Pumps							
G1310B Iso Pump G1311B Quat Pump G1311C Quat Pump VL G1312B Bin Pump G1312C Bin Pump VL 1376A Cap Pump G2226A Nano Pump G5611A Bio-inert Quat Pump	2	Yes	No	Yes	1	Yes	
G4220A/B Bin Pump	2	No	Yes	Yes	No	Yes	
G1361A Prep Pump	2	Yes	No	Yes	No	Yes	CAN-DC- OUT for CAN slaves
Samplers							
G1329B ALS G2260A Prep ALS	2	Yes	No	Yes	No	Yes	THERMOSTAT for G1330B
G1364B FC-PS G1364C FC-AS G1364D FC- μ S G1367E HiP ALS G1377A HiP micro ALS G2258A DL ALS G5664A Bio-inert FC-AS G5667A Bio-inert Autosampler	2	Yes	No	Yes	No	Yes	THERMOSTAT for G1330B CAN-DC- OUT for CAN slaves
G4226A ALS	2	Yes	No	Yes	No	Yes	

Table 9 Agilent 1200 Infinity Series Interfaces

Module	CAN	LAN/BCD (optional)	LAN (on-board)	RS-232	Analog	APG Remote	Special
Detectors							
G1314B VWD VL G1314C VWD VL+	2	Yes	No	Yes	1	Yes	
G1314E/F VWD	2	No	Yes	Yes	1	Yes	
G4212A/B DAD	2	No	Yes	Yes	1	Yes	
G1315C DAD VL+ G1365C MWD G1315D DAD VL G1365D MWD VL	2	No	Yes	Yes	2	Yes	
G1321B FLD G1362A RID	2	Yes	No	Yes	1	Yes	
G4280A ELSD	No	No	No	Yes	Yes	Yes	EXT Contact AUTOZERO
Others							
G1170A Valve Drive	2	No	No	No	No	No	Requires a HOST module with on-board LAN (e.g. G4212A or G4220A with minimum firmware B.06.40 or C.06.40) or with additional G1369C LAN Card
G1316A/C TCC	2	No	No	Yes	No	Yes	
G1322A DEG	No	No	No	No	No	Yes	AUX
G1379B DEG	No	No	No	Yes	No	No	AUX
G4227A Flex Cube	2	No	No	No	No	No	
G4240A CHIP CUBE	2	Yes	No	Yes	No	Yes	CAN-DC- OUT for CAN slaves THERMOSTAT for G1330A/B (NOT USED)

NOTE

The detector (DAD/MWD/FLD/VWD/RID) is the preferred access point for control via LAN. The inter-module communication is done via CAN.

- CAN connectors as interface to other modules
- LAN connector as interface to the control software
- RS-232C as interface to a computer
- REMOTE connector as interface to other Agilent products
- Analog output connector(s) for signal output

Overview Interfaces

CAN

The CAN is inter-module communication interface. It is a 2-wire serial bus system supporting high speed data communication and real-time requirement.

LAN

The modules have either an interface slot for an LAN card (e.g. Agilent G1369A/B LAN Interface) or they have an on-board LAN interface (e.g. detectors G1315C/D DAD and G1365C/D MWD). This interface allows the control of the module/system via a connected PC with the appropriate control software.

NOTE

If an Agilent detector (DAD/MWD/FLD/VWD/RID) is in the system, the LAN should be connected to the DAD/MWD/FLD/VWD/RID (due to higher data load). If no Agilent detector is part of the system, the LAN interface should be installed in the pump or autosampler.

RS-232C (Serial)

The RS-232C connector is used to control the module from a computer through RS-232C connection, using the appropriate software. This connector can be configured with the configuration switch module at the rear of the module. Refer to *Communication Settings for RS-232C*.

NOTE

There is no configuration possible on main boards with on-board LAN. These are pre-configured for

- 19200 baud,
- 8 data bit with no parity and
- one start bit and one stop bit are always used (not selectable).

The RS-232C is designed as DCE (data communication equipment) with a 9-pin male SUB-D type connector. The pins are defined as:

Table 10 RS-232C Connection Table

Pin	Direction	Function
1	In	DCD
2	In	RxD
3	Out	TxD
4	Out	DTR
5		Ground
6	In	DSR
7	Out	RTS
8	In	CTS
9	In	RI

12 Hardware Information

Interfaces

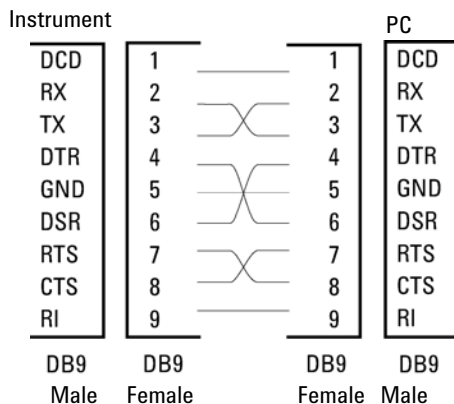


Figure 19 RS-232 Cable

Analog Signal Output

The analog signal output can be distributed to a recording device. For details refer to the description of the module's main board.

APG Remote

The APG Remote connector may be used in combination with other analytical instruments from Agilent Technologies if you want to use features as common shut down, prepare, and so on.

Remote control allows easy connection between single instruments or systems to ensure coordinated analysis with simple coupling requirements.

The subminiature D connector is used. The module provides one remote connector which is inputs/outputs (wired- or technique).

To provide maximum safety within a distributed analysis system, one line is dedicated to **SHUT DOWN** the system's critical parts in case any module detects a serious problem. To detect whether all participating modules are switched on or properly powered, one line is defined to summarize the **POWER ON** state of all connected modules. Control of analysis is maintained by signal readiness **READY** for next analysis, followed by **START** of run and optional **STOP** of run triggered on the respective lines. In addition **PREPARE** and **START REQUEST** may be issued. The signal levels are defined as:

- standard TTL levels (0 V is logic true, + 5.0 V is false),
- fan-out is 10,
- input load is 2.2 kOhm against + 5.0 V, and
- output are open collector type, inputs/outputs (wired- or technique).

NOTE

All common TTL circuits operate with a 5 V power supply. A TTL signal is defined as "low" or L when between 0 V and 0.8 V and "high" or H when between 2.0 V and 5.0 V (with respect to the ground terminal).

Table 11 Remote Signal Distribution

Pin	Signal	Description
1	DGND	Digital ground
2	PREPARE	(L) Request to prepare for analysis (for example, calibration, detector lamp on). Receiver is any module performing pre-analysis activities.
3	START	(L) Request to start run / timetable. Receiver is any module performing run-time controlled activities.
4	SHUT DOWN	(L) System has serious problem (for example, leak: stops pump). Receiver is any module capable to reduce safety risk.
5		Not used
6	POWER ON	(H) All modules connected to system are switched on. Receiver is any module relying on operation of others.
7	READY	(H) System is ready for next analysis. Receiver is any sequence controller.
8	STOP	(L) Request to reach system ready state as soon as possible (for example, stop run, abort or finish and stop injection). Receiver is any module performing run-time controlled activities.
9	START REQUEST	(L) Request to start injection cycle (for example, by start key on any module). Receiver is the autosampler.

Special Interfaces

Some modules have module specific interfaces/connectors. They are described in the module documentation.

Setting the 8-bit Configuration Switch (without On-board) LAN

The 8-bit configuration switch is located at the rear of the module.

Modules that do not have their own LAN interface (e.g. the TCC) can be controlled through the LAN interface of another module and a CAN connection to that module.

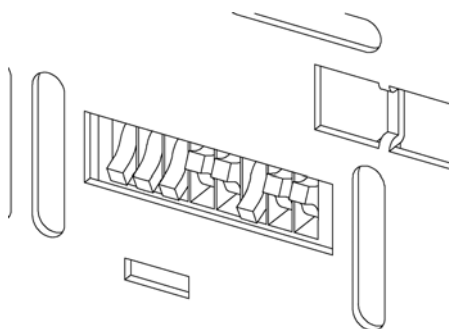


Figure 20 Configuration switch (settings depend on configured mode)

All modules without on-board LAN:

- default should be ALL DIPS DOWN (= best settings)
 - Bootp mode for LAN and
 - 19200 baud, 8 data bit / 1 stop bit with no parity for RS-232
- DIP 1 DOWN and DIP 2 UP allows special RS-232 settings
- for boot/test modes DIPS 1+2 must be UP plus required mode

NOTE

For normal operation use the default (best) settings.

Switch settings provide configuration parameters for serial communication protocol and instrument specific initialization procedures.

NOTE

With the introduction of the Agilent 1260 Infinity, all GPIB interfaces have been removed. The preferred communication is LAN.

NOTE

The following tables represent the configuration switch settings for the modules without on-board LAN only.

Table 12 8-bit Configuration Switch (without on-board LAN)

Mode Select	1	2	3	4	5	6	7	8
RS-232C	0	1	Baudrate			Data Bits	Parity	
Reserved	1	0	Reserved					
TEST/BOOT	1	1	RSVD	SYS		RSVD	RSVD	FC

NOTE

The LAN settings are done on the LAN Interface Card G1369A/B. Refer to the documentation provided with the card.

Communication Settings for RS-232C

The communication protocol used in the column compartment supports only hardware handshake (CTS/RTR).

Switches 1 in down and 2 in up position define that the RS-232C parameters will be changed. Once the change has been completed, the column instrument must be powered up again in order to store the values in the non-volatile memory.

Table 13 Communication Settings for RS-232C Communication (without on-board LAN)

Mode Select	1	2	3	4	5	6	7	8
RS-232C	0	1	Baudrate			Data Bits	Parity	

Use the following tables for selecting the setting which you want to use for RS-232C communication. The number 0 means that the switch is down and 1 means that the switch is up.

12 Hardware Information

Setting the 8-bit Configuration Switch (without On-board) LAN

Table 14 Baudrate Settings (without on-board LAN)

Switches			Baud Rate	Switches			Baud Rate
3	4	5		3	4	5	
0	0	0	9600	1	0	0	9600
0	0	1	1200	1	0	1	14400
0	1	0	2400	1	1	0	19200
0	1	1	4800	1	1	1	38400

Table 15 Data Bit Settings (without on-board LAN)

Switch 6	Data Word Size
0	7 Bit Communication
1	8 Bit Communication

Table 16 Parity Settings (without on-board LAN)

Switches		Parity
7	8	
0	0	No Parity
0	1	Odd Parity
1	1	Even Parity

One start bit and one stop bit are always used (not selectable).

Per default, the module will turn into 19200 baud, 8 data bit with no parity.

Special Settings

The special settings are required for specific actions (normally in a service case).

Boot-Resident

Firmware update procedures may require this mode in case of firmware loading errors (main firmware part).

If you use the following switch settings and power the instrument up again, the instrument firmware stays in the resident mode. It is not operable as a module. It only uses basic functions of the operating system for example, for communication. In this mode the main firmware can be loaded (using update utilities).

Table 17 Boot Resident Settings (without on-board LAN)

	Mode Select	SW1	SW2	SW3	SW4	SW5	SW6	SW7	SW8
No LAN	TEST/BOOT	1	1	0	0	1	0	0	0

Forced Cold Start

A forced cold start can be used to bring the module into a defined mode with default parameter settings.

CAUTION

Loss of data

Forced cold start erases all methods and data stored in the non-volatile memory. Exceptions are diagnosis and repair log books which will not be erased.

→ Save your methods and data before executing a forced cold start.

If you use the following switch settings and power the instrument up again, a forced cold start has been completed.

Table 18 Forced Cold Start Settings (without on-board LAN)

	Mode Select	SW1	SW2	SW3	SW4	SW5	SW6	SW7	SW8
No LAN	TEST/BOOT	1	1	0	0	1	0	0	1

12 Hardware Information

Setting the 8-bit Configuration Switch (without On-board) LAN



13 Appendix

General Safety Information	182
The Waste Electrical and Electronic Equipment Directive	185
Batteries Information	186
Radio Interference	187
Sound Emission	188
Agilent Technologies on Internet	189

This chapter provides addition information on safety, legal and web.



General Safety Information

General Safety Information

The following general safety precautions must be observed during all phases of operation, service, and repair of this instrument. Failure to comply with these precautions or with specific warnings elsewhere in this manual violates safety standards of design, manufacture, and intended use of the instrument. Agilent Technologies assumes no liability for the customer's failure to comply with these requirements.

WARNING

Ensure the proper usage of the equipment.

The protection provided by the equipment may be impaired.

→ The operator of this instrument is advised to use the equipment in a manner as specified in this manual.

Safety Standards

This is a Safety Class I instrument (provided with terminal for protective earthing) and has been manufactured and tested according to international safety standards.

Operation

Before applying power, comply with the installation section. Additionally the following must be observed.

Do not remove instrument covers when operating. Before the instrument is switched on, all protective earth terminals, extension cords, auto-transformers, and devices connected to it must be connected to a protective earth via a ground socket. Any interruption of the protective earth grounding will cause a potential shock hazard that could result in serious personal injury. Whenever it is likely that the protection has been impaired,

the instrument must be made inoperative and be secured against any intended operation.

Make sure that only fuses with the required rated current and of the specified type (normal blow, time delay, and so on) are used for replacement. The use of repaired fuses and the short-circuiting of fuse holders must be avoided.

Some adjustments described in the manual, are made with power supplied to the instrument, and protective covers removed. Energy available at many points may, if contacted, result in personal injury.

Any adjustment, maintenance, and repair of the opened instrument under voltage should be avoided whenever possible. When inevitable, this has to be carried out by a skilled person who is aware of the hazard involved. Do not attempt internal service or adjustment unless another person, capable of rendering first aid and resuscitation, is present. Do not replace components with power cable connected.

Do not operate the instrument in the presence of flammable gases or fumes. Operation of any electrical instrument in such an environment constitutes a definite safety hazard.






Do not install substitute parts or make any unauthorized modification to the instrument.

Capacitors inside the instrument may still be charged, even though the instrument has been disconnected from its source of supply. Dangerous voltages, capable of causing serious personal injury, are present in this instrument. Use extreme caution when handling, testing and adjusting.

When working with solvents please observe appropriate safety procedures (e.g. goggles, safety gloves and protective clothing) as described in the material handling and safety data sheet by the solvent vendor, especially when toxic or hazardous solvents are used.

Safety Symbols

Table 19 Safety Symbols

Symbol	Description
	The apparatus is marked with this symbol when the user should refer to the instruction manual in order to protect risk of harm to the operator and to protect the apparatus against damage.
	Indicates dangerous voltages.
	Indicates a protected ground terminal.
	Indicates eye damage may result from directly viewing the light produced by the deuterium lamp used in this product.
	The apparatus is marked with this symbol when hot surfaces are available and the user should not touch it when heated up.

WARNING

A WARNING

alerts you to situations that could cause physical injury or death.

- Do not proceed beyond a warning until you have fully understood and met the indicated conditions.

CAUTION

A CAUTION

alerts you to situations that could cause loss of data, or damage of equipment.

- Do not proceed beyond a caution until you have fully understood and met the indicated conditions.

The Waste Electrical and Electronic Equipment Directive

Abstract

The Waste Electrical and Electronic Equipment (WEEE) Directive (2002/96/EC), adopted by EU Commission on 13 February 2003, is introducing producer responsibility on all electric and electronic appliances starting with 13 August 2005.

NOTE

This product complies with the WEEE Directive (2002/96/EC) marking requirements. The affixed label indicates that you must not discard this electrical/electronic product in domestic household waste.

Product Category:

With reference to the equipment types in the WEEE Directive Annex I, this product is classed as a Monitoring and Control Instrumentation product.



NOTE

Do not dispose off in domestic household waste

To return unwanted products, contact your local Agilent office, or see www.agilent.com for more information.

Batteries Information

WARNING

Lithium batteries may not be disposed-off into the domestic waste. Transportation of discharged Lithium batteries through carriers regulated by IATA/ICAO, ADR, RID, IMDG is not allowed.

Danger of explosion if battery is incorrectly replaced.

- Discharged Lithium batteries shall be disposed off locally according to national waste disposal regulations for batteries.
 - Replace only with the same or equivalent type recommended by the equipment manufacturer.
-



WARNING

Lithiumbatteri - Eksplosionsfare ved fejlagtig håndtering.

Udskiftning må kun ske med batteri af samme fabrikat og type.

- Lever det brugte batteri tilbage til leverandøren.
-

WARNING

Lithiumbatteri - Eksplosionsfare.

Ved udskiftning benyttes kun batteri som anbefalt av apparatfabrikanten.

- Brukt batteri returneres apparatleverandøren.
-

NOTE

Bij dit apparaat zijn batterijen geleverd. Wanneer deze leeg zijn, moet u ze niet weggooien maar inleveren als KCA.

Radio Interference

Cables supplied by Agilent Technologies are screened to provide optimized protection against radio interference. All cables are in compliance with safety or EMC regulations.

Test and Measurement

If test and measurement equipment is operated with unscreened cables, or used for measurements on open set-ups, the user has to assure that under operating conditions the radio interference limits are still met within the premises.

Sound Emission

Manufacturer's Declaration

This statement is provided to comply with the requirements of the German Sound Emission Directive of 18 January 1991.

This product has a sound pressure emission (at the operator position) < 70 dB.

- Sound Pressure $L_p < 70$ dB (A)
- At Operator Position
- Normal Operation
- According to ISO 7779:1988/EN 27779/1991 (Type Test)

Agilent Technologies on Internet

For the latest information on products and services visit our worldwide web site on the Internet at:

<http://www.agilent.com>

Select Products/Chemical Analysis

It will provide also the latest firmware of the modules for download.

Index

8

8-bit configuration switch
without On-Board LAN 176

A

accessory kit 29, 144
active inlet valve cartridge 114
active inlet valve 114, 142
active seal wash 10, 58
adapter 114
Agilent Diagnostic software 69
Agilent Lab Advisor software 69
Agilent Lab Advisor 69
Agilent
on internet 189
algae growth 49
alternative seal material 59
ambient non-operating temperature 25
ambient operating temperature 25
analog signal 174
analog
cable 156
apg remote 174
AUTO mode 16

B

ball-screw drive 13
battery
safety information 186
BCD
cable 161
bench space 24
bio-inert

active inlet valve 114
materials 17
multi channel gradient valve
(MCGV) 131
outlet valve 117
pump head 126
purge valve 119

blockage 86
bottle head assembly 148
buffer application 48
buffer solution 10

C

cable
analog 156
BCD 161
CAN 163
external contact 164
interface 35
LAN 163
remote 158
RS-232 165

cables
analog 154
BCD 154
CAN 155
external contact 155
LAN 155
overview 154
remote 154
RS-232 155

CAN
cable 163
changing solvents 44
cleaning 112

Communication settings
RS-232C 177
compensation sensor open 78
compensation sensor short 78
compressibility compensation 15, 60
condensation 24
configuration
one stack 30
connections, flow 38
control software 37

D

damaged parts 28
defect on arrival 28
degasser cannot read signal 94
degasser limit not reached 94
degasser, operational hints 56
degasser, when to use 56
Diagnostic software 69
dimensions 25
dual-piston in-series design 11

E

electrical connections
descriptions of 168
electronic waste 185
electrostatic discharge (ESD) 134
EMF
early maintenance feedback 19
encoder missing 87
error messages, degasser cannot read
signal 94
error messages, degasser limit not
reached 94

Index

- error messages
 - pump error 80
- error messages
 - compensation sensor open 78
 - compensation sensor short 78
 - encoder missing 87
 - fan failed 79
 - ignition without cover 79, 79
 - index adjustment 90
 - index limit 90
 - index missing 91
 - initialization failed 92
 - inlet valve fuse 84
 - inlet-valve missing 87
 - leak sensor open 77
 - leak sensor short 77
 - leak 76
 - missing pressure reading 83
 - motor drive power 86
 - pressure above upper limit 81
 - pressure below lower limit 82
 - pump configuration 83
 - pump head missing 89
 - remote timeout 75
 - selection valve failed 82
 - servo restart failed 89
 - shut-down 74
 - solvent zero counter 80
 - stroke length 91
 - synchronization lost 76
 - temperature limit exceeded 88
 - temperature out of range 88
 - timeout 74
 - valve failed 85
 - valve fuse 84
 - wait timeout 93
- exchanging
 - active inlet valve 114
 - active inlet valve cartridge 114
 - inlet valve 111
- multi channel gradient valve (MCGV) 131
- optional interface board 134
- outlet valve 117, 111
- pistons 111
- pump seals 111
- purge valve frit 119, 111
- purge valve 119, 111
- external contact
 - cable 164
- F**
- fan failed 79
- firmware
 - updates 136, 136
 - upgrade/downgrade 136
 - upgrade/downgrade 136
- flow connections 38
- frequency range 25
- G**
- general error messages 74
- gradient valve 131
- H**
- hints for successful use 48
- humidity 25
- hydraulic path 149
- I**
- index limit 90
- index adjustment 90
- index missing 91
- initialization failed 92
- initialization 15
- inlet-valve fuse 84
- inlet-valve missing 87
- installation, pump module 33
- installation
 - bench space 24
- instrument layout 20
- interface cables 35
- interfaces 170
- internet 189
- introduction to the pump 10
- L**
- LAN
 - cable 163
- leak rate test 102
 - running the test 104
- leak rate test
 - evaluating results 104
- leak sensor open 77
- leak sensor short 77
- leak 76
- line frequency 25
- line voltage 25
- lithium batteries 186
- M**
- maintenance
 - feedback 19
 - overview 111
 - pump head 126
 - replacing firmware 136, 136
- materials
 - bio-inert 17
- message
 - ignition without cover 79, 79
 - remote timeout 75
- missing pressure reading 83
- missing parts 28
- motor-drive power 86
- multi channel gradient valve (MCGV) 131

Index

N

- non-operating altitude 25
- non-operating temperature 25

O

- operating Altitude 25
- operating temperature 25
- operational hints, MCGV 57
- optimization
 - stack configuration 30
- optional interface board 134
- outlet valve 117, 140
- overview, pump 11

P

- packaging
 - damaged 28
- parts identification
 - cables 153
- parts
 - accessory kit 144
 - active inlet valve 142
 - bottle head assembly 148
 - damaged 28
 - hydraulic path 149
 - missing 28
 - outlet valve 140
 - pump head assembly 138
 - purge valve 141
 - solvent cabinet 146
 - starter kit 145
 - system tool kit 151
- physical specifications 25
- piston chamber 11
- piston 13, 48
- power supply indicator 66
- power consideration 22
- power consumption 25

- power cords 23
- power switch 34
- pressure above upper limit 81
- pressure below lower limit 82
- pressure pulsation 16, 60
- pressure range 59
- priming
 - with a pump 43, 56
 - with a syringe 56
- proportioning valve, high-speed 11
- PTFE frit 119
- pump error messages 80
- pump head assembly 122, 138
- pump head missing 89
- pump configuration 83
- pump head 126
- pump piston 48
- purge valve frit 48
- purge valve 119, 141

R

- radio interference 187
- reassembling the pump head 129
- remote
 - cable 158
- removing
 - pump head assembly 122
- repairs
 - replacing firmware 136, 136
- restart without cover 80
- RS-232C
 - cable 165
 - communication settings 177

S

- safety class I 182
- safety information
 - lithium batteries 186

- safety
 - general information 182
 - standards 25
 - symbols 184
- sapphire piston 13
- seal wash option 138
- seal wash 10, 11, 58
 - when to use 58
- seal, alternative material 59
- selection valve failed 82
- servo restart failed 89
- shut-down 74
- simple repairs 108
- site requirements
 - power cords 23
- snap fasteners 39
- solvent inlet filters 48
- solvent zero counter 80
- solvent cabinet 48, 146
- solvent filter
 - checking 113
 - replacing 113
- solvent filters
 - prevent blocking 49
- solvent information 51
- solvents, change 44
- sound emission 188
- special interfaces 175
- special settings
 - boot-resident 179
 - forced cold start 179
- specification
 - physical 25
- starter kit 145
- status indicator 67
- stroke length 91
- stroke volume 13, 16
- synchronization lost 76

Index

- system pressure test, evaluating results 99
- system pressure test 65, 97
 - running the test 99
- system setup and installation
 - optimizing stack configuration 30
- system tool kit 151
- WEEE directive 185
- weight 25

T

- temperature limit exceeded 88
- temperature out of range 88
- temperature sensor 76
- test functions 64
- test function
 - introduction 96
- timeout 74
- troubleshooting
 - error messages 73, 64
 - status indicators 64, 66

U

- unpacking the pump 28
- user interfaces 68

V

- vacuum degasser 10, 39
- valve failed 85
- valve frit 119
- valve fuse 84
- variable reluctance motor 13
- variable stroke volume 16
- voltage range 25

W

- wait timeout 93
- waste
 - electrical and electronic equipment 185

In This Book

This manual contains technical information about the Agilent 1260 Infinity Bio-inert Quaternary Pump G5611A. The manual describes the following:

- introduction,
- site requirements and specifications,
- installing the pump,
- using the pump,
- optimizing performance,
- troubleshooting and diagnostics,
- maintenance,
- parts and materials for maintenance,
- identifying cables,
- hardware information,
- appendix.

© Agilent Technologies 2011-2012

Printed in Germany
05/12



G5611-90000 Rev. C



Agilent Technologies