

# Air, water, CO<sub>2</sub> and ethylene

## Application Note

Energy & Fuels

### Authors

Agilent Technologies, Inc.

### Introduction

Gas chromatography with an Agilent PoraPLOT U column and Agilent 490 Micro GC separates air, water, carbon dioxide and ethylene in 60 seconds.



**Agilent Technologies**


## Conditions

Technique : Micro-GC  
GC-Channel : Agilent PoraPLOT U, 0.32 mm x 10 m  
Temperature : 50 °C  
Carrier Gas : Helium, 150 kPa  
Injector : 100 ms, 110 °C  
BackFlush : -  
Detector : chip TCD

Courtesy : Jim Luong and Rhonda Gras, Dow Chemical Canada

## Peak identification

1. air
2. water
3. carbon dioxide
4. ethylene

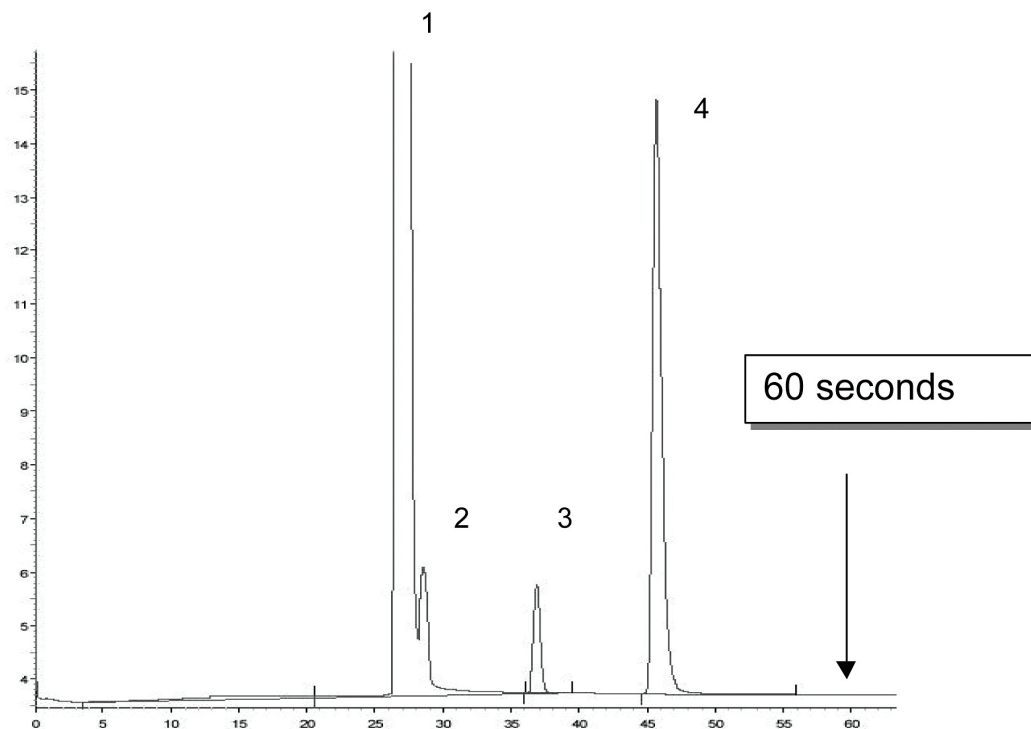

[www.agilent.com/chem](http://www.agilent.com/chem)

This information is subject to change without notice.

© Agilent Technologies, Inc. 2011

Printed in the USA

31 October, 2011

First published prior to 11 May, 2010

A02066



**Agilent Technologies**