

## Analysing Quaternary Ammonium Salts (QAS) in Seawater by Difficult Matrix Introduction (DMI)

Diane Nicholas

- **No sample preparation necessary**
- **Removes water under controlled conditions prior to analysis**
- **Retains the salt and involatiles within the microvial**

### Instrumentation

- ATAS Optic 2-200 programmable injector
- Agilent 5890 GC with 5971 MSD

### Principles

- Inject 2-3 uL of seawater into a microvial and place in fritted liner in Optic injector
- Vent the water, heating the injector and column oven to 100 °C
- Thermally desorb the sample at high temperature to degrade the QAS, under static flow conditions
- Transfer the analytes directly onto the GC column and analyse by GC-MS

### Chromatogram

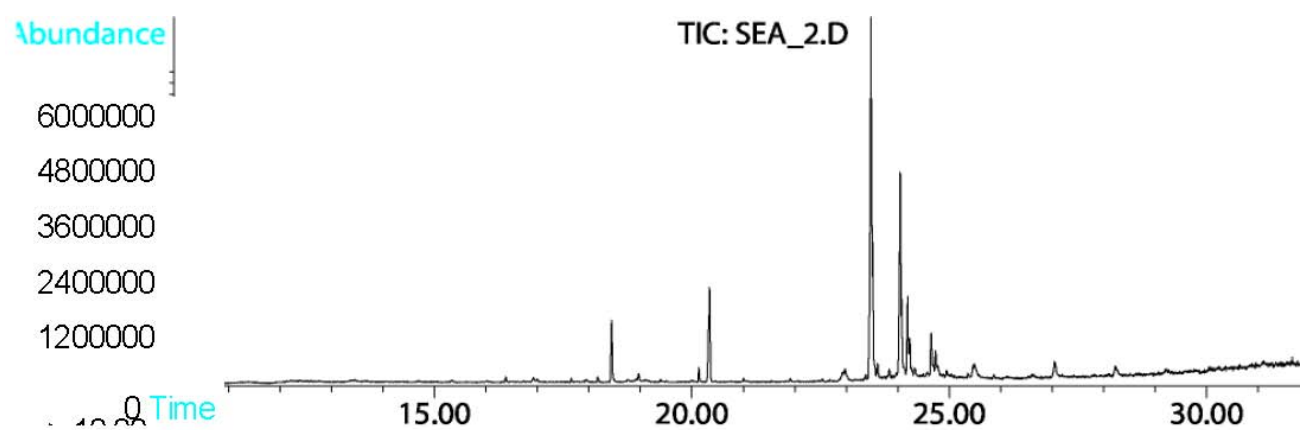


Figure 1: TIC of analysis of seawater by DMI

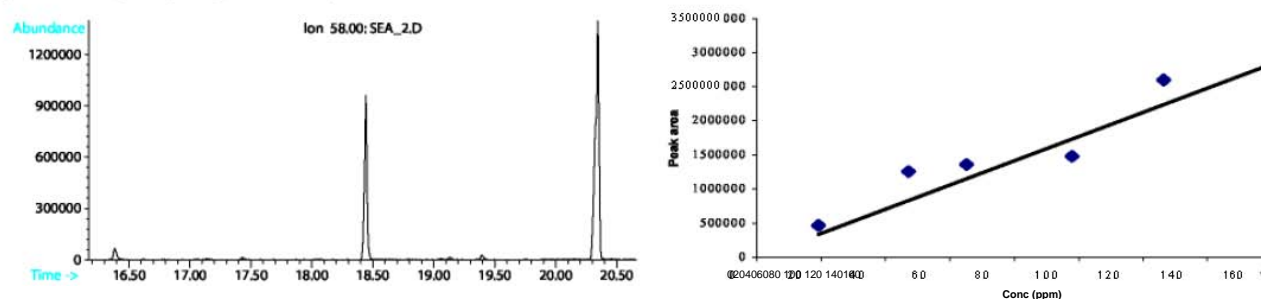


Figure 2: Selected ion for the quaternary ammonium salt Figure 3: Calibration with manual injection using the un-optimised method

We would like to thank John Sellers and Andy Dowding from AkzoNobel, UK for providing these samples

ATAS GL INTERNATIONAL B.V.  
De Sleutel 9, 5652 AS, Eindhoven, The Netherlands  
Tel. +31 (0)40 254 95 31 Fax. +31 (0)40 254 97 79  
E-mail: [info@atasgl.com](mailto:info@atasgl.com) Internet: [www.atasgl.com](http://www.atasgl.com)  
Web shop: [www.atasgl.com/shop](http://www.atasgl.com/shop)

Liner: ATAS Fritted  
Microvial: 6 mm  
Mode: Expert  
Gas Flows: Vent: 370 ml/min  
                  Split: 50 ml/min  
Initial temperature: 40 °C  
Ramp rate 1: 4 °C/s  
Final temperature 1: 100 °C  
Hold time 1: 1.5 mins  
Ramp rate: 16 °C/s  
Final temperature: 350 °C  
End time: 37.5 mins  
Vent pressure: 8 psi  
Vent time: 1.5 mins  
Split open time: 1.5 mins  
Desorption pressure: 0 psi  
Desorption time: 5 mins  
Transfer pressure: 7.4 psi  
Transfer time: 0.5 mins  
Initial pressure: 7.4 psi  
Final pressure: 22 psi

**GC Parameters:**


Column: DB5-MS 30m x 0.25mm i.d. x 0.25um film  
Initial temperature: 45 °C  
Ramp rate 1:  
Final temperature 1: 20 °C/min  
Hold time 1: 100 °C 4.5  
Ramp rate 2: mins 10  
Final temperature 2: °C/min 350  
°C

**MS Parameters:**

Acquisition mode: Scan  
Low mass: High 50 m/z  
mass: Sampling 400 m/z 2  
number: Threshold: 500  
Transfer line: 280 °C  
Solvent delay: 10 mins

**Appendix**

**Optic Parameters:**

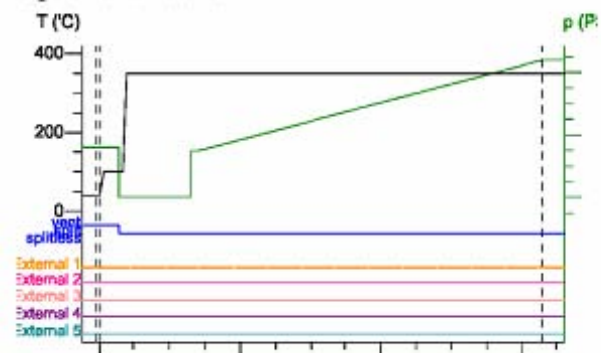

ATAS GL INTERNATIONAL B.V.  
De Sleutel 9, 5652 AS, Eindhoven, The Netherlands  
Tel. +31 (0)40 254 95 31 Fax. +31 (0)40 254 97 79  
E-mail: [info@atasgl.com](mailto:info@atasgl.com) Internet: [www.atasgl.com](http://www.atasgl.com)  
Web shop: [www.atasgl.com/shop](http://www.atasgl.com/shop)