

Clarity2Go

Analysis monitoring for smartphones

Clarity2Go is a mobile companion to Clarity Chromatography Station.

It enables to monitor selected parameters of running Sequence or Single Analysis conveniently from your mobile. Its use is **free** for all customers with Clarity station.

The monitoring works only if both (smartphone and PC with Clarity) have access to the internet.

The whole solution consists of three parts:

- Clarity Software (Full, minimal version 4.0.4) sending the data to an online server.
- Clarity2Go Server available at <http://clarity2go.dataapex.com>
- Clarity2Go application for smartphone or tablet.


Mobile application is currently available for Android (Google Play) and iOS (Appstore).



How to use Clarity2Go

In Clarity software

In Clarity Software on your PC use the *System / Clarity2Go* command to invoke the *Clarity2Go Configuration* dialog.

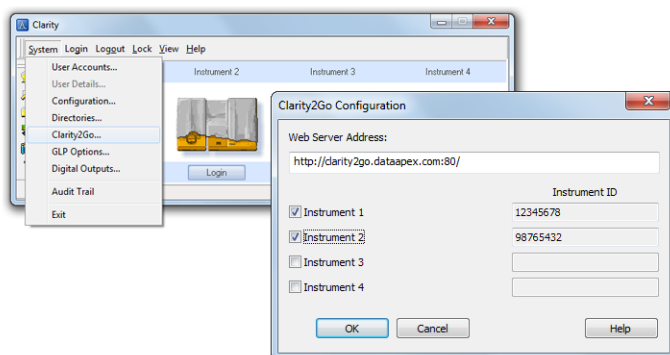


Select the instruments for monitoring, in the following example we have selected Instrument 1 and Instrument 2.

After checking the checkbox the Instrument is automatically registered on our server and you will receive unique ID for it. Use these Instrument IDs in Clarity2Go application on your smartphone/tablet.

In your smartphone

In application's Settings, enter the Instrument ID into "ID" field. Any number of Instruments can be added, not only from a single Clarity station but from different stations as well. Every Instrument can be enabled or disabled. Check "Show demo version" checkbox to display only simulated data. Switch between Instruments by swiping the screen left or right.

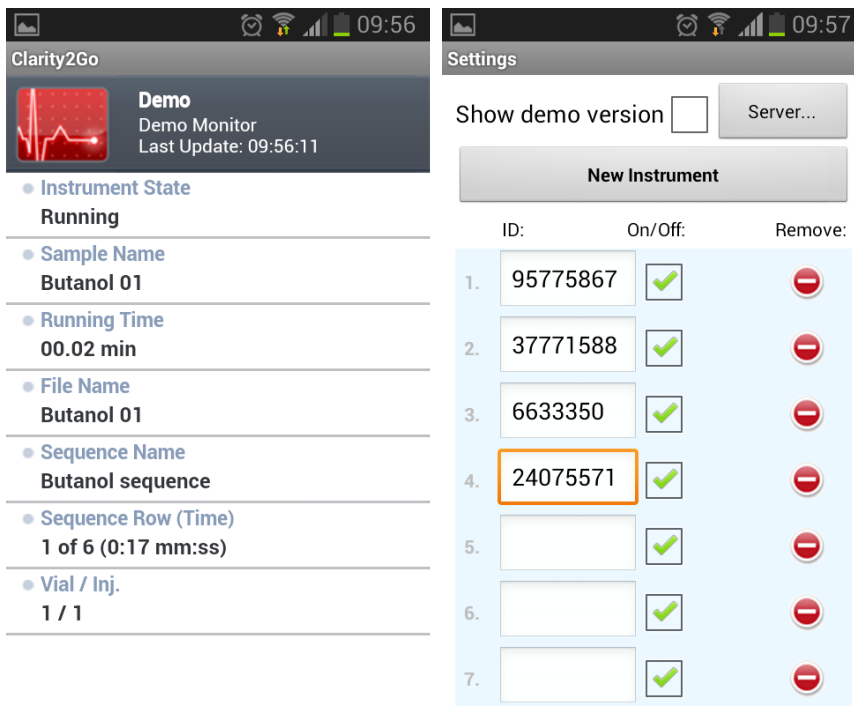

Clarity2Go was released with only few features and further development will follow the needs of Clarity users . If you have any suggestions or encounter some inconvenience, let us know at support@dataapex.com.

Help us make Clarity2Go better and give us your feedback!