

## Agilent G2709AA MassHunter StreamSelect B.06.00– Site Preparation Checklist

Thank you for purchasing an Agilent instrument. To get you started and to assure a successful and timely installation of your Agilent **software**, please refer to this specification or set of requirements.

Correct site preparation is the key first step in ensuring that your instruments and software systems operate reliably over an extended lifetime. This document is an **information guide and checklist** prepared for you that outlines the supplies, consumables, space and utility requirements for your equipment.

### Customer Responsibilities

**Make sure your site meets the following prior specifications before the installation date. For details, see specific sections within this checklist, including:**

- The computing environment and the necessary space is made available
- The number & location of electrical outlets for your computer systems and peripherals are planned
- That your site meets the software, hardware and networking specifications below
- Locate your sales order information, software authorization codes and/or software licenses/certificates
- The necessary software media, disks etc are available including upgrade/update disks
- That a suitable backup solution is identified for your software
- Availability of a system/network administrator as needed to connect to your intranet
- Please consult Other Requirements section below for other product-specific information.
- Complete Final Check: Software Site Preparation Tool. Please visit the following Agilent website to download the Software Installation Site Preparation Tool.

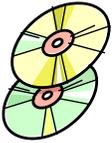
<http://www.chem.agilent.com/en-US/Support/Downloads/Utilities/SWSitePrepTool/Pages/Default.aspx>

This easy-to-use tool verifies that your PC meets the necessary hardware and software pre-requisites before installing the Agilent software.

**If Agilent is delivering installation and familiarization services, users of the instrument should be present throughout these services; otherwise, they will miss important operational, maintenance and safety information.**

### Important Customer Information

1. If you have questions or problems in providing anything described as a Customer Responsibility above, please contact your local Agilent or partner support/service organization for assistance prior to delivery. In addition, Agilent and/or its partners reserve the right to reschedule the installation dependent upon the readiness of your laboratory.
2. Should your site not be ready for whatever reasons, please contact Agilent as soon as possible to re-arrange any services that have been purchased.
3. Other optional services such as additional training, operational qualification (OQ) and consultation for user-specific applications may also be provided at the time of installation when ordered with the system, but should be contracted separately.

**Agilent G2709AA MassHunter StreamSelect B.06.00–  
Site Preparation Checklist**

**Software Specifications for  
Workstations, Clients, and Servers**
**Special Notes**

1. MassHunter StreamSelect software is an Add-On program that requires MassHunter Acquisition for 6400 Series QQQ software.

Specification Description	Minimum	Recommended (if applicable)	Comments (if applicable)
Operating system type(s), versions	Windows 7 - 64 Bit Professional		Requirements commensurate with QQQ Acquisition B.06.00
O/S .NET & other add-ons			Requirements commensurate with QQQ Acquisition B.06.00
Language settings/compatibility	U.S. English		Requirements commensurate with QQQ Acquisition B.06.00
Regional settings/compatibility	U.S. English		Requirements commensurate with QQQ Acquisition B.06.00
Account settings/privileges			Requirements commensurate with QQQ Acquisition B.06.00
Specific drivers			Requirements commensurate with QQQ Acquisition B.06.00
MassHunter Acquisition for 6400 Series QQQ	B.06.00 SP3	B.06.00 SP3	


**Computer Hardware Specifications for  
Workstations, Clients, and Servers**
**Special Notes**

1. Author to add special considerations or notes.

Specification Description	Minimum	Recommended (if applicable)	Comments (if applicable)
Processor type & speed	Intel Pentium Xeon W3520		Requirements commensurate with QQQ Acquisition B.06.00
Memory	4 GB	8 GB	Requirements commensurate with QQQ

**Agilent G2709AA MassHunter StreamSelect B.06.00-  
Site Preparation Checklist**

			Acquisition B.06.00
Internal Storage/devices/media			Requirements commensurate with QQQ Acquisition B.06.00
External Storage/devices/media			Requirements commensurate with QQQ Acquisition B.06.00
Video devices	1280 x 1024	1920 x 1200	Requirements commensurate with QQQ Acquisition B.06.00
Audio devices			Requirements commensurate with QQQ Acquisition B.06.00
Printing devices		HP LJ P3005D	Requirements commensurate with QQQ Acquisition B.06.00
Pointing devices			Requirements commensurate with QQQ Acquisition B.06.00



## Networking Specifications

### Special Notes

1. Author to add special considerations or notes.

Specification Description	Minimum	Recommended (if applicable)	Comments (if applicable)
Network type, bandwidth, speed, protocol etc	100 Mbps	1000 Mbps (1 Gbps)	Requirements commensurate with QQQ Acquisition B.06.00
Additional network or instrument devices/cards requirements			Requirements commensurate with QQQ Acquisition B.06.00



## Other Requirements

1. Important web links  
Microsoft Hardware Compatibility Lists  
<http://www.microsoft.com/whdc/hcl/default.msp>

**Agilent G2709AA MassHunter StreamSelect B.06.00-  
Site Preparation Checklist**

Links to specific O/S fixes, updates needed

<http://support.microsoft.com/>

2. Refer to the **WINDOWS 7 PROFESSIONAL (64-BIT) FOR MASSHUNTER WORKSTATION SETUP GUIDE** (p/n G3336-90016) for details on MassHunter specific Windows Configuration requirements:

[http://www.chem.agilent.com/Library/usermanuals/Public/G3336-90016\\_Windows7\\_Setup.pdf](http://www.chem.agilent.com/Library/usermanuals/Public/G3336-90016_Windows7_Setup.pdf)

## Agilent G2709AA MassHunter StreamSelect B.06.00- Site Preparation Checklist

### Important Customer Web Links

- For additional information about our solutions, please visit our web site at <http://www.chem.agilent.com/en-US/Pages/HomePage.aspx>
- Need to get information on your product?  
Literature Library - <http://www.agilent.com/chem/library>
- Need to know more?  
Customer Education - <http://www.agilent.com/chem/education>
- Need technical support, FAQs? - <http://www.agilent.com/chem/techsupp>
- Need supplies? - <http://www.agilent.com/chem/supplies>

Document part number: G2709-90002