



Controlling the Agilent 1260 Infinity/1290 Infinity II Multisampler (G7167A/B) in Waters Empower 3 environment using the Agilent Instrument Control Framework

Technical Note

Technical Guide for the configuration and use of the Agilent 1290 Infinity II Multisampler (G7167B) and the Agilent 1260 Infinity Multisampler (G7167A) with Waters Empower 3 using the Agilent Instrument Control Framework (ICF).

Contents

Introduction 2

Prerequisites 2

Defining and Assigning Plates in Waters Empower 3

Multisampler Method Setup 8

Start an Analysis in Empower 3 11

Appendices 13

A - Configuring an Agilent LC in Empower 13

B - Importing Agilent Well Plate Definitions 19

C - Manual Set Up of Well Plates and Example Configuration Layouts for Standard Agilent Well Plates 23

D - Optional Multisampler Method Setup Parameters 26



Agilent Technologies

Introduction

The Agilent Instrument Control Framework (ICF) is a software component that manages the communication between the chromatography data system (CDS) and the instrument driver. This component makes it easier and faster for CDS software vendors to implement control of Agilent LC, CE and GC instruments in their data systems or workstations. Based on new standard instrument drivers from Agilent, ICF eliminates much of the delay and effort of using low-level instrument control codes and the need of software developers to write their own native drivers.

Waters Instrument Control Software (ICS) is Waters Corporation adoption of the Agilent Instrument Control Framework for their data systems.

This guide is designed to assist you in setting up a Multisampler based method. Additional appendices explain the process to perform additional, sometimes optional, configuration steps to use the Multisampler.

NOTE

For more information on support, refer to the Waters Release Note.

Prerequisites

Instrument Prerequisites

- Agilent 1290 Infinity II Multisampler (G7167B) or Agilent 1260 Infinity Multisampler (G7167A) with Firmware Rev. D.06.60 or higher.
- Other Agilent LC modules in the LC system *must meet or exceed* the minimum firmware requirements specified by the 3rd-party CDS software vendor and meets Agilent's firmware set/firmware interoperability requirements.
http://www.chem.agilent.com/_layouts/agilent/downloadFirmware.aspx?whid=79809

Software Prerequisites:

The operating system and Empower prerequisites are documented in the Waters Empower Instrument Control Release Notes for Agilent Instrument Control.

At minimum the following components are required:

- Waters Instrument Control version (ICS): ICS 2.1 Hotfix #716004653 including
 - Agilent ICF Version A.02.03 DU1 HF2
 - Agilent ICF-LC Driver Version A.02.11 SP1 [72]

NOTE

Ensure that all prerequisites are met before proceeding to the next step. The component versions can be verified in the operating systems Control Panel/Program and Features.

Account set up

- For installation, please ensure the account being used has Administrator rights and privileges.


Defining and Assigning Plates in Waters Empower

NOTE


In order to use Agilent well plates with Empower and the Multisamplers, well plate definitions for these plates must exist already in Empower. If they are not present, then you will need to import at least one Agilent plate definition file for the first time. See Appendix “[B - Importing Agilent Well Plate Definitions](#)” on page 19 for importing Agilent well plate definitions.

Proceed with the following steps to define the Agilent well plate within Empower.

- 1 Go to the main Edit menu then click **Plates**.

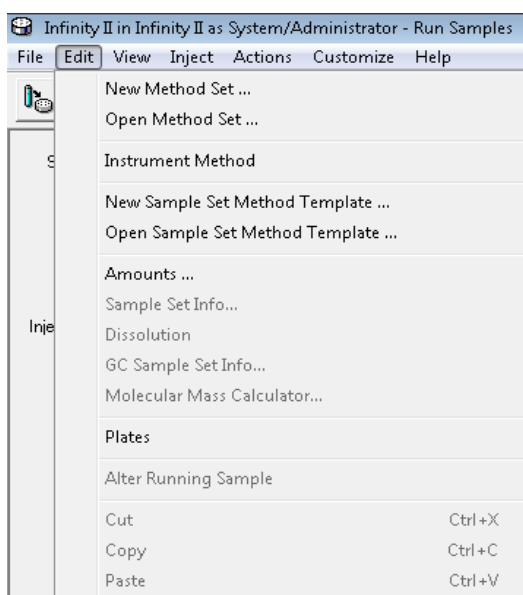

- 2 Within a row of the Plate Type Name column click the drop down arrow and select the plate of interest. If the plate type is not present it need either to be imported (Appendix “B - Importing Agilent Well Plate Definitions” on page 19) or to be created manually (Appendix “C - Manual Set Up of Well Plates and Example Configuration Layouts for Standard Agilent Well Plates” on page 23).

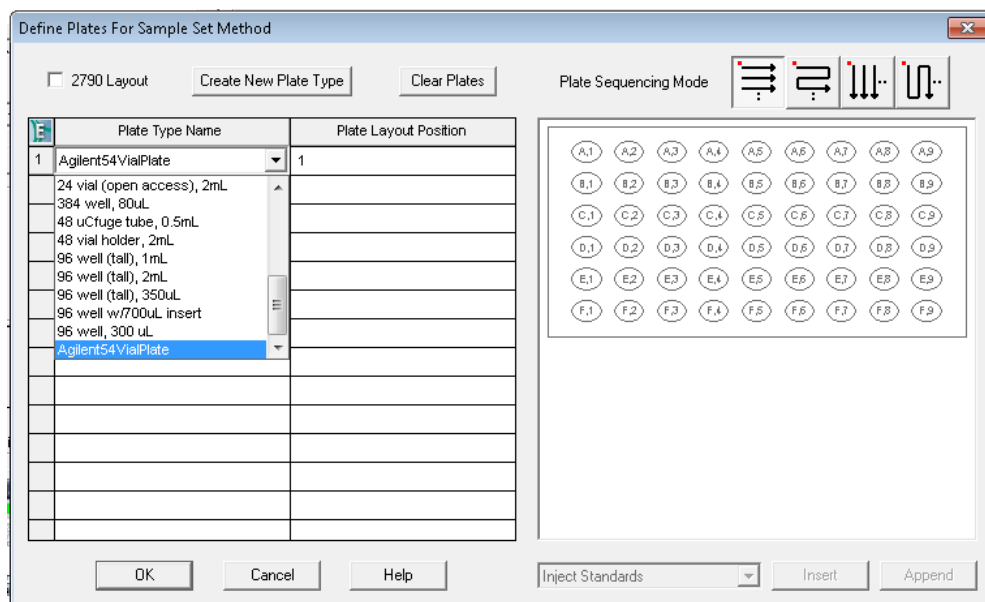

NOTE

The selected plates defined in the table above must be the same type and position as defined in the Tray and Plate Configuration dialog of the Multisampler graphical user interface (refer to step 8 on page 6). Otherwise the mismatch prevents starting the analysis.


- 3 Ensure that the just defined well/vial plates for Empower matching the assigned well/vial plates for the Multisamplers.
- 4 Go to the **Instrument Status** display.




- 5 Right mouse click on the **Instrument Status** of the Multisampler to display the context-sensitive menu.

NOTE

Perform these steps only if you have changed the drawer configuration from last time.

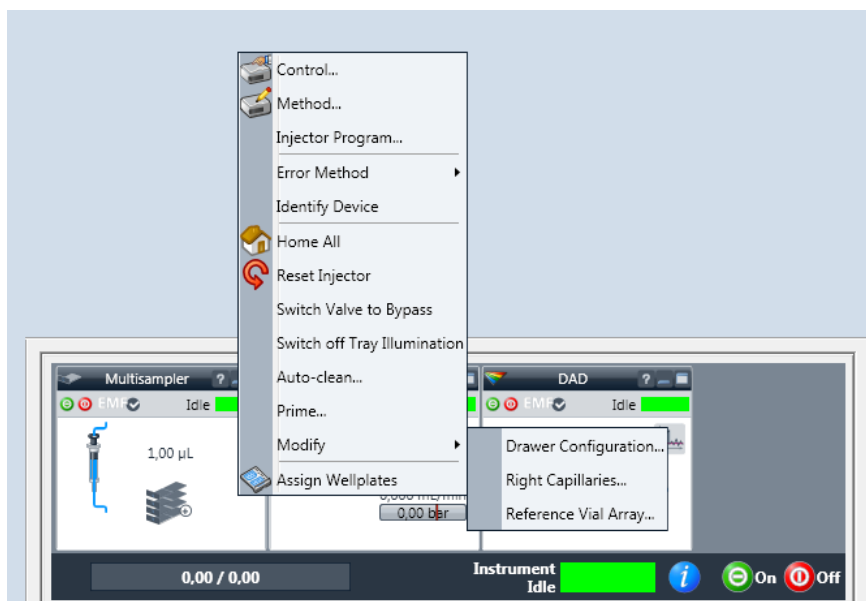


- 6 Click **Modify** from the drop-down menu list and select the **Drawer Configuration** option.
- 7 Click the **Start** button.

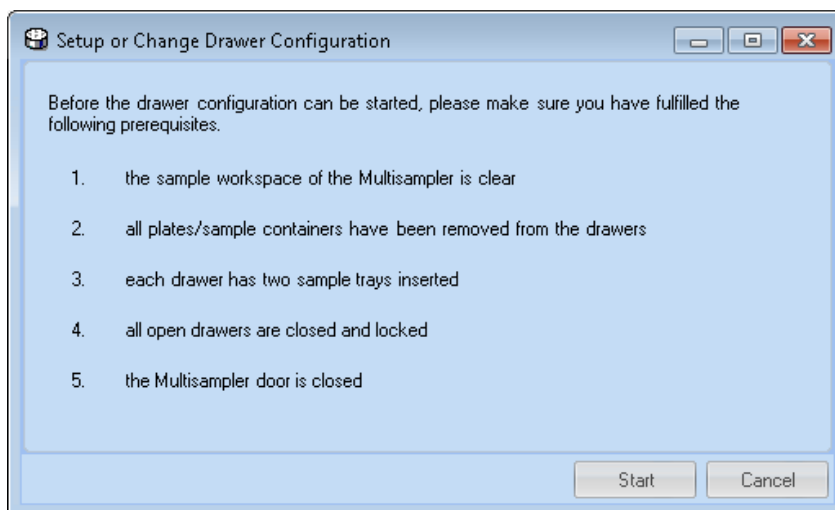

The Multisampler will run through its configuration procedure – this may take some minutes.

- 8 Right mouse click on Multisampler **Instrument Status** and click **Assign Wellplates**.

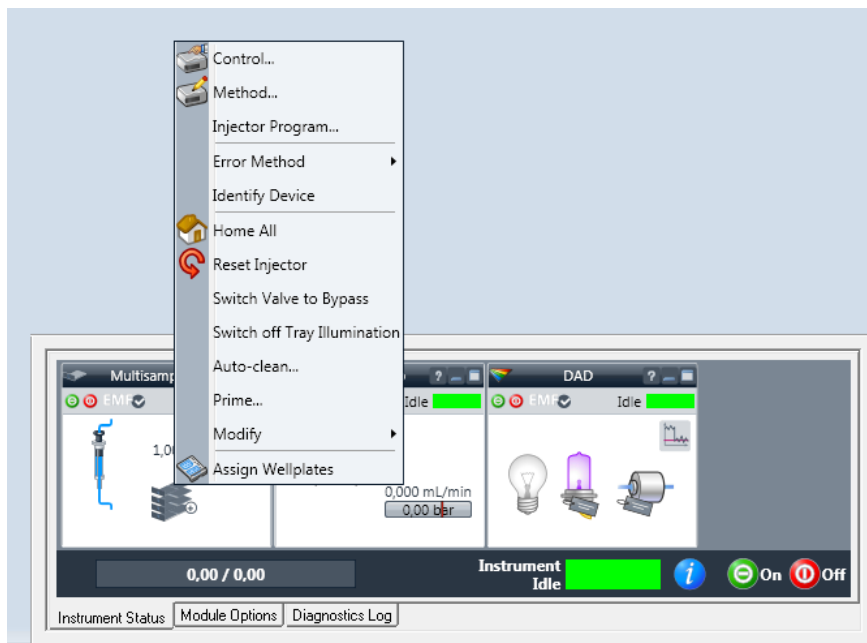

This accesses the Tray and Plate Configuration of the Multisampler.


- 9 Select Tray and Plate Configuration of P1 up to PX (depending of the configured drawers) and select the well/vial plate matching the type previously set in defined plate within Empower. Then click **OK**.

NOTE


If the correct drawer configuration is not directly displayed in the LC Status Window after the **Drawer Configuration Change** either scan the instrument again or reboot the LAC/E box.

NOTE

The selected plates here must be the same type as in **Define Plates** configured within Empower.

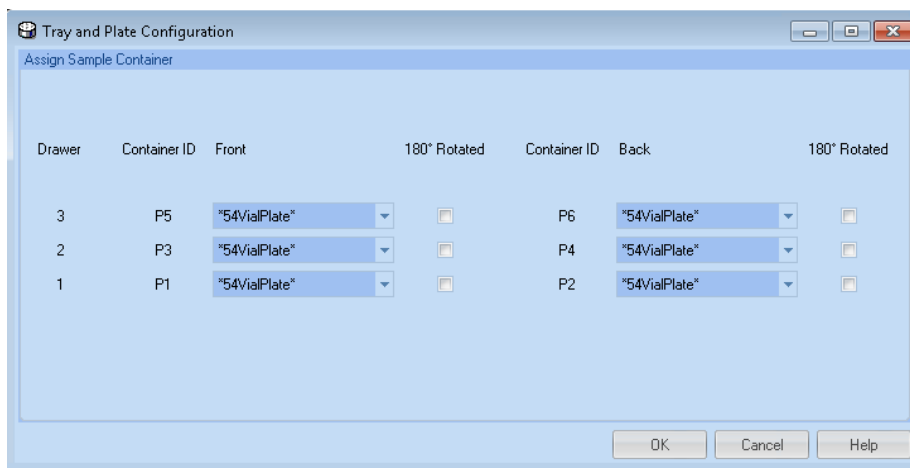

- 10 Return to the Multisampler **Instrument Status**, click on **Modify**, click on **Right Capillaries**, enter the appropriate settings for the Loop Capillary and the Seat Capillary, and then click **Assign**.

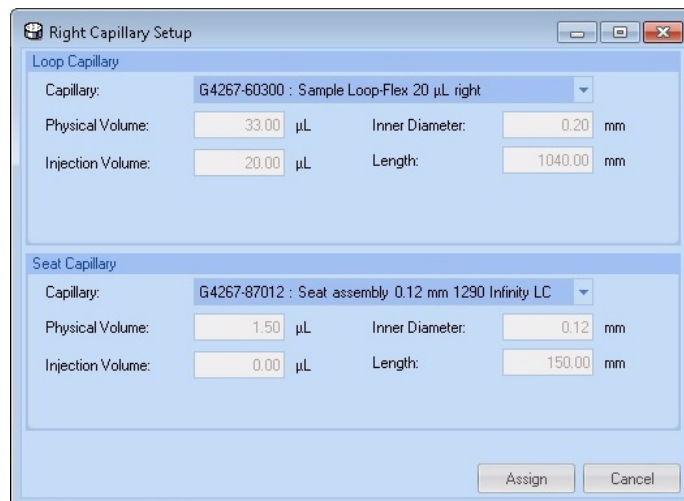

The image shows a software dialog box titled "Right Capillary Setup". It contains two sections: "Loop Capillary" and "Seat Capillary". Each section has a dropdown menu for "Capillary" and three input fields for "Physical Volume", "Inner Diameter", and "Length". The "Loop Capillary" section is currently selected.

Section	Capillary	Physical Volume (μL)	Inner Diameter (mm)	Length (mm)
Loop Capillary	G4267-60300 : Sample Loop-Flex 20 μL right	33.00	0.20	1040.00
Seat Capillary	G4267-87012 : Seat assembly 0.12 mm 1290 Infinity LC	1.50	0.12	150.00

At the bottom right of the dialog box are two buttons: "Assign" and "Cancel".


Multisampler Method Setup

This section describes the setup of the Multisamplers method and its various parameters once the LC system has been setup and configured in Empower.

NOTE

It assumes you have completed the necessary steps to connect the modules and configured them within Empower. See Appendix [“A - Configuring an Agilent LC in Empower”](#) on page 13 for how to configure a Multisampler system.

It also assumes you have already imported, defined and assigned the required Agilent well plate definition. See the previous section.

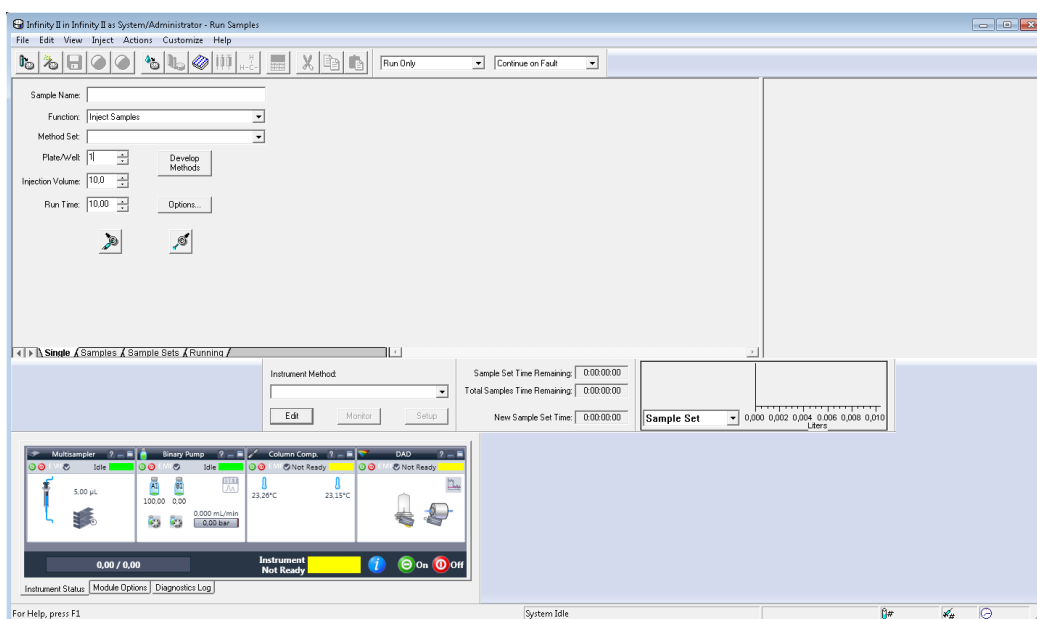


The Agilent Instrument Status screen and ICF software under Empower is shown above with the Multisampler configured at the bottom of the screen.

The instrument method and the method set can be created in Empower by accessing the Empower Run Samples screen. The Instrument method is set up and saved through the Edit/Instrument Method dialog.


The Multisampler dialog screens for setting appropriate method parameters are shown below ([“Edit Instrument Method”](#) on page 9). Additional/optional Multisampler method setup steps are provided in Appendix [“D - Optional Multisampler Method Setup Parameters”](#) on page 26.

Edit Instrument Method


- 1 Go to **Edit** and click **Instrument Method**.

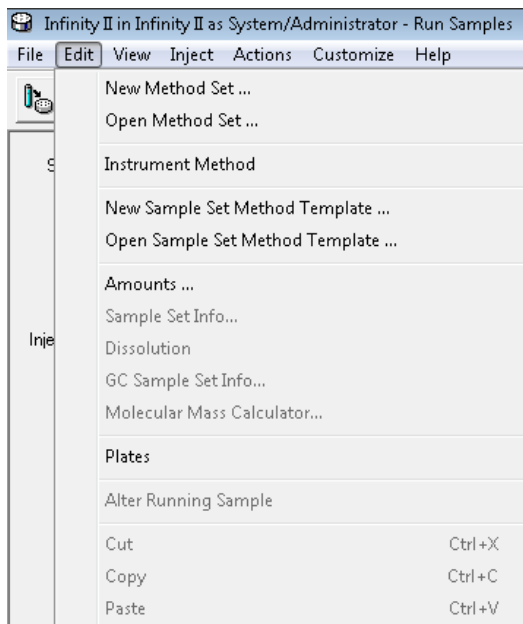

- 2 Click the Multisampler tab. The Instrument Method Editor is displayed with the default settings and the available Advanced features. Enter the appropriate parameters in these fields and click the save icon to save the method.

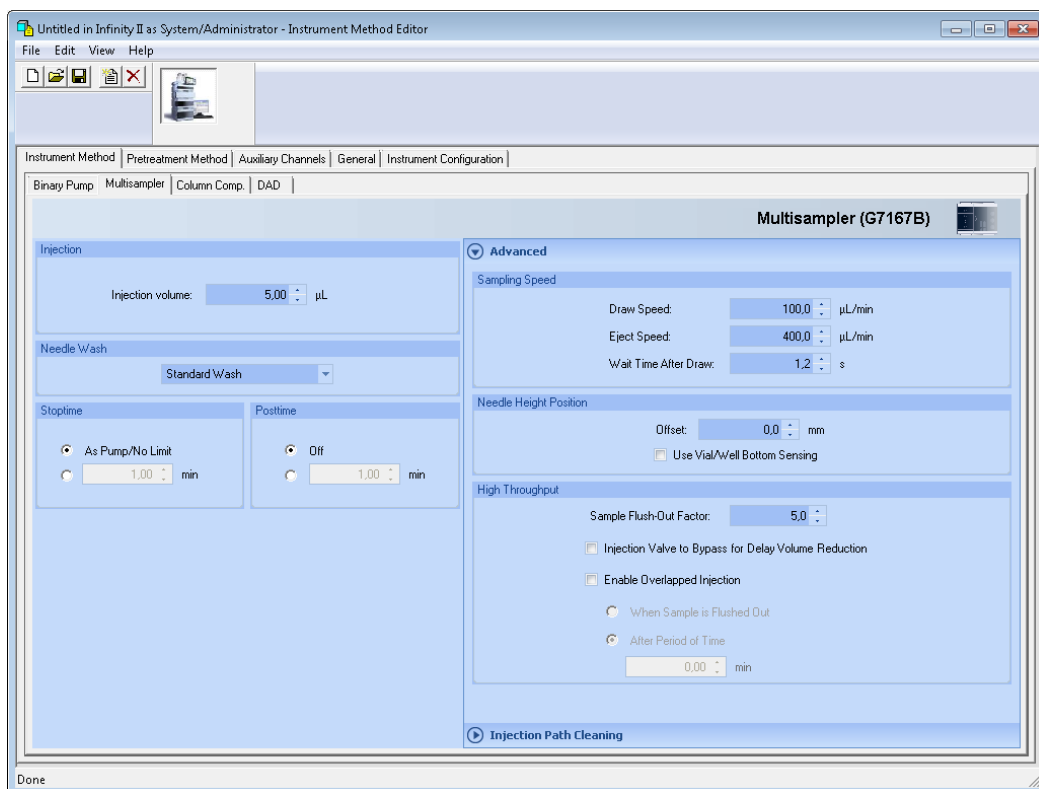

- 3 Enter the Name, Method Comments (optional), click **Save** and close the Instrument Method Editor.

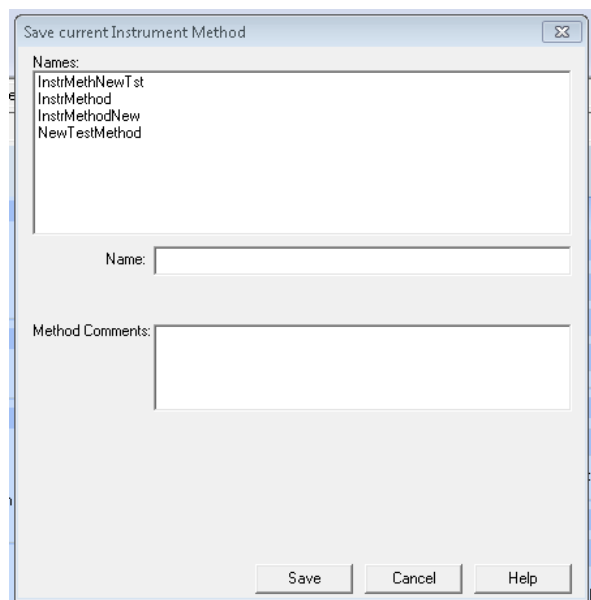

The image shows a dialog box titled "Save current Instrument Method" with a close button (X) in the top right corner. Inside the dialog, there is a section labeled "Names:" containing a list box with four entries: "InstrMethNewTst", "InstrMethod", "InstrMethodNew", and "NewTestMethod". Below the list box, there is a "Name:" label followed by a text input field. Underneath that, there is a "Method Comments:" label followed by a larger text input area. At the bottom of the dialog, there are three buttons: "Save", "Cancel", and "Help".


Start an Analysis in Empower 3

NOTE

Ensure the Agilent well plate is positioned in the lowest drawer position of the Multisampler! All positions in the Sample Hotel must be filled either with dummies or drawers. The drawers must be installed from bottom to top.

Create a Single Injection/Run

To create a single injection/run, follow the steps below for creating a Method Set. This can be performed from the main Empower Run Samples screen.

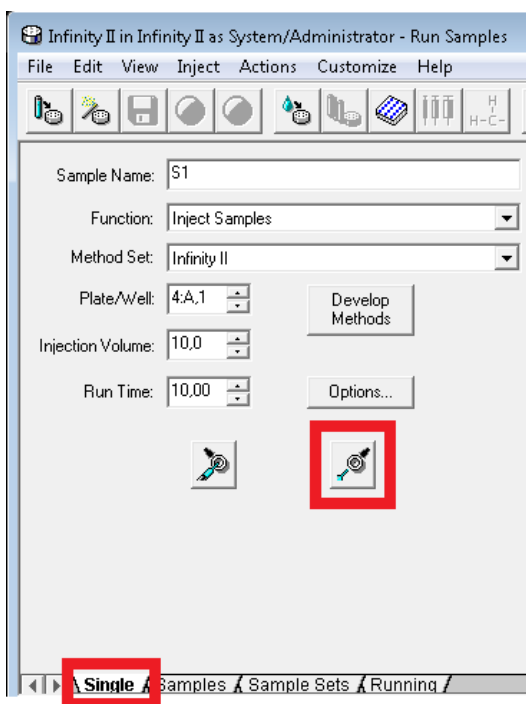

- 1 Ensure you are on the Single tab as outlined above and enter appropriate values including Sample Name, Function, Plate/Well, Injection Volume and Run Time.
- 2 When entering the info for the Plate/Well, select the appropriate position of the sample or vial. For example "4:A,1" refers to a vial in sample position A1 being used in Plate 4 (P4).


NOTE

In Empower 3, "4:A,1" corresponds to "P4-A-1" in the Multisampler user interface. For more information on plate configurations, refer to the Appendix "[C - Manual Set Up of Well Plates and Example Configuration Layouts for Standard Agilent Well Plates](#)" on page 23 and the Agilent 1200 Infinity Series Multisamplers User Manual.

- 3 Click the **Inject** Icon (outlined above) to start the analysis.

Create a Series of Injections

To create a series of injections, follow the steps below for creating a Method Set. This can be performed from the main Empower Run Samples screen.

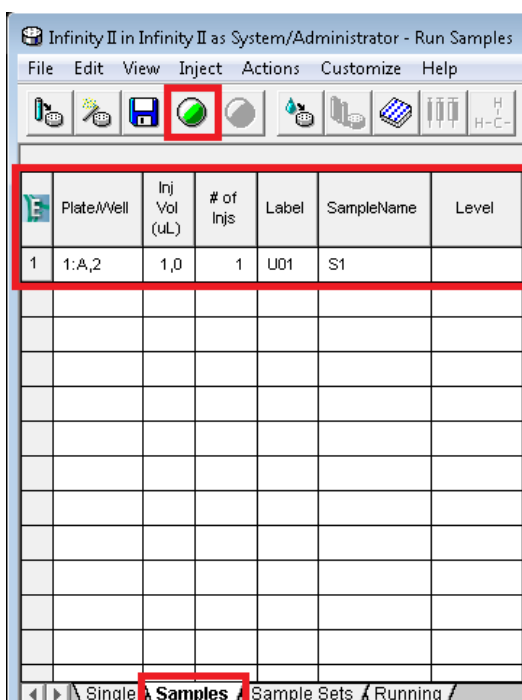

- 1 Select File/New Sample Set Method/Using Sample Set Wizard from the top level menu and follow the steps provided on screen. In the example above in the Samples tab, the Sample Set Method table contains a series of parameters for making multiple injections.
- 2 Ensure there are values for Plate/Vial position, injection volume, the number of injections per vial, Label and Sample Name in the Sample Set Method table. In this example above, "1:A,2" refers to sample position A2 being used in Plate 1 (P1).
- 3 Click on the green **Run** icon (outlined above) to start the analysis.

NOTE

In Empower 3, “1:A,2” corresponds to “P1-A-2” in the Multisampler user interface. For more information on plate configurations, refer to the Appendix **“C - Manual Set Up of Well Plates and Example Configuration Layouts for Standard Agilent Well Plates”** on page 23 and the Agilent 1200 Infinity Series Multisamplers User manual.

Appendices

A - Configuring an Agilent LC in Empower


NOTE

These configuration steps have to be followed whenever a new module is added or removed. Previously, the old configuration has to be deleted from the DHCP server configuration. Then the Empower software has to be shut down and the LAN connection to the module has to be switched off and on again. When the LAN connection has been restored the new configuration procedure can be started.

HINT


Ensure that the module hosting the LAN connection is enabled to use DHCP. This can be achieved by setting up the DIP switches of the module in a specific order. Please refer to the manual of the module hosting the communication.

- 1 Access the Configuration Manager as System Administrator as shown below.

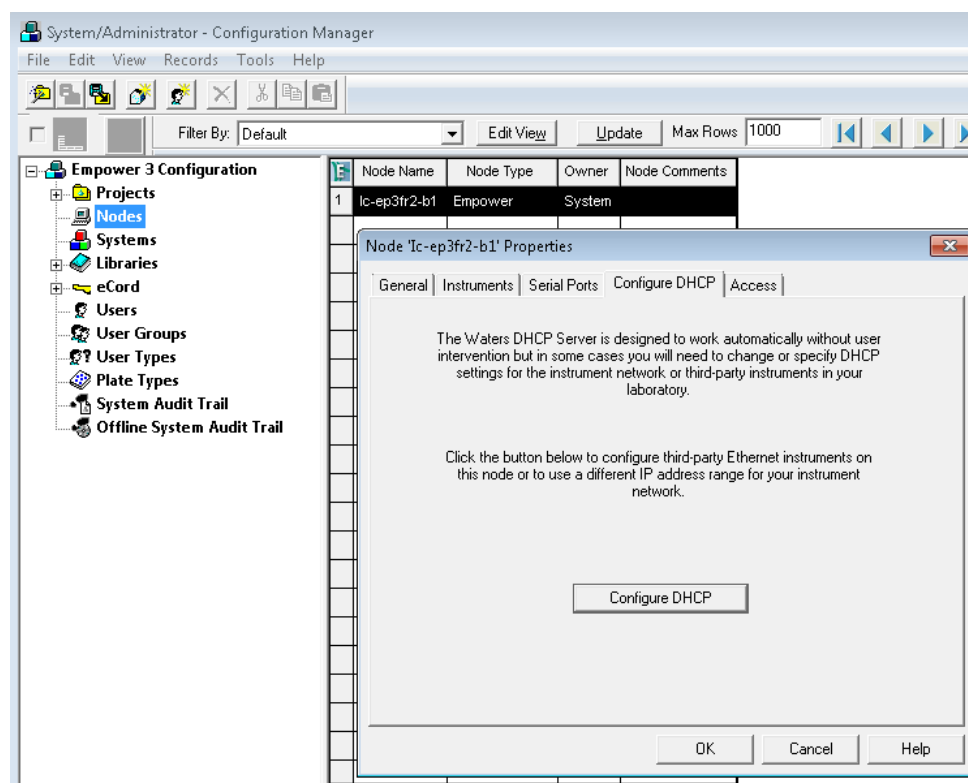


- 2 Go to Properties of the Empower Nodes.
- 3 Click **Configure DHCP** button on the Configure DHCP tab.

- 4 Check the IP Address and the MAC Address of the instrument. If the parameters are correct click **OK**. To add a new instrument, Click **Add...**

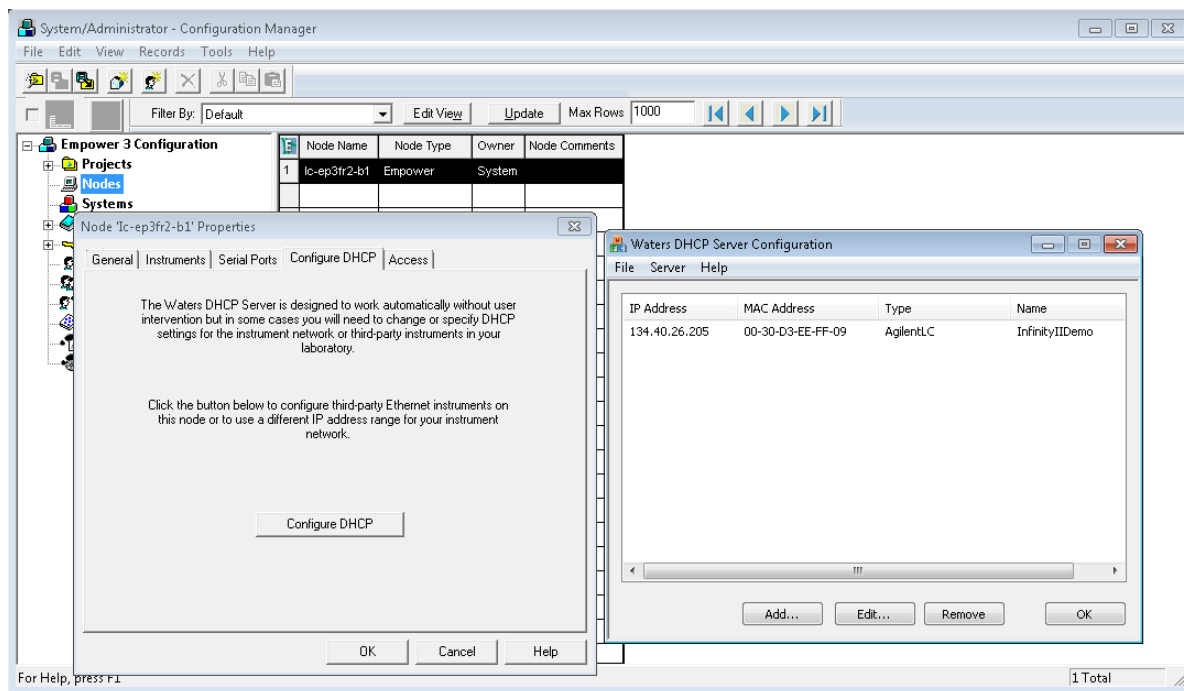


- 5 Use **Add IP Address** screen to set the Instrument Type (select **Agilent LC** from the drop-down list) and enter a unique name for the system in the Serial Number/Unique Name field then click **OK**.

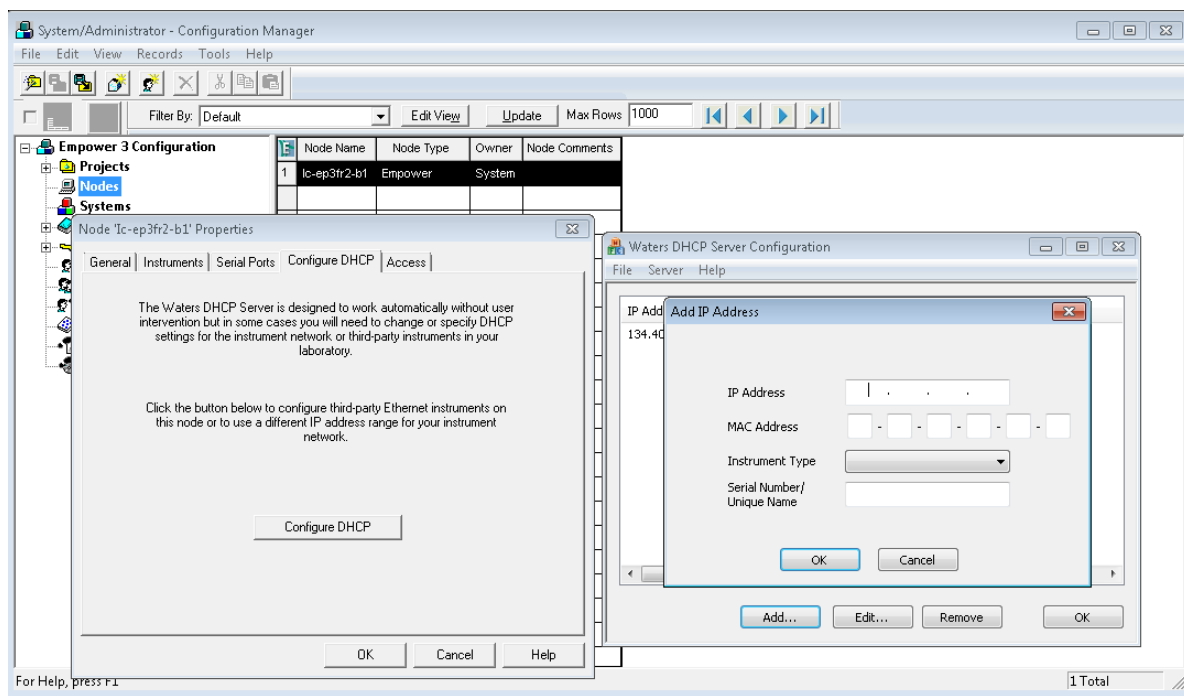

The unique name defined here will then be used as the address in the Node/Properties/Instrument tab.

- 6 From the main menu, navigate to the **Node/Properties/Instrument** tab and click the **Scan Instruments** button to check you have a connected instrument.

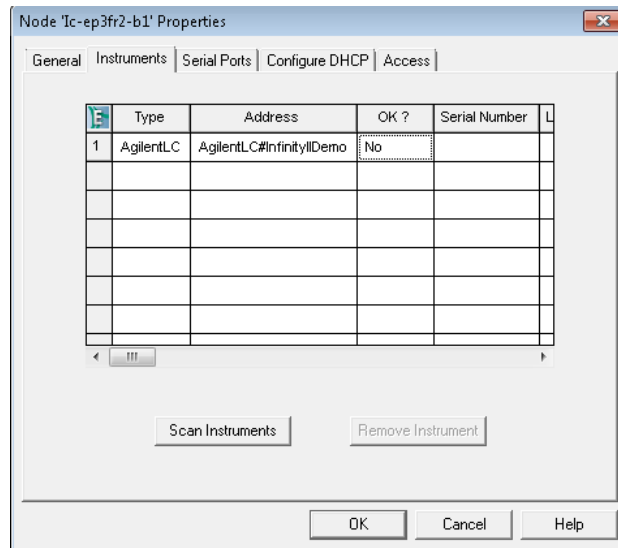


From the example LC Status Windows information view (accessible via  in the LC status window), you can review the list of connected LC modules and related information as shown.

Vendor	Name	Part Number	Serialnumber	Firmware Revision	Connection Info	Driver Version	Additional Information
Agilent	Quat. Pump	G7104A	DEBAX00158	B.06.73 [0002]	134.40.26.205	A.02.11 SI 517.05	
Agilent	Binary Pump	G7120A	DEBA200149	B.06.73 [0001]	134.40.26.205	A.02.11 SI 517.05	
Agilent	Multisampler	G71678	DEBAS00205	D.06.72 [0002]	134.40.26.205	A.02.11 SI 517.05	
Agilent	Column Comp.	G71168	DEBA400197	D.06.72 [0002]	134.40.26.205	A.02.11 SI 517.05	Slave Firmware: C.06.72 [0001]
Agilent	RID	G1362A	DEAA601801	A.06.54 [002]	134.40.26.205	A.02.11 SI 517.05	
Agilent	DAD	G71178	DEBAW00171	D.06.72 [0002]	134.40.26.205	A.02.11 SI 517.05	Access Point


Configuration as of last online instrument connection

The next series of steps enables the connected system to be configured and visible in Empower.


- 1 Click **File, New** and click **Chromatographic System** to make the new LC system accessible to Empower.

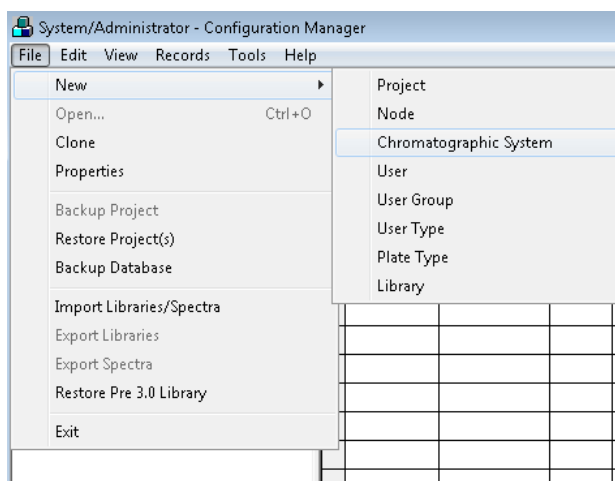


- 2 Select **Create New System** to initiate the wizard and click **Next >**.

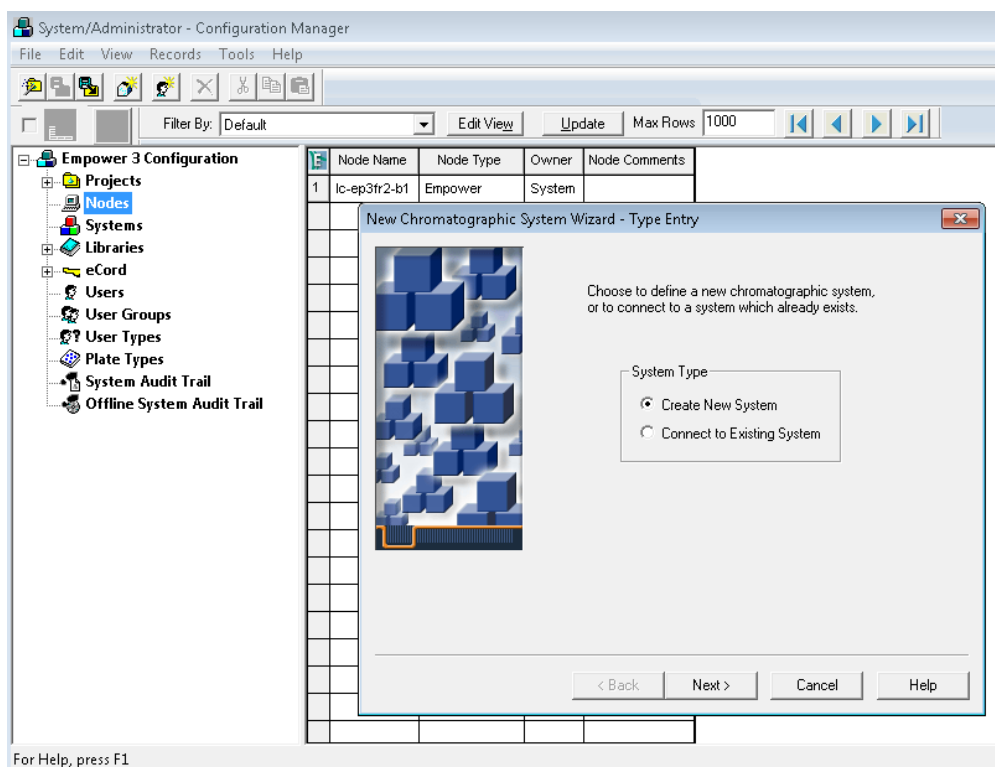


- 3 Drag and drop the desired instrument from the Available Instruments list to New System Instruments list and click **Next >**.

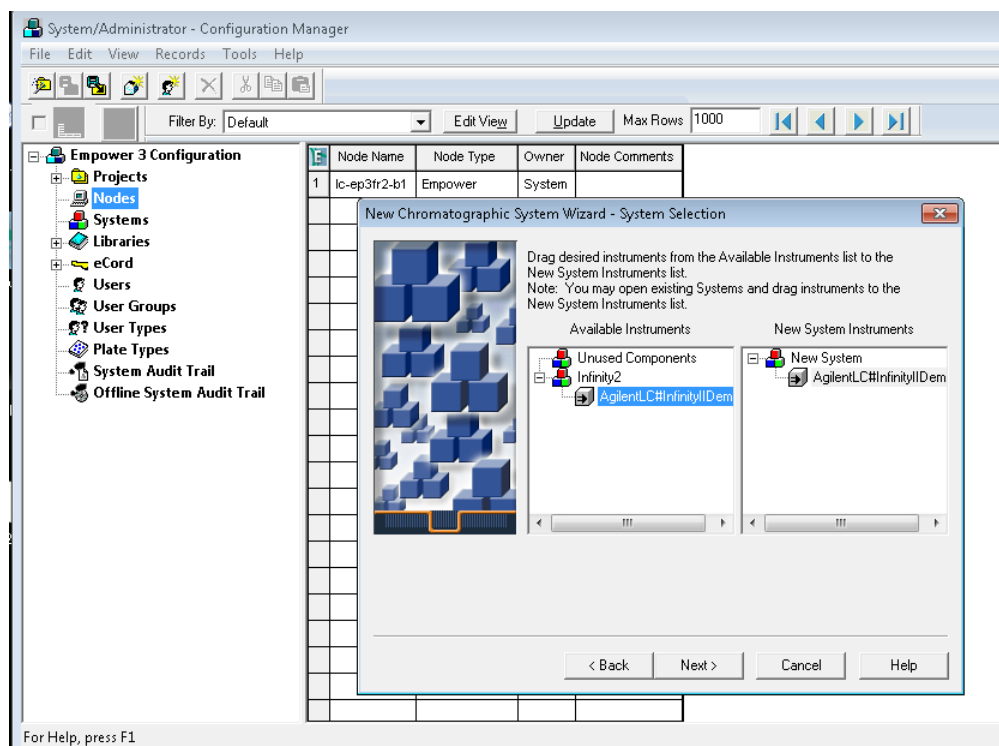


- 4 In the Allowed Access dialog box, select the radio button **Owner, Group and World**. Click **Next >**.

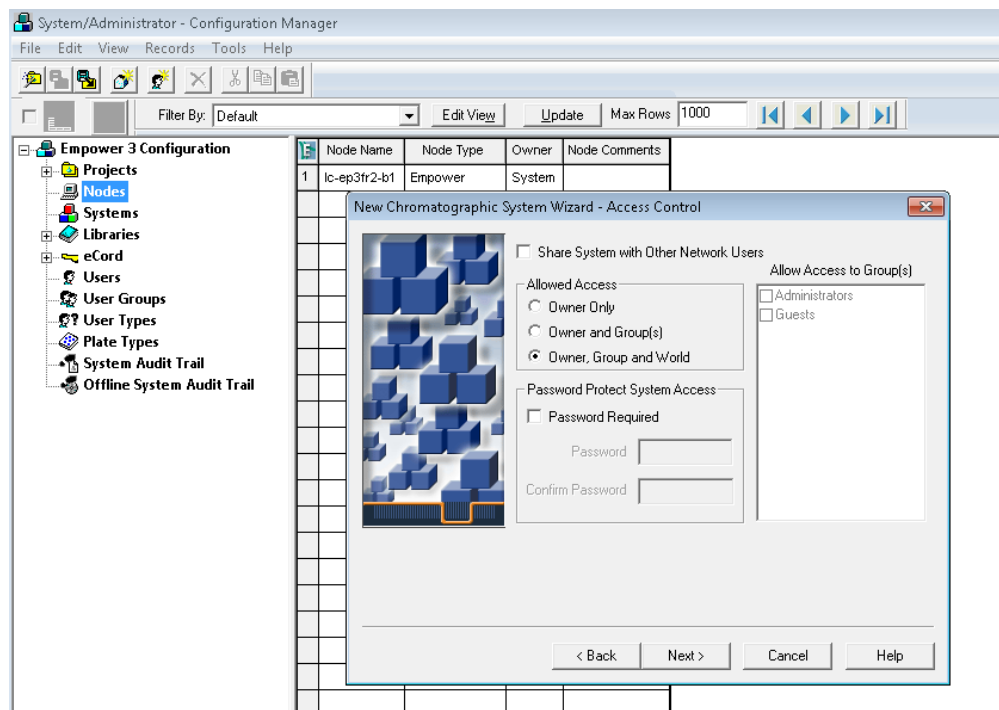

- 5 Enter the **System Name** and **System Comment** (optional) then click **Finish**.

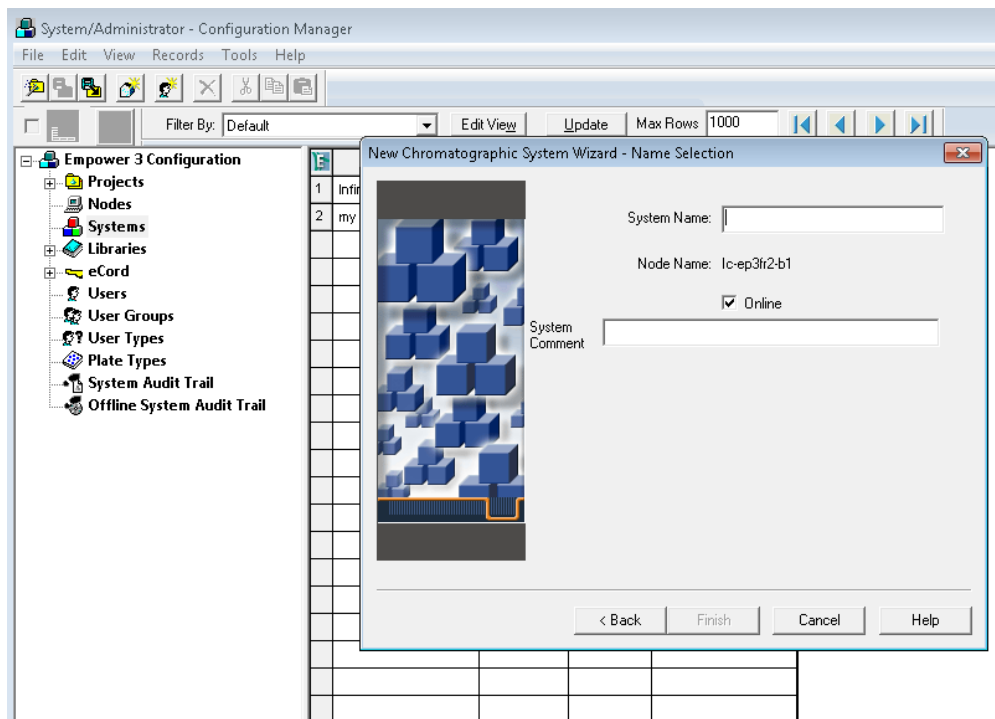

After the configuration procedure, the LC system should appear in the list of systems as shown in the example below.

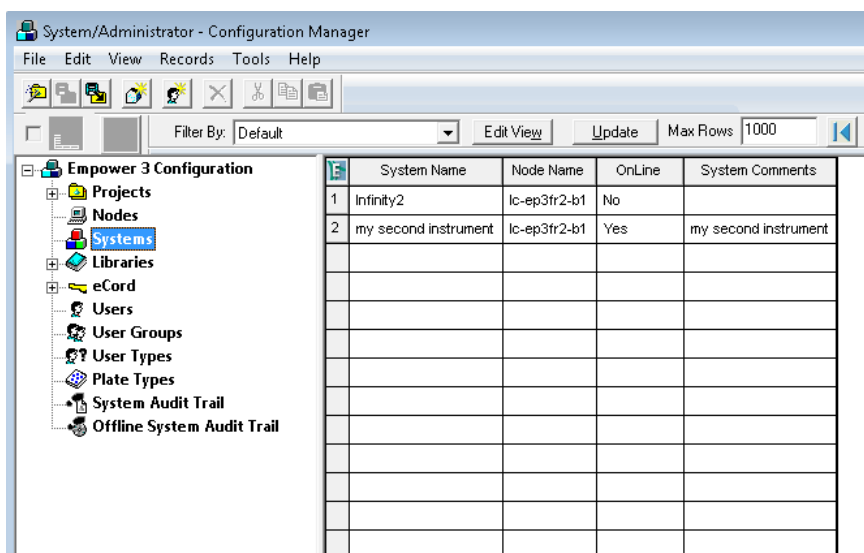

NOTE

You can only run an analysis when OnLine is set as Yes in the table above.


B - Importing Agilent Well Plate Definitions

In order to use Agilent well plates in conjunction with Empower 3 and a Multisampler, well plate definitions must exist already. These steps outline the procedure for importing new Agilent well plate definition files into Empower.

NOTE

If the Agilent well plate types are not present in the list, it will be necessary to search and import at least the Agilent 54-well plate definition .txt file as this is not installed automatically by Empower 3. Also, importing of plate definition .txt files can only be performed one at a time.


From the Configuration Manager window, select **Plate Types** in the left navigation pane to display the current list of well plate types available that can be used within Empower. Determine if the Agilent well plate types are present. In the example, the plates do not exist so follow the steps below.

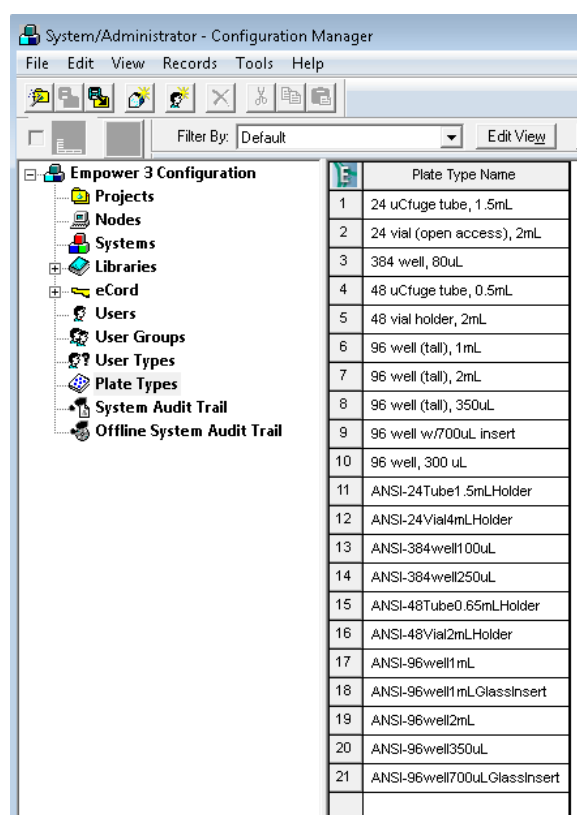


Appendices

B - Importing Agilent Well Plate Definitions


- 1 Right click on an empty row of the **Plate Type Name** table to select the list of plate menu functions.

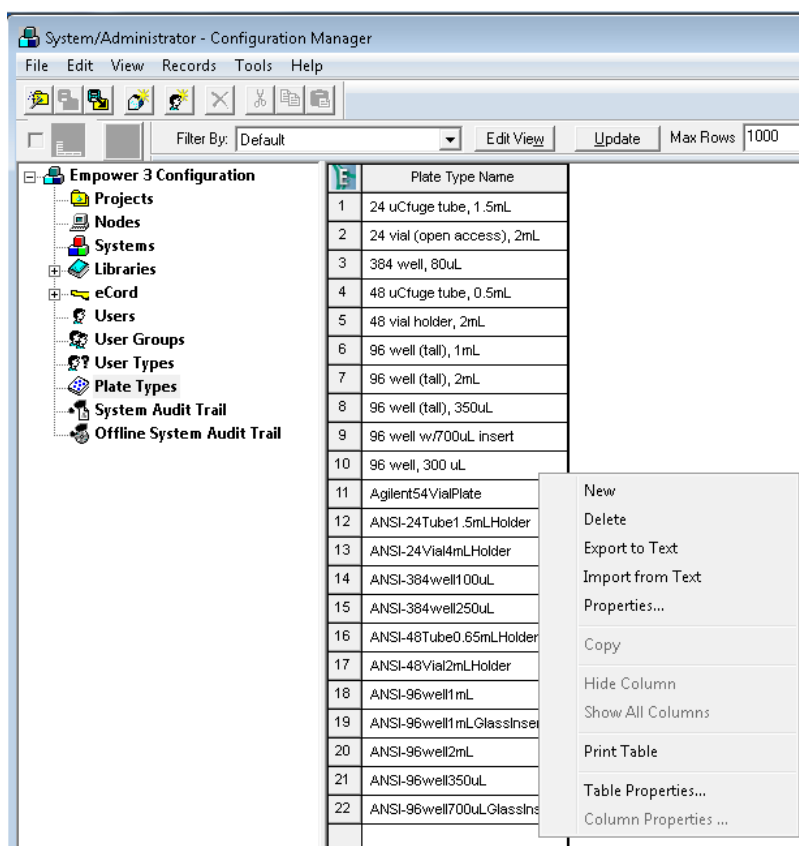


- 2 Select **Import from Text** from the menu list.

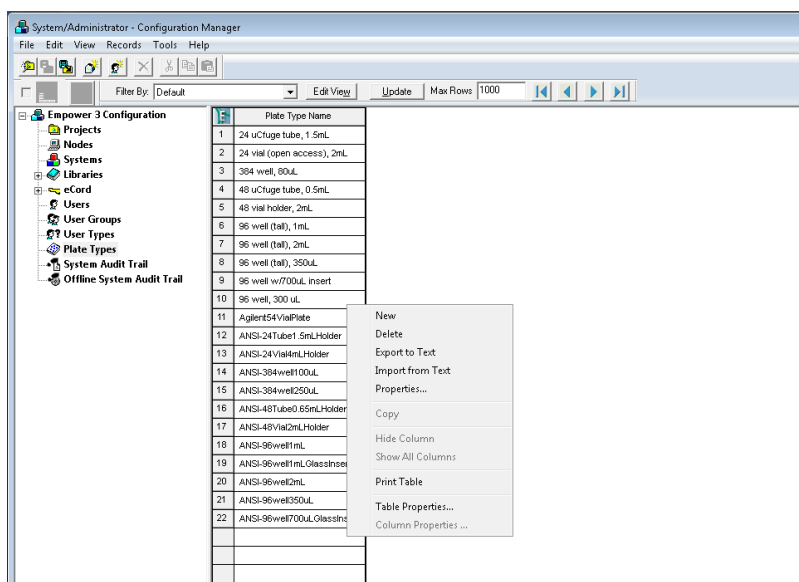


- 3 The **Import Plate Type From Text File** dialog appears. Click the **Browse** button and navigate to the **Agilent Plates for Import** folder as described in the next step.

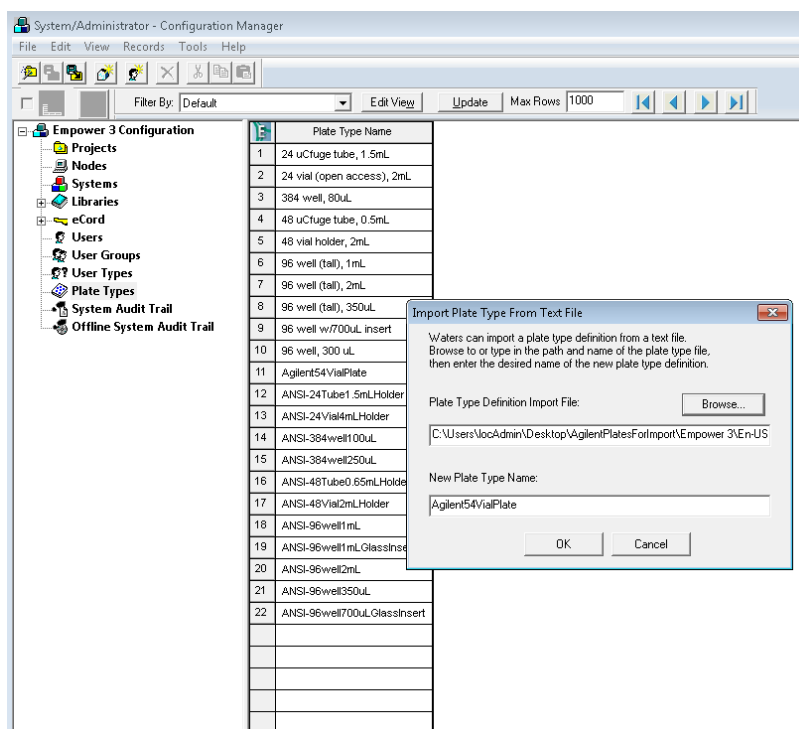

- 4 The location of the Agilent well plate import files are part of the Waters Empower Driver pack currently located on the Empower Driver Pack CD in folder **Agilent ICF/Agilent ICF Support v 2.1 Hotfix 1/AgilentPlatesForImport/Empower 3/En-US** and take it from here. Navigate to this folder using Explorer.

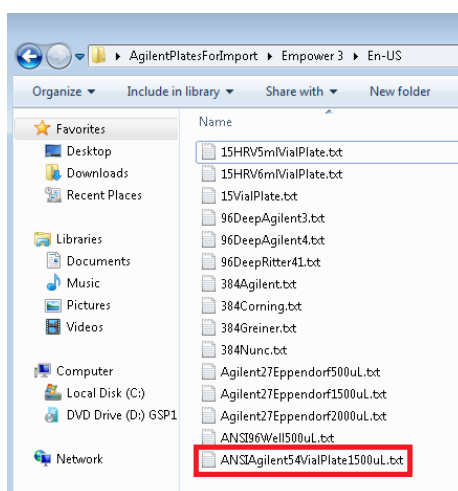

- 5 In this example, an Agilent 54-well plate will be selected. Select the **Agilent 54VialPlate** text file and click **OK** to import.

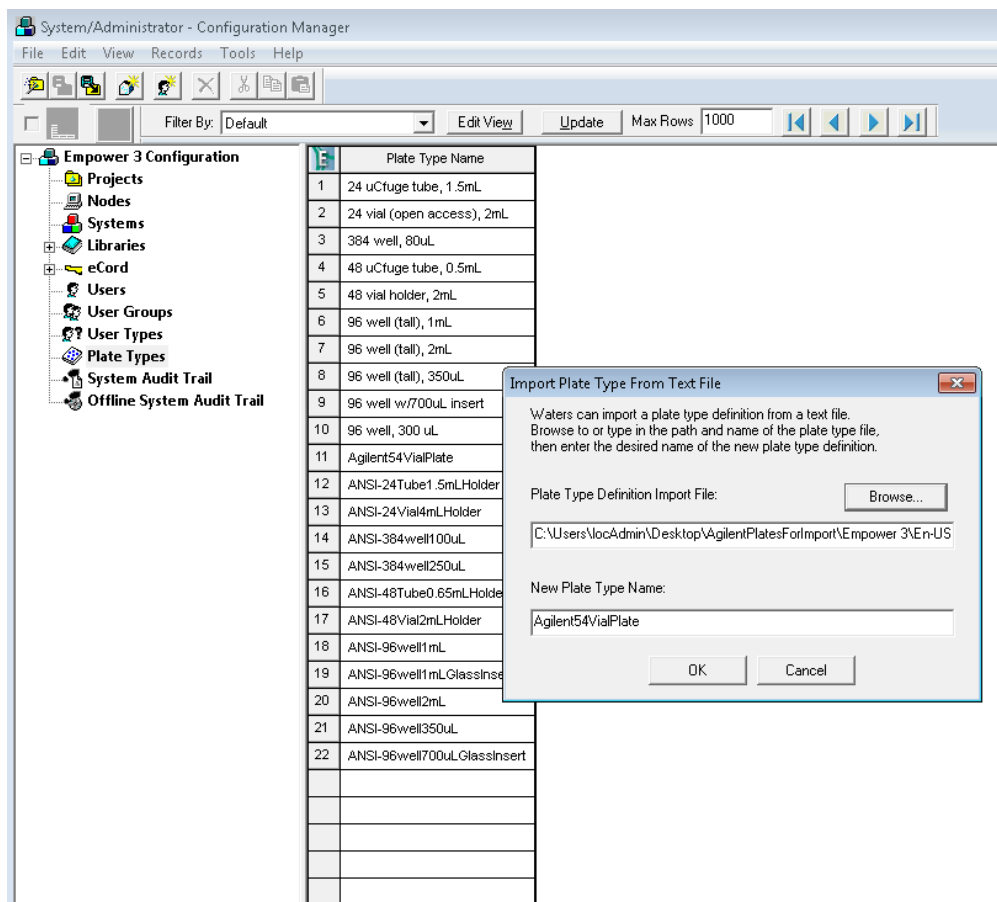

C - Manual Set Up of Well Plates and Example Configuration Layouts for Standard Agilent Well Plates

If the well/vial plate of interest is neither present in Empower nor part of the available well/vial plate definitions for import, the well/vial plate need to be manually defined within Empower.

As all dimensions are required to successfully define the well/plates you find the parameters for the some of the Agilent well/plates below. If you know the dimensions, proceed to the next step – manual set up.

The 3 screens shown above refer to Agilent's 54 vial plates.

The 3 screens shown above refer to Agilent's 96 well plates.

Appendices

C - Manual Set Up of Well Plates and Example Configuration Layouts for Standard Agilent Well Plates

The image displays three screenshots of the Agilent Well Plate Setup dialog box, showing the configuration for an Agilent 384 well plate. The dialog box has three tabs: Plate, Rows and Columns, and Referencing.

Plate Tab:

- Plate Type: Agilent 384 well plate
- Format: XY
- Plate Dimensions - mm:
 - X: 127.80
 - Y: 85.60
 - Height: 14.40
- Well Dimensions - mm:
 - Top Left Well Location:
 - X: 3.70
 - Y: 3.70
 - Well Size:
 - Diameter: 3.70
 - Depth: 10.20

Rows and Columns Tab:

- Row and Column Dimensions - mm:
 - Rows: 16
 - Columns: 24
 - Spacing: 4.50
- Row and Column Offsets - mm:
 - Row Offset:
 - None (selected)
 - Odd
 - Even
 - Offset: 0.00
 - Column Offset:
 - None (selected)
 - Odd
 - Even
 - Offset - mm: 0.00

Referencing Tab:

- Origin:
 - Top Left (selected)
 - Top Right
 - Bottom Left
 - Bottom Right
- Scheme:
 - Referencing: XY (selected), Sequential
 - Horizontal: ABC... (selected), 123...
 - Vertical: ABC... (selected), 123...
 - Sequential Continuous (unchecked)
 - Horizontal First Priority (unchecked)
- Terminology:
 - Plate: Plate
 - Well: well

The 3 screens shown above refer to Agilent 384 well plates.

Manual Set Up of Well/Vial Plates

- 1 To manually define plates go to the main **Edit** menu then click **Plates**.
- 2 In the upcoming dialog select **Create New Plate Type**.

Define Plates For Sample Set Method

☐ 2790 Layout Plate Sequencing Mode:

	Plate Type Name	Plate Layout Position
1	Agilent54	1
2	Agilent54	2
3	Agilent54	3
4	Agilent54	4
5	Agilent54	5
6	Agilent54	6

Visual representation of a 96-well plate layout (A1-A9, B1-B9, C1-C9, D1-D9, E1-E9, F1-F9).

Inject Standards

NOTE

The selected plates defined in the table above must be the same type and position as defined in the **Tray and Plate Configuration** dialog of the Multisampler graphical user interface (refer to step 8 on page 6) otherwise a mismatch will result when starting the analysis.

New Plate Type

Plate Rows and Columns Referencing

Row and Column Dimensions - mm

	Number	Spacing
Rows	8	9,00
Columns	12	9,00

Row and Column Offsets - mm

Row Offset: ☐ None ☐ Odd ☐ Even Offset: 0,00

Column Offset: ☐ None ☐ Odd ☐ Even Offset - mm: 0,00

D - Optional Multisampler Method Setup Parameters

The information provided below are additional/optional steps in setting up a Multisampler method.

- 1 This set of parameters can be used to reduce the amount of potential carryover in the system. Click **Injection Path Cleaning** and enter the corresponding values in the fields provided.

Instrument Method | Pretreatment Method | Auxiliary Channels | General | Instrument Configuration |

Binary Pump | Multisampler | Column Comp. | DAD |

Multisampler (G7167B)

Injection

Injection volume: 5.00 µL

Needle Wash

Standard Wash

Stoptime

☒ As Pump/No Limit
☐ 1.00 min

Posttime

☒ Off
☐ 1.00 min

Advanced

Injection Path Cleaning

Standard Wash


Mode: Flush Port
Time: 3 s
Location:
Repeat: 3

Multi-wash


Step	Solvent	Time [s]	Seat Back Flush	Needle Wash	Comment
1	S1	3	☒	☒	
2	S2	3	☒	☒	
3	S3	3	☒	☒	
Start Cond.	S1		☒	☒	

Done

- 2 Right mouse click on Multisampler **Instrument Status** and click **Control**. Select the parameters as shown in the example and click **OK**.

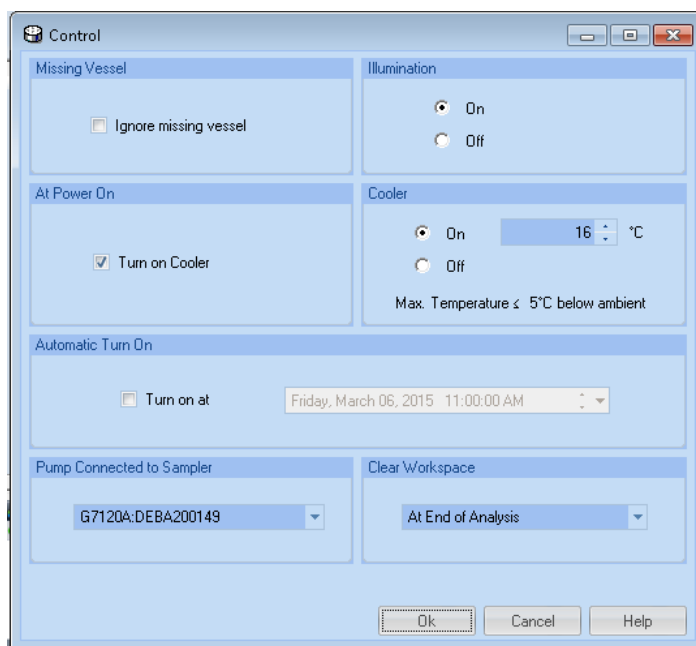


- 3 Right mouse click on Multisampler **Instrument Status** and click **Auto-clean**.

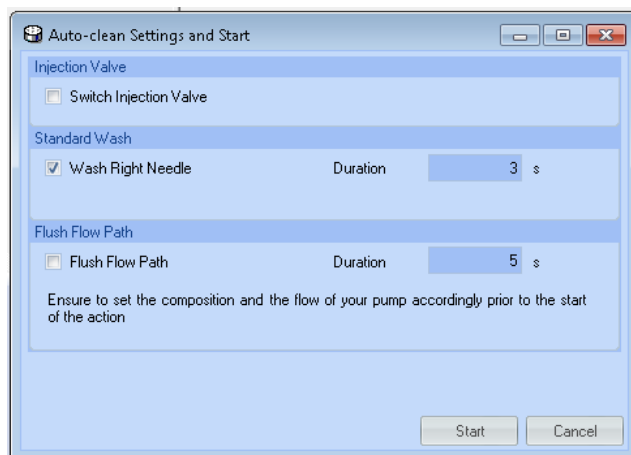

- 4 Select the **Duration** and click the **Start** button.

- 5 Return to the Multisampler **Instrument Status**, click on **Prime**, enter the Duration period and then click **Start**.

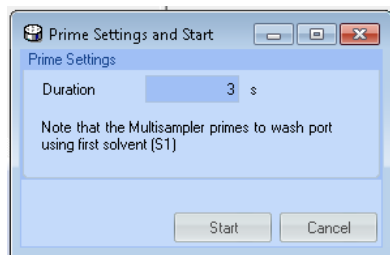

NOTE

The use of the reference vial array is currently not supported.

It is not possible to enter the reference vial position (1-5) for the sample set.

NOTE

For more information on the available functions, please consult the Agilent 1200 Infinity Series Multisamplers User Manual.



Part Number: G7167-90150

Edition: 10/2015
Printed in Germany

© Agilent Technologies, Inc 2015

Agilent Technologies, Inc
Hewlett-Packard-Strasse 8
76337 Waldbronn
Germany