


# Column Test Mix

On GraceSmart™ HPLC Column

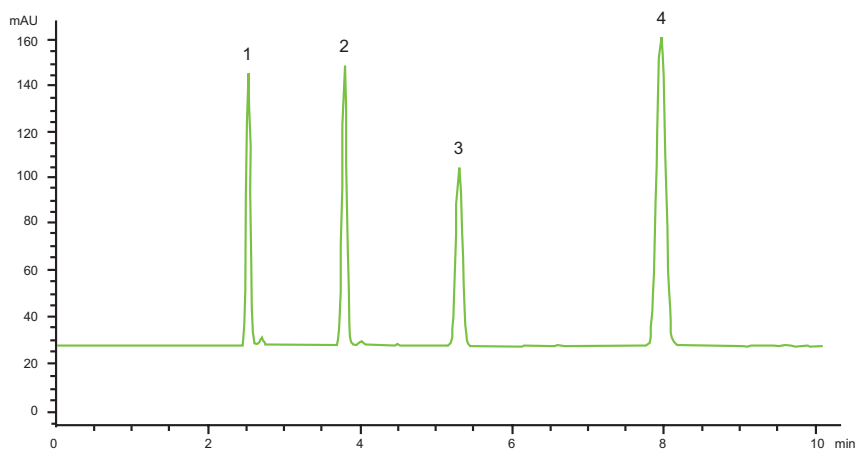

1. Uracil
2. Phenol
3. Diethy-ethyl toluamide
4. Toluene

	Tailing Factor	Theoretical plates	Plates/meter
Toluene	1.008	25847	103388

### Test condition

**Sample:** C of A. Test Mix

**Mobile Phase:** ACN:H<sub>2</sub>O 58:42

**Column:** GraceSmart™ C18, 120Å, 5µm, 4.6 x 250mm (Part #: 5138810)

**Flow:** 1ml/min

**Detection:** 254nm

**Column temp:** 30°C

**Injection Volume:** 5µl

GRACE®, GRACESMART™, and GRACE DAVISON DISCOVERY SCIENCES™ are trademarks and/or registered trademarks of W. R. Grace & Co.-Conn. The trademarks designated with the ® symbol are registered in the United States and/or other countries. This trademark list has been compiled using available published information as of the publication date of this brochure and may not accurately reflect current trademark ownership. Grace Davison Discovery Sciences is a product group of W. R. Grace & Co.-Conn. Alltech Associates, Inc. is a wholly owned subsidiary of W. R. Grace & Co.-Conn. © Copyright 2006 Alltech Associates, Inc. All rights reserved. The information presented herein is derived from our testing and experience. It is offered for your consideration and verification. Since operating conditions vary significantly, and are not under our control, we disclaim all warranties on the results that may be obtained from the use of our products. Grace reserves the right to change prices and/or specifications without prior notification. Printed in the USA.

11/20/2006, M178

### Grace Davison Discovery Sciences™ Regional Headquarters:

**In the Americas:**  
 2051 Waukegan Rd.  
 Deerfield, IL 60015  
 Tel: 847-948-8600  
 Email: [contact.alltech@grace.com](mailto:contact.alltech@grace.com)

**In Europe:**  
 Brandstraat 12  
 B-9160 Lokeren, Belgium  
 Tel: 09-340-65 65  
 Email: [info-lokeren.alltech@grace.com](mailto:info-lokeren.alltech@grace.com)

**In Asia:**  
 19th Floor, K.Wah Center  
 1010 Hua Hai Zhong Road  
 Shanghai 200031 P.R.C  
 Tel: 86-21-5467 4678  
 Email: [dsbiz.asia@grace.com](mailto:dsbiz.asia@grace.com)

**In Australia/New Zealand:**  
 30 Brookhollow Avenue  
 Baulkham Hills  
 New South Wales 2153, Australia  
 Tel: 1300 36 24 12  
 Email: [custcare-aust.discoverysciences@grace.com](mailto:custcare-aust.discoverysciences@grace.com)