

# Agilent MassHunter Optimizer

## *Automated MS Method Development Software*

### Quick Start Guide

What is MassHunter Optimizer?	2
Startup	3
User Interface	4
Menus and Toolbar	5
Compound Setup Tab	8
Precursor Ion Selection Tab	8f
Product Ion Selection Tab	8
Optimizer Setup Tab	7
Database Browser Window	8
Ion Breakdown Profile Dialog Box	9
Getting Started	10
Using Projects	12
Setting Optimization Parameters	14
Setting Up Compounds and Working with Databases	18
Setting Up Sequences	30
Optimization and Results	37

## Where to find more information

Use the online help for in-depth information not included in this *Quick Start Guide*.

Display online help by pressing the **F1** key to get more information about a window, tab, or dialog box.



**Agilent Technologies**

# What is MassHunter Optimizer?

**Automated MS Method Development Software** MassHunter Optimizer software lets you automatically optimize the data acquisition parameters for MRM mode (multiple-reaction monitoring) on a triple quad mass spectrometer instrument for each individual compound analyzed. Specifically, it automates the selection of the best precursor ions, the optimization of the fragmentor voltage for each precursor ion, selection of the best product ions, and optimization of collision energy values for each transition for a list of compounds you specify.

**Automated Optimization** Automation of the following optimization tasks results in a significant time saving over manual optimization procedures:

- Selection of optimum precursor ion via rules, such as M+H, M+Na, M+K, and M+NH<sub>4</sub>.
- Optimization of fragmentor voltage to maximize precursor ion intensity for up to 4 precursor ions (adducts and/or charge states) for each polarity in MS SIM
- Selection of multiple best product ions (MS/MS full scan)
- Optimization of collision energy to maximize product ion intensity (MS/MS full scan)
- Optionally perform fine fragmentor optimization for maximum sensitivity.

**Additional Features**

- Stores optimum precursor and product ions and parameters in a compound database.
- Imports optimization results into Agilent MassHunter Workstation triple quad Data Acquisition software.
- Lets you import compounds for optimization and update compound retention times from Agilent MassHunter Workstation triple quad Data Acquisition methods
- Supports sample introduction via manual infusion (syringe pump), automatic infusion (via loop injection at lower flows), and injection analysis with column (to maximize optimization yield).
- Data that is acquired during the optimization run can be displayed with Agilent MassHunter Qualitative Analysis software.

**NOTE**

You can optimize other source and ion funnel parameters using the "ion funnel and source optimization" program. For more information, see *online Help* and the *Familiarization Guide* for MassHunter Workstation Data Acquisition software.

## Startup

MassHunter Optimizer software is automatically installed with Agilent MassHunter Acquisition software. Two versions of MassHunter Optimizer software are installed on your system: one for small molecules and one for peptides.


Open MassHunter Optimizer software in *either* of the following ways:



- Double-click the **Optimizer** icon on the desktop, *or*
- Select **Programs> Agilent> MassHunter Workstation> Optimizer** from the Windows Start menu.

## To optimize peptides

A version of MassHunter Optimizer software is available for optimizing peptides. Open Optimizer for Peptides software in *either* of the following ways:

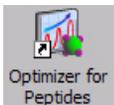

- Double-click the **Optimizer for Peptides** icon on the desktop, *or*
- Select **Programs> Agilent> MassHunter Workstation> Optimizer for Peptides** from the Windows Start menu.

### NOTE

Projects are specific to the version of Optimizer software that you use to create them. In other words, projects created with Optimizer for Peptides software *cannot* be opened in MassHunter Optimizer software and vice versa.

## To remove the software

- Click **Programs and Features** in Windows 7 Control Panel.
- Select **Agilent MassHunter Workstation Acquisition Software** in the list of programs, then click **Uninstall**.

Note that Optimizer can only be removed as part of Agilent MassHunter Workstation Acquisition software, not separately.

# User Interface

The MassHunter Optimizer software has the following UI elements that let you enter parameters, set options, view status, and perform functions:

- “Menus and Toolbar” on page 5
- “Status Bar” on page 7

### Tabs

- “Optimizer Setup Tab” on page 7
- “Precursor Ion Selection Tab” on page 8
- “Product Ion Selection Tab” on page 8
- “Compound Setup Tab” on page 8

### Other Windows and Dialog Boxes

- “Peptide Fragment Ion Predictor Dialog Box” on page 8
- “Database Browser Window” on page 8
- “Ion Breakdown Profile Dialog Box” on page 9

## Menus and Toolbar

The following menu items and toolbar buttons are available in MassHunter Optimizer software main window.

### File menu



**New Project** Clears the current settings and lets you set up a new project for use in MassHunter Optimizer software. See “To create a project” on page 12.



**Open Project** Lets you open a project. See “To open a project” on page 12.



**Save Project** Lets you save changes to the current project.



**Save As Project** Lets you save the current project under a new name.



**Save Compounds** Lets you save compound information to a database, for use with other projects. See “To save compounds to a database” on page 25.



**Delete Projects** Lets you select projects and remove them.

**Exit** Closes the MassHunter Optimizer software window.

### Edit menu

**Copy** Copies the selected rows, which can then be pasted into another position in the table.

**Cut** Cuts the selected rows from the table, which can then be pasted into another position in the table.

**Paste** Pastes the most recently copied or cut rows into the selected position in the table.

**Delete** Removes the selected rows from the table

**Select All** Selects all rows in the table.

**Fill** Copies the value in the top cell down to the rest of the cells of a highlighted block or to the entire column.



**Clear Results** Lets you clear all fragmentor values, all collision energy values, or all results, including methods and all precursor and product ions for all compounds in the selected row and below in the table.

### View menu

**Show/Hide Columns** Lets you select which columns of information to display in the table. At least one column in every row must be displayed.

**Expand/Collapse All Rows** Lets you display all rows in the table or collapse the table to show only Compound rows. This affects the currently selected row and all rows below it in the table.

### Import/Export menu



**Import from Database** Opens the Database Browser window, which lets you select compounds for optimization from the database. See “Database Browser Window” on page 8.



**Import from Excel** Lets you import compounds from an Excel file. See “To import compounds from an Excel file” on page 27.



**Import from Acquisition Methods** Lets you import all compounds from Acquisition methods into the current project.



**Export to Excel** Lets you export the current compound list for use in Excel. See “To export compound information to an Excel file” on page 27.

### Optimization menu



**Start Optimization** Starts the optimization process with the current parameters. The results are saved to the database and displayed in the Compound Setup tab. A report is automatically printed on completion. See “To start optimization” on page 38.



**Ion Breakdown Profile** Runs optimization with the current parameters and displays the results in the Ion Breakdown Profile dialog box when optimization is complete. Collision energy values can be selected from the graphs, instead of using those determined by optimization. See “To start optimization and display ion breakdown profiles” on page 39.



**Stop Optimization** Lets you terminate optimization before completion.

## Tools menu



**Update RT** Lets you update compound retention times from Data Acquisition methods. See “To update retention times from a MassHunter Acquisition method” on page 29.

*The following items only appear in Optimizer for Peptides software.*



**Extract Peptide** Opens the Peptide Fragment Ion Predictor dialog box, which lets you extract peptide sequences for optimization from a selected MassHunter Qualitative Analysis method. See “Peptide Fragment Ion Predictor Dialog Box” on page 8.



**Convert .csv to .psq** Lets you convert multiple peptide sequences with or without modification into multiple **.psq** files that can be imported into MassHunter Qualitative Analysis software. See “[To convert multiple peptide sequences for import into MassHunter Qualitative Analysis](#)” on page 36.

## Help menu

**Help Topics** Opens the online help window with the **Contents** tab displayed.

**About Optimizer** Displays version information for the software.

## Status Bar

The following items appear in the status bar at the bottom of the MassHunter Optimizer software main window: Project Name, Optimizer Status, Current Record progress bar, and All Records progress bar.

## Optimizer Setup Tab

Lets you set the acquisition parameters used for optimization, including sample introduction method, Acquisition methods, and ranges for optimization of fragmentor and collision energy voltages. See “To set Optimizer acquisition parameters” on page 14.

## Precursor Ion Selection Tab

Lets you select adduct species to include in the search for precursor ions. The default is **+H** for positive and **-H** for negative. See “To select precursor ions” on page 15.

## Product Ion Selection Tab

Lets you set criteria to use in the search for product ions. See “To select product ions” on page 16.

## Compound Setup Tab

Displays the list of compounds to optimize, and lets you enter precursor and product ion values, if known. It also displays the results of optimization, when you mark the **Show results summary** check box. Sequences rather than compounds are used for optimization in Optimizer for Peptides software.

## Peptide Fragment Ion Predictor Dialog Box

*Available in Optimizer for Peptides software only.*



This lets you select peptide sequences for optimization from a MassHunter Qualitative Analysis method. The selected sequences are added to the Compound Setup tab. After you click the **Extract Peptide** toolbar button shown at the left, you are prompted to select the method of interest. See “To import peptide sequences from a MassHunter Qualitative Analysis method” on page 32.

## Database Browser Window



This window appears when you click the **Import from Database** toolbar button. It lets you select compounds for optimization from the database. The selected compounds are added to the Compound Setup tab. See “To import compounds from a database” on page 22.

The Database Browser window also lets you open a different database, save the current database with a new name, and make the current database the default database.

## Ion Breakdown Profile Dialog Box



This dialog box automatically appears after optimization is complete when you click the **Ion Breakdown Profile** toolbar button. The product ion profiles determined by optimization are displayed and you can manually select collision energy values for each ion to replace the results of optimization. See “[To start optimization and display ion breakdown profiles](#)” and “[To manually set collision energy voltages](#)” on page 39.

## Getting Started

The following topics will help you get started using MassHunter Optimizer software.

- Using Projects**
  - “To create a project” on page 12
  - “To open a project” on page 12
  - “To edit a project” on page 12
  - “To save parameters to the current project” on page 13
  - “To save parameters to a new project” on page 13
- Setting Optimization Parameters**
  - “To set Optimizer acquisition parameters” on page 14
  - “To select data path and methods to use for optimization” on page 15
  - “To select precursor ions” on page 15
  - “To select product ions” on page 16
  - “To set Optimizer acquisition parameters for peptides” on page 16
- Setting Up Compounds and Working with Databases**
  - “To set up a compound list” on page 19
  - “To use known ions for individual compounds” on page 21
  - “To open a database and make it the default database” on page 22
  - “To import compounds from a database” on page 22
  - “To select transitions and triggers for Dynamic and Triggered MRM” on page 24
  - “To save compounds to a database” on page 25
  - “To filter compounds in Database Browser” on page 26
  - “To search compounds in Database Browser” on page 26
  - “To import compounds from an Excel file” on page 27
  - “To export compound information to an Excel file” on page 27
  - “To import compounds from a MassHunter Data Acquisition method” on page 27
  - “To update retention times from a MassHunter Acquisition method” on page 29
- Setting Up Sequences**
  - “To set up a sequence list for peptides” on page 30

**Optimization and  
Results**

- “To import peptide sequences from a MassHunter Qualitative Analysis method” on page 32
- “To convert multiple peptide sequences for import into MassHunter Qualitative Analysis” on page 36
- “To start optimization” on page 38
- “To review optimization results” on page 38
- “To start optimization and display ion breakdown profiles” on page 39
- “To manually set collision energy voltages” on page 39
- “To import optimization results to Acquisition for dMRM/tMRM” on page 41
- “To import optimization results to Acquisition for MRM time segments” on page 44

# Using Projects

## To create a project

Use this procedure to store all the values and settings used for optimization in a project.



- 1 Click the **New Project** button on the toolbar.

If current parameters have not already been saved, you are prompted to save the project. The current parameters are cleared.

- 2 Make the desired changes to parameters on the following tabs:

- Optimizer Setup tab
- Precursor and Product Ion Selection tabs
- Compound Setup tab



- 3 Click the **Save Project** or **Save As Project** button and enter a name for the project.

## To open a project

Use this procedure to use optimization values and settings from an existing project.



- 1 Click the **Open Project** button on the toolbar.
- 2 Click to select the project on the Select Project Name dialog box.
- 3 Click the **Open** button.

## To edit a project

- 1 Open the project to edit as described above.
- 2 Make the desired changes to parameters on the following tabs:
  - Optimizer Setup tab
  - Precursor and Product Ion Selection tabs
  - Compound Setup tab
- 3 Save the parameters as described below.

## To save parameters to the current project



- 1 Click the **Save Project** toolbar button.

## To save parameters to a new project



- 1 Click the **Save As Project** toolbar button.
- 2 Enter a name for the project and click the **Save** button.

## Setting Optimization Parameters

To set Optimizer acquisition parameters

# Setting Optimization Parameters

Use the following procedures to set up the optimization parameters for MassHunter Optimizer software.

“To set Optimizer acquisition parameters” *below*

“To select data path and methods to use for optimization” *below*

“To select precursor ions” on page 15

“To select product ions” on page 16

### Workflow for Peptides

“To set Optimizer acquisition parameters for peptides” on page 16

“To select data path and methods to use for optimization” *below*

## To set Optimizer acquisition parameters

- 1 Click the **Optimizer Setup** tab.
- 2 Set the desired **Sample introduction** method (**Injection**, **Automatic infusion**, or **Manual infusion**).
- 3 Set the Fragmentor ramp parameters as follows:  
*(Skip this step if your instrument is a model 6490 LC/MS. The fragmentor parameters are not displayed, since fragmentor voltage is not optimized for that instrument.)*
  - a Set the range for ramping the **Fragmentor** values.
  - b *(optional)* Mark the **Fragmentor Fine** check box and enter a **Step** size for Fine ramping. Fine optimization refines the coarse ramping values and provides better optimization but takes longer to run.
- 4 Set the range for ramping the **Collision Energy** values.
- 5 Continue to select data path and methods to use for optimization as described below.

## To select data path and methods to use for optimization

- 1 If not already displayed, click the **Optimizer Setup** tab.
- 2 Select a **Path for data files** to store the optimization run data. The data can be displayed later with Agilent MassHunter Qualitative Analysis software.
- 3 In the table on the right side of the screen, select the acquisition methods to use for optimization as follows. *These methods will be used to optimize all compounds, except for those that have specific methods selected on the Compound Setup tab.*
  - a Right-click in the table and select **Add Method** from the shortcut menu. This adds a row to the end of the table for a new method.
  - b Click the button on the right side of the **Acq Method** cell to open the Open Method dialog box.
  - c Select the method of interest and click **OK**. The **Polarity** and **Ion Source** will be filled in from the values set in the selected method.
  - d Check to make sure that the **Ion Source** from the method matches the physical configuration of your instrument.
  - e (*optional*) Repeat steps 3a - d to select additional methods for optimization.

**Tip** Additional options are available by right-clicking in the method table, such as **Delete Methods**, **Copy**, and **Fill**.

## To select precursor ions

Use this procedure to specify precursor ions to use to optimize the fragmentor voltage. The information entered here will be used for all compounds, except for those that have precursor ions specified on the Compound Setup tab.

- 1 Click the **Precursor Ion Selection** tab.
- 2 Select the **Positive** and **Negative** adduct ions to include.  
If necessary, add Positive or Negative adduct ions to the list by typing the value in the box and clicking the  button.
- 3 Set the search priority of the precursor ions in *one* of the following ways:
  - **Mark** the **Use most abundant precursor ion** option to use the most abundant precursor ion.

## Setting Optimization Parameters

### To select product ions

- *Clear* the **Use most abundant precursor ion** option and use the **Up** and **Down** arrow buttons to set the search order (ions at the top of the list are given more priority).
- 4 To exclude certain masses from consideration, click **Exclude masses** at the bottom of the screen. Enter the **m/z values** to exclude separated by commas and/or enter a **Minimum abundance** value in counts.

### To select product ions

Use this procedure to specify product ions to use to optimize the collision energy voltage. The information entered here will be used for all compounds, except for those that have product ions specified on the Compound Setup tab.

- 1 Click the **Product Ion Selection** tab.
- 2 Enter a **Low mass cut-off** value. Select whether to specify the value as **Mass (m/z)** or as a **% Precursor mass (m/z)** and enter the cut-off value.
- 3 To exclude certain masses from consideration, click the **Exclude masses** option at the bottom of the screen. Enter the **m/z values** to exclude separated by commas and/or enter a **Minimum abundance** value in counts.
- 4 If desired, you can also enter **Neutral Losses** to exclude, e.g. water (H<sub>2</sub>O). Enter a formula in the box and click the  button to add it to the list.

### To set Optimizer acquisition parameters for peptides

- 1 Click the **Optimizer Setup** tab in Optimizer for Peptides software.
- 2 Set the desired **Sample introduction** method (**Injection**, **Automatic infusion**, or **Manual infusion**).
- 3 Set the Fragmentor Ramp Settings parameters as follows. *Skip this step if your instrument is a model 6490 LC/MS, since fragmentor voltage is not optimized for that instrument.*
  - a In the Fragmentor area, mark the **Fixed Fragmentor** check box and enter the fragmentor voltage to use for optimization. This value will be

used for all compounds, unless you enter specific values for individual compounds on the Compound Setup tab.

- Tip** Using a fixed fragmentor voltage saves time by skipping the SIM run to determine the optimal fragmentor voltage. However, if multiple adducts are specified, then a SIM run is made to determine the precursor ion to use.
- b** Mark the **Overwrite existing** check box to use this value for all compounds, *even if* values are set for individual compounds on the Compound Setup tab.
  - 4** In the Collision Energy area, mark the **Start from predicted value** check box, which lets the Optimizer software set a collision energy value, using the formula shown. In the formula, **m/z** is the mass of the precursor ion.
  - 5** (optional) Mark the **Put instrument into standby mode at end** check box to put the instrument into standby mode at the end of the optimization run.
- See** “To select data path and methods to use for optimization” on page 15.

# Setting Up Compounds and Working with Databases

Use the following procedures to set up compounds for use in MassHunter Optimizer software.

## Setting up compounds

- “To set up a compound list” on page 19
- “To use known ions for individual compounds” on page 21

## Using compound databases

- “To open a database and make it the default database” on page 22
- “To import compounds from a database” on page 22
- “To select transitions and triggers for Dynamic and Triggered MRM” on page 24
- “To save compounds to a database” on page 25
- “To filter compounds in Database Browser” on page 26
- “To search compounds in Database Browser” on page 26

## Importing and exporting compound information from Excel

- “To import compounds from an Excel file” on page 27
- “To export compound information to an Excel file” on page 27

## Using compound information from MassHunter Acquisition methods

- “To import compounds from a MassHunter Data Acquisition method” on page 27
- “To update retention times from a MassHunter Acquisition method” on page 29

## To set up a compound list

Use this procedure to set up the list of compounds to optimize and enter precursor and product ion values, if known.

- 1 Click the **Compound Setup** tab.
- 2 (*optional*) Import compounds as described in “To import compounds from an Excel file” on page 27 or “To import compounds from a database” on page 22.
- 3 Clear the **Show results summary** check box above the table while you set up the compound list.
- 4 To add a compound:
  - a Right-click in the table and select **Add Compound** from the shortcut menu to add a row to the end of the table.
  - b Enter the **Compound Name** for the new compound.
  - c (*optional*) If you want to be able to find groups of compounds in the database such as sulfa drugs or pesticides, assign a group name in the **Groups** column. You can then specify the desired group names when you filter compounds to import from the database as described on page 22.
  - d Enter the **Formula** of the compound. The **Mass** will be automatically calculated from the formula. If you don't know the formula, then enter the monoisotopic mass in the **Mass** column.
  - e Enter the **Sample Position** for the new compound.
  - f (*optional*) Adjust the **Optimization dwell time** value to set a shorter or longer cycle time if desired; otherwise a default value is used.
  - g (*optional*) To speed up optimization, specify precursor and product ions for the compound as described in “To use known ions for individual compounds” below. If no methods or ions are specified here, then optimization for the compound uses the methods from the Optimizer Setup tab and information from the Precursor Ion Selection and Product Ion Selection tabs to generate the ions.
  - h Repeat the steps above to add more compounds to the table.

### NOTE

Compounds are global to all projects. Compound information such as name, group, formula, and mass in one project will be reflected in the entire database.

## Setting Up Compounds and Working with Databases

### To set up a compound list

- Tip** To cut or copy and paste compounds (rows) in the table:
- a** Select the rows of interest in the table.
  - b** Right-click on the selected rows and select **Cut** or **Copy** from the shortcut menu.
  - c** Right-click on the new location for the cut or copied rows and select **Paste** from the shortcut menu.
- 5** Mark the **Select** columns for the compounds (rows) to use for optimization.
  - 6** Save the compound list as described in “To save compounds to a database” on page 25 or “To save parameters to the current project” on page 13.
  - 7** (*optional*) Export the compound list as described in “To export compound information to an Excel file” on page 27.

## To use known ions for individual compounds

Use this procedure to optionally specify methods, precursor ions, and product ions for individual compounds on the Compound Setup tab. Using known ions speeds up the optimization process.

- 1 Add compounds to the Compound List as described in “[To set up a compound list](#)” on page 19.
- 2 To add methods for the current compound:
  - a Right-click in the **Compound** row of interest and select **Add Method** from the shortcut menu. This adds a Method row below the selected compound.
  - b Click the down arrow button on the right end of the **Acq Method** cell and select an acquisition method to use for the selected compound. If you don't see the method of interest, add it to the Method Table on the Optimizer Setup tab as described in “[To set Optimizer acquisition parameters](#)” on page 14.
  - c Repeat the above steps to add more methods for the current compound.
- 3 To specify a known precursor ion for the current compound:
  - a Right-click in the **Method** row for the compound of interest and select **Add Precursor Ion** from the shortcut menu. This adds a Precursor Ion row for the current compound.
  - b Enter a value for the **Precursor Ion**.
- 4 To specify product ions for the current compound:
  - a Right-click in the **Precursor Ion** row for the compound of interest and select **Add Product Ion** from the shortcut menu. This adds a Product Ion row for the current compound.
  - b Enter a value for the **Product Ion**.
  - c Repeat the above steps to add the product ions for the current compound and precursor ion.
- 5 Repeat steps 2 - 4 to specify methods and ions for other compounds in the Compound Table.

## To open a database and make it the default database

Use this procedure when you want to use a database other than the default optimizer database.



**1** If the Database Browser window is not already open, click the **Import from Database** button on the toolbar to open it.



**2** Click the **Open Database** button on the Database Browser toolbar.

**3** Select the database of interest and click **OK**.

**4** You are asked if you want to make this database the default database.

### NOTE

If you are importing additional compounds into an existing project, click **No**.

Otherwise, unsaved changes made to the current project will be lost.

All saved compound information in the previous project will still exist in the previous default database, but it won't be available for use with the new default database.

### Tip



If you decide later to make this database the default database, click the **Set Default Database** button on the Database Browser toolbar.

## To import compounds from a database

Use this procedure to import compounds that have previously been saved as described in “[To save compounds to a database](#)” on page 25.



**1** Click the **Import from Database** toolbar button to open the Database Browser.

**2** If not already open, open the database of interest as described on [page 22](#).

**3** Select compounds for optimization from the database.

- If desired, filter compounds as described on [page 26](#).
- If desired, search compounds as described on [page 26](#).
- To display all compounds in the database, mark the **Show All Records** check box.

- 4 When the compounds from step 3 appear in the Compound table in the lower half of the Database Browser window, select the transition ions to import in *one* of the following ways:
  - Individually mark the check boxes in the first column of the table for the transition ions you want to import.
  - Click the **Select top ranked transitions** option, then enter how many top transitions to import for each compound, and click the **Select Transitions** button. You can rank the top transitions by either **Abundance** or **Response Factor**.
  - To select the primary transitions, click the **Primary Transitions** option, then click the **Select Transitions** button.
  - To select the secondary transitions, click the **Secondary Transitions** option, then click the **Select Transitions** button.

**NOTE**

To set up the primary and trigger transition ions to import, see “To select transitions and triggers for Dynamic and Triggered MRM” on page 24.

- 5 Do one of the following:
  - If the compounds in the table are the only ones you want to import at this time, click **Import**. This closes the Database Browser window and adds the compounds to the Compound Setup tab.
  - If you want to do additional search and filter operations on the current database or a different database, click **Add to Import List**.
- 6 Do one of the following:
  - To find more compounds in the *same* database, repeat steps 3 - 5.
  - To select compounds from a *different* database, repeat steps 2 - 5.
- 7 Click the **Import List** tab to review the list of compounds.
  - Right-click and select **Remove** from the shortcut menu to remove any compounds you don't want to import.
- 8 Click the **Import** button.

The imported compounds will appear on the Compound Setup tab.

## To select transitions and triggers for Dynamic and Triggered MRM

- 1 If not already open, open the database of interest as described on [page 22](#).
- 2 In the Database Browser window, display compounds for optimization from the database.
  - If desired, filter compounds as described on [page 26](#).
  - If desired, search compounds as described on [page 26](#).
  - To display *all* compounds in the database, mark the **Show All Records** check box.
- 3 Deselect any unwanted compounds by clearing the check box next to the **Compound Name** in the first column of the Compound table.

### Triggered MRM method

- 4 Select transitions for a *triggered MRM method* as follows:
  - a In the **Select top ranked transitions** field, enter the maximum number of transitions per compound that you want to show in the import list. You can have two primary transitions, plus additional secondary transitions (*see step c below*).
  - b Click the **Select Transitions** button to display the transitions in the Compound table.
  - c In the **Set primary and trigger flags** area, specify the number of **Primary** transitions to use. The rest are assigned as Secondaries. Typically, two primaries are used.
  - d Click **Set Primaries and Trigger**. The most abundant transition is automatically assigned as the Trigger.
  - e Review the compound list and edit as needed.
    - A very abundant low  $m/z$  ion may be unsuitable as a Primary transition.
    - You can manually select a different transition for a compound as the Primary transition, and define a low  $m/z$  ion as a Secondary.
    - You may want to select two Primaries as triggers for a compound.

### Dynamic MRM method

- 5 Select transitions for a *dynamic MRM method* as follows:
  - a Click **Primary transitions**.
  - b Click the **Select Transitions** button to display the transitions in the Compound table.
    - At this point, all transitions are displayed as primaries.

- Typically, the transitions are ranked by abundance.
- c** Review the compound list and edit as needed.
- Deselect any unwanted transitions.
- 6** When the list of transitions is set, click the **Add to Import List** button below the Compound table.
- 7** Do one of the following:
- To find more compounds in the *same* database, repeat steps 2- 6.
  - To select compounds from a *different* database, repeat steps 1- 6.
- 8** Click the **Import List** tab to review the list of compounds.
- Right-click and select **Remove** from the shortcut menu to remove any compounds you don't want to import.
- 9** Click the **Import** button.
- The imported compounds will appear on the Compound Setup tab.

## To save compounds to a database

Do one of the following to save compounds to a database:



- To save compounds to the current database:
  - Click the **Save Compounds** button on the main toolbar, or



- Click the **Save Database** button on the Database Browser toolbar.



- To save compounds to a *new* database, click **Save As Database** button on the Database Browser toolbar, enter a new name, and click the **Save** button.

You can import the compounds into other projects at a later time as described in “[To import compounds from a database](#)” on page 22.

## To filter compounds in Database Browser

- 1 If not already open, open the database of interest as described on [page 22](#).
- 2 Enter the Filter criteria in the upper left area of the Database Browser Search/Filter tab. Note that filter criteria are unavailable if **Show All Records** is selected.
  - Mark the **Enable Filters** check box.
  - Mark the check boxes for the filter criteria to use, such as **Optimized Compounds, Date, Group Name, Project Name, Polarity, Model, and Method**. Enter values for each selected criteria.
- 3 Click the **Search/Filter** button.

The compounds that pass the filter are displayed in the Compound table in the lower part of the screen.

## To search compounds in Database Browser

- 1 If not already open, open the database of interest as described on [page 22](#).
- 2 Enter the **Search Text** in the upper right area of the Database Browser Search/Filter tab, such as **Compound Name** and **Formula**. This can be either a partial or complete entry as described in step 4.

### NOTE

The values you enter here for **Product** or **Precursor** ions are treated as *numeric values* rather than text. Matches are returned whose values fall between 0.7 amu below the entered mass and 0.3 amu above the specified mass.

- 3 Mark the columns to search or click the top entry in the list (Select Columns) to search *all* columns.
- 4 Do one of the following:
  - To find compounds that match only a partial name or formula, *clear* the **Match entire word for each string** check box.
  - To find only compounds that exactly match the text you entered in step 2, *mark* the **Match entire word for each string** check box.
- 5 To ignore all filters for this search, *clear* the **Enable filters** check box.

**NOTE**

If you want to display *all items* in the database, mark the **Show All Records** check box. This automatically disables all the search and filter options.

- 6 Click the **Search/Filter** button.
  - The compounds that meet the search criteria are displayed in the Compound Table in the lower part of the screen.
  - Search text that was *not found* in the search is shown in red.

## To import compounds from an Excel file



- 1 Click the **Import from Excel** button on the main toolbar.
- 2 Select the **.xls**, **.xlsx**, or **.csv** file to import, then click **Open**. All compounds from the file will be imported and listed on the Compound Setup tab.  
The Select Method for Linking Compounds dialog box may be displayed to assign methods to the compounds.

## To export compound information to an Excel file



- 1 Click the **Export to Excel** button on the main toolbar.
- 2 Enter a name to export the compound list in **.xls** format for use in Excel, click **Save**.

## To import compounds from a MassHunter Data Acquisition method

- 1 Create MassHunter Acquisition methods for the compounds you want to import.



- 2 Click the **Import from Acquisition Methods** toolbar button in MassHunter optimizer software.
- 3 Select the MassHunter Acquisition methods from step 1 that contain the compounds of interest, then click the **Open** button.

## Setting Up Compounds and Working with Databases

To import compounds from a MassHunter Data Acquisition method

All compounds in the data acquisition methods will be added to the current project and can be seen on the Compound Setup tab.

## To update retention times from a MassHunter Acquisition method

- 1 Create MassHunter Acquisition methods for the compounds of interest.
- 2 Be sure that compounds whose retention times you want to update are part of a project. Only retention times for compounds that are part of a project will be updated. Retention times for compounds that are not in a project will not be updated, even if those compounds are in the database.



- 3 Click the **Update RTs** button on the MassHunter optimizer software main toolbar.
- 4 Select the MassHunter Acquisition methods from step 1 that contain the compounds of interest, then click the **Open** button.
- 5 When the Update Retention Time Options dialog box appears:
  - a Select whether to **Update in current project** or to **Search database for eligible projects**.
  - b To update retention time windows, mark the **Update RT Windows** check box. If marked, the retention time windows in the acquisition method will be used to update the retention time windows in the database.
  - c If transition ions are found that do not match existing transition ions, you can add them by marking the **Add new transitions found** check box. Otherwise these transitions will be ignored.
  - d Click **OK**.
- 6 Several scenarios may result:

**Projects are found that contain the same methods and transition ions** In this case, select the project to update, and click **OK**.

*or*

**No project matches for update** In this case, you are prompted to select whether to add the transition ions to the **Current Project** or a **New Project**. Click the desired option, then click **OK**.

All transition ions of the same compound will be updated with the first retention time in the acquisition method. The updated compound information can be seen on the Compound Setup tab.

- 7 Be sure to save the project to save the updated retention times.

# Setting Up Sequences

Use the following procedures to set up peptide sequences for use in Optimizer for Peptides software.

- “To set up a sequence list for peptides” below
- “To import peptide sequences from a MassHunter Qualitative Analysis method” on page 32
- “To convert multiple peptide sequences for import into MassHunter Qualitative Analysis” on page 36

## To set up a sequence list for peptides

Use this procedure to set up the list of peptides to optimize in Optimizer for Peptides software.

- 1 Click the **Compound Setup** tab.
- 2 Clear the **Show results summary** check box above the table while you set up the sequence list.
- 3 Import sequences as described in “To import peptide sequences from a MassHunter Qualitative Analysis method” on page 32.
- 4 (*optional*) If you want to be able to find groups of sequences in the database, assign a group name in the **Groups** column. You can then specify the desired group name when you filter sequences to import from the database.
- 5 Leave the **Formula** column blank. The **Mass** is imported from the method and displayed in the table automatically.
- 6 Enter the **Sample Position** for the sequence.
- 7 Mark the **Select** columns for the sequences (rows) that you want to use for optimization.
- 8 Mark the **Perform multi-compound run** check box. This lets you perform collision energy optimization when all the fragment ions are known, a typical case in peptides using ion prediction.
- 9 (*optional*) Adjust the **Optimization dwell time** value to set a shorter or longer cycle time if desired.

- 10** Click the **Fill Values** button to calculate the Cycle Time parameters and the peak width necessary to obtain ten data points required for optimization when using Fixed Fragmentor or CE value formula on the Optimizer Setup tab. See “[To set Optimizer acquisition parameters for peptides](#)” on page 16.
- 11** Save the sequence list as described in “[To save compounds to a database](#)” on page 25 or “[To save parameters to the current project](#)” on page 13.
- 12** (*optional*) Export the sequence list as described in “[To export compound information to an Excel file](#)” on page 27.


# To import peptide sequences from a MassHunter Qualitative Analysis method

*Applies to Optimizer for Peptides software only.*

- 1 Create a MassHunter Qualitative Analysis method that contains the peptide sequences you want to import.
  - If desired, create the **.psq** files for MassHunter Qualitative Analysis as described in “[To convert multiple peptide sequences for import into MassHunter Qualitative Analysis](#)” on page 36.
- 2 If you need to add custom modifications:
  - a Select **Open Chemical Data Dictionary Editor** from the Tools menu.
  - b Click the + button on the **Modifications** tab to add a custom modification.
  - c Type in a name for a new modification, then enter its **Chemistry** values.

### Example

Assume you wish to add a custom modification “heavy Proline”, where the carbons and nitrogen in Proline are replaced with 13-C and 15-N. A Proline residue has five Carbons and one Nitrogen. Assuming complete labeling, specify **Losses** of C5 N1 and **Gains** of Cx5 Nx1 for a Delta mass of 6.01381. Mark the check box for Proline and clear all others to specify this custom modification for Proline only.

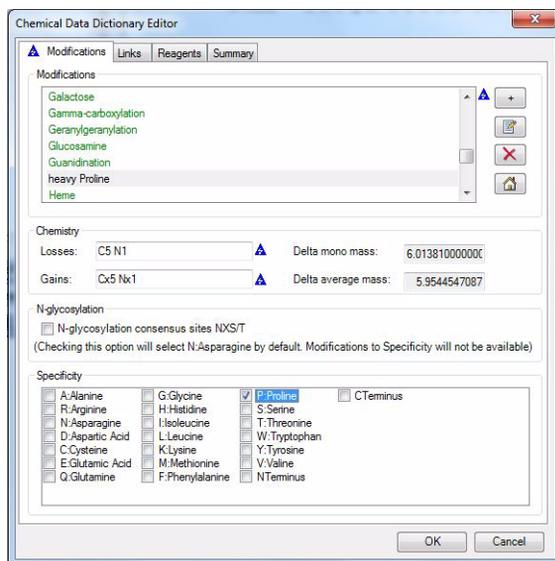

- d Click **OK** to close the Chemical Data Dictionary Editor.
- 3 Apply the custom modification to the sequences of interest as follows:
  - a Open the **Sequence Editor** window if it is not already open, and select the sequence of interest.
  - b Using the example above, select the Proline amino acid in the sequence, then apply the custom modification created in Step 2.
  - c Save the method, which contains the modified sequences.
- 4 Be sure to observe the following caveats if you modified the Chemical Data Dictionary as described in Step 2 above:
  - Close the MassHunter Qualitative Analysis software window. Otherwise, the custom dictionary will not be modified.
  - If Optimizer was open while the custom modifications were added, restart Optimizer to make sure that it reads the new custom dictionary.
  - If you plan to use this method on a different computer, you need to add the same custom modifications to the Chemical Data Dictionary on that system. Otherwise Optimizer for Peptides won't recognize the custom modifications and an error will result.

## Setting Up Sequences

To import peptide sequences from a MassHunter Qualitative Analysis method

5 Import the sequences into MassHunter Optimizer for Peptides software as follows:

a Review the charge states in the **Precursor Ion Selection** tab. Make sure that the desired charge states are listed, and add them if they are not.

### Tip

For peptide analysis, charge state precursors 2+ and 3+ are usually preferred, and singly-charged precursors are ignored.



b Click **Extract Peptide** on the toolbar in Optimizer for Peptides software.

c Select the MassHunter Qualitative Analysis method from step 1 that contains the peptide sequences of interest, then click the **Open** button. All sequences from the selected method are displayed in the “**Peptide Fragment Ion Predictor Dialog Box**” on page 8.

d Right-click in the table and select **Expand/Collapse All** from the shortcut menu to see the Precursor Ions and Charge states for each sequence.

e Mark the sequences that you want to use for optimization.

### Tip

To make it easier to edit long transitions lists, mark the **Show selected ions** option above the table to display only the selected ions.

- Within each peptide sequence, mark the precursor charge state to use for that peptide and *deselect* all charge states that you don't want to use. For example, you may want to examine only the 2+ charge state precursor because it produces abundant product ions, but the 3+ charge state precursor does not.
- If you decide to examine multiple charge states for a given peptide, then select all of the charge states of interest.
- Select the product ions of interest that result from each peptide precursor. Use the following check boxes in combination with selecting individual product ions in the table to make your selections.
  - Show selected ions
  - Show only y-ions
  - Show only b-ions
  - Show ions > than m/z value
  - Show ions > than % of precursor

f Click the **Add** button. The selected peptide sequences from the method are imported and listed on the “**Compound Setup Tab**” on page 8.

- 6 Review the list of precursors and product ions.
  - a Right-click in the table and select **Expand/Collapse All Rows** from the shortcut menu to show all precursors and product ions.
  - b To edit the list, right-click on a product ion row, and select Delete Product Ions or Add Product Ion from the shortcut menu.
- 7 If you want to use only a single charge state for each peptide to optimize, then edit the charge states on the **Precursor Ion Selection** tab to make sure that only one charge state is selected for each peptide. That is, remove the charge state selection from all but one charge state. For example, clear charge state selections except for Charge State 2.

**NOTE**

If more than one charge state is active, then during optimization, the software tries to determine which charge state has the highest abundance.

---

- 8 (*optional*) Save the compounds.

## Setting Up Sequences

To convert multiple peptide sequences for import into MassHunter Qualitative Analysis

# To convert multiple peptide sequences for import into MassHunter Qualitative Analysis


*Applies to Optimizer for Peptides software only.*

Use this procedure to convert multiple peptide sequences with or without modification into multiple **.psq** files for import into MassHunter Qualitative Analysis software.

- 1 Create a comma separated (**.csv**) file that has one sequence per line, with the following format:

**NameString, SequenceString, Modification1String, Location1String, Modification2String, Location2String**

### Example

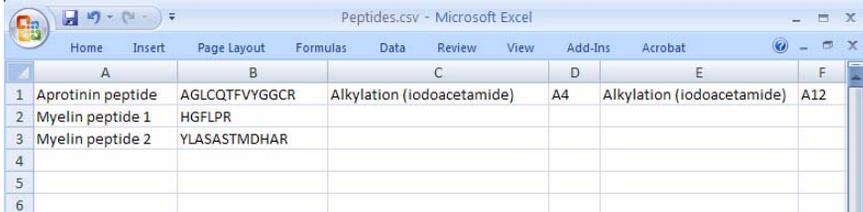

	A	B	C	D	E	F
1	Aprotinin peptide	AGLCQTFVYGGCR	Alkylation (iodoacetamide)	A4	Alkylation (iodoacetamide)	A12
2	Myelin peptide 1	HGFLPR				
3	Myelin peptide 2	YLASASTMDHAR				
4						
5						
6						



- 2 Click the **Convert .csv to .psq** toolbar button in Optimizer for Peptides software.

- 3 Select the **.csv** file created in step 1, and click the **Open** button.

The file is converted into multiple **.psq** files, one for each sequence, which are saved in a folder named **OutputPSQFiles** in the same folder that contains the original **.csv** file.

- 4 Import the **.psq** files into MassHunter Qualitative Analysis software and save them in a Data Analysis method.
- 5 Import the sequences for use in Optimizer for Peptides software as described in “[To import peptide sequences from a MassHunter Qualitative Analysis method](#)” on page 32.

## Optimization and Results

Use the following procedures to run the optimization procedure, evaluate the results, and import them into MassHunter Acquisition software.

“To start optimization” on page 38

“To review optimization results” on page 38

“To start optimization and display ion breakdown profiles” on page 39

“To manually set collision energy voltages” on page 39

“To import optimization results to Acquisition for dMRM/tMRM” on page 41

“To import optimization results to Acquisition for MRM time segments” on page 44

## To start optimization

- 1 Set the optimization parameters as described on [page 14](#).
- 2 Set up the compound list as described in “[To set up a compound list](#)” on [page 19](#) or “[To set up a sequence list for peptides](#)” on [page 30](#).
- 3 Start the optimization process in *either* of the following ways:



- Click the **Start Optimization** button on the toolbar.



- Click the **Ion Breakdown Profile** button on the toolbar if you want to display the product ion profiles determined by optimization before they are saved to the database. For more information, see “[To start optimization and display ion breakdown profiles](#)” on [page 39](#).



- 4 If necessary, you can stop the optimization process by clicking the **Stop Optimization** button on the toolbar.
- 5 Review results as described below.
- 6 (*optional*) Use Agilent MassHunter Qualitative Analysis software to look at the data acquired during the optimization run.

## To review optimization results

- 1 Run the optimization procedure as described above.
- 2 Click the **Compound Setup** tab.
- 3 Mark the **Show results summary** check box above the table.
- 4 Review the following values for each transition in the Compound table:

**Fragmentor**—The optimum voltage to apply to the exit end of the capillary for the compound, determined by MassHunter Optimizer software. (*Does not apply to instrument model 6490 LC/MS.*)

**Collision Energy**—The optimum voltage to apply to the ions in the collision cell for the compound, determined by MassHunter Optimizer software.
- 5 Review the printed optimization report.
- 6 See “[To import optimization results to Acquisition for MRM time segments](#)” on [page 44](#), to use optimization values for data acquisition.

## To start optimization and display ion breakdown profiles

Use this procedure to display the product ion profiles determined by optimization before they are saved to the database and to manually select collision energy values for each ion if desired.

1 Set the optimization parameters as described on [page 14](#).

2 Set up the compound list as described on [page 19](#).



3 Click the **Ion Breakdown Profile** toolbar button to start optimization.

4 When optimization is complete, the Ion Breakdown Profile dialog box is displayed. Review the graphs, which plot collision energy steps on the x-axis and abundance values on the y-axis for each product ion. The values shown in the table are the collision energy values determined by optimization. See an example of Ion Breakdown Profile results in [Figure 1](#) on the next page.

5 If desired, you can override the collision energy values as described below. Otherwise, click **Save & Exit** to close the Ion Breakdown Profile dialog box.

## To manually set collision energy voltages

1 Set optimization parameters and start optimization as described above.

2 Click on the Ion Breakdown Profile graph to select the desired collision energy value. See an example Ion Breakdown Profile graph in [Figure 1](#) on the next page.

3 Apply the selected value to either **All ions** or only the **Selected ion** by clicking the appropriate option for **Apply manual selection to**.

4 If desired, you can reload the collision energy values set by optimization at any time by clicking the **Load Optimizer Defaults** button. Doing this overwrites the manual selections you made above.

5 Click the **Save & Exit** button to close the Ion Breakdown Profile dialog box.

## Optimization and Results

To manually set collision energy voltages

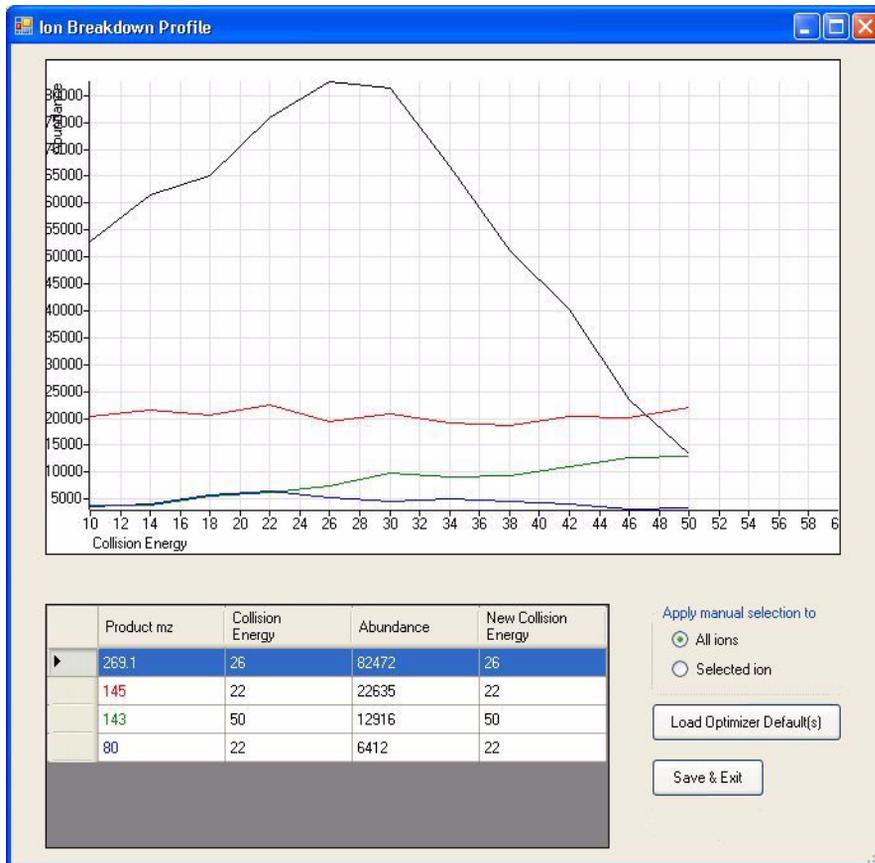


Figure 1 Ion Breakdown Profile Results

## To import optimization results to Acquisition for dMRM/tMRM

- 1 Display the **Method pane** in MassHunter Workstation Data Acquisition software.
- 2 Create or open the method of interest.
- 3 Click the **MS QQQ** tab in the Method Editor pane.
- 4 Click the **Acquisition** tab on the right side of the screen.

Time segments							
#	Start Time	Scan Type	Div Valve	Delta EMV (+)	Delta EMV (-)	Stored	
▶ 1	0	Dynamic MRM	To MS	200	0	<input checked="" type="checkbox"/>	

- 5 Set **Scan Type** to **Dynamic MRM** as shown above.

- 6 For triggered MRM:
  - a Mark the **Triggered** check box in the lower right area of the Acquisition tab.



- b Notice that additional columns are now available in the Scan segments table, including **Primary**, **Trigger**, **Trigger Window**, **Trigger MRM Threshold**, and **Trigger Entrance Delay**.
- 7 Use the expected retention time and RT window parameters to set the timetable for monitoring transitions. Usually only a single Time segment is used for Dynamic MRM and Triggered MRM. Note that the Cycle time remains constant.
    - As transitions are added to the list in real time, the Dwell time is reduced for each transition that is active to keep the Cycle time constant.
    - The reverse happens for transitions whose time has elapsed, and dwell times are *increased* for the remaining active transitions.
    - For a single compound on the list, one transition serves as the quantitation response. Other transitions for the same compound are used as qualifiers.

## Optimization and Results

### To import optimization results to Acquisition for dMRM/tMRM

- 8** Right-click in the **Acquisition** tab and select **Import from Optimizer** from the shortcut menu.
- 9** When the Database Browser window appears:
  - a** If desired, filter the compounds as described on [page 26](#).
  - b** If desired, search compounds as described on [page 26](#).
  - c** Mark the transition ions of interest from the Compound table in the lower half of the Database Browser window in *one* of the following ways:
    - Manually select the transition ions you want to import for each compound, *or*
    - Enter the top number of transitions to import for each compound, then click the **Select** button. You can select the top transitions based on either Abundance or Response Factor. Deselect any transitions that you don't want to use by clearing their check box.
- 10** If the transition ions selected in the table are the only ones you want to import at this time, click **Import**. This closes the Data Browser window and adds the transition ions into the current time segment.
  - Otherwise, to do additional search and filter operations on the current database or a different database, click **Add to Import List**.
- 11** To select additional compounds:
  - a** To select compounds from a different database, open the database of interest as described on [page 22](#).

#### NOTE

When asked if you want to make this database the default database, click **No**. Otherwise, unsaved changes made to the current project will be lost. All saved compound information in the previous project will still exist in the previous default database, but won't be available for use with the new default database.

- b** Repeat steps 10a - c to select the transition ions of interest.
  - c** Click **Add to Import List**.
- 12** Click the **Import List** tab to review the list of transition ions. To remove any transition ions you don't want to import, right-click and select **Remove** from the shortcut menu.
- 13** Click the **Import** button. The selected transitions are imported into the current time segment.

- 14 Review the timings to make sure that not too many transitions are being monitored concurrently:
  - a Right-click in the Scan segments table and select **Edit dMRM method** from the shortcut menu.
  - b When the Dynamic MRM Viewer appears, review how many transitions are active at any point in time during the run.

**NOTE**

When too many transitions are active at the same time, the dwell time per transition may become too short (1-2 msec per transition or less). Under these conditions, the signal/noise ratio is low, and the quantitation accuracy and precision may suffer.

- c If necessary, change the retention time windows or delete transitions to increase the average dwell time.
  - d You can also use the **Split Method** feature in the **Options and Parameters** pane of the Dynamic MRM Viewer to divide the transitions from the current method across two or more methods.
    - Select the criteria for splitting the method from one of the following: **Minimum dwell time**, **Maximum concurrent MRM**, or **Number of new methods**.
    - Note that all transitions for the same compound are assigned to the same method.
- 15 (*optional*) Save the method.

## To import optimization results to Acquisition for MRM time segments

- 1 Display the Method pane in MassHunter Workstation Data Acquisition software.
- 2 Create or open the method of interest.
- 3 Click the **MS QQQ** tab in the Method Editor pane.
- 4 Click the **Acquisition** tab on the right side of the screen.
- 5 Select the time segment of interest or create a new one.

Time segments						
#	Start Time ▲	Scan Type	Polarity	Div Valve	Delta EMV	Stored
1	0	MRM	Positive	To MS	200	<input checked="" type="checkbox"/>

- 6 Set **Scan Type** to **MRM**.
- 7 Right-click in the **Acquisition** tab and select **Import from Optimizer** from the shortcut menu.
- 8 When the Database Browser window appears:
  - a If desired, filter the compounds as described on [page 26](#).
  - b If desired, search compounds as described on [page 26](#).
  - c Mark the transition ions of interest from the Compound table in the lower half of the Database Browser window in *one* of the following ways:
    - Manually select the transition ions you want to import for each compound, *or*
    - Enter the top number of transitions to import for each compound and click the **Select** button. You can select the top transitions based on either Abundance or Response Factor. Deselect any transitions that you don't want to use by clearing their check box.
- 9 If the transition ions selected in the table are the only ones you want to import at this time, click **Import**. This closes the Data Browser window and adds the transition ions into the current time segment.
  - Otherwise, to do additional search and filter operations on the current database or a different database, click **Add to Import List**.

**10** To select additional compounds:

- a** To select compounds from a different database, open the database of interest as described on [page 22](#).

**NOTE**

When asked if you want to make this database the default database, click **No**. Otherwise, unsaved changes made to the current project will be lost. All saved compound information in the previous project will still exist in the previous default database, but won't be available for use with the new default database.

---

- Repeat steps 8a - c to select the transition ions of interest.
- Click **Add to Import List**.
- b** Click the **Import List** tab to review the list of transition ions. To remove any transition ions you don't want to import, right-click and select **Remove** from the shortcut menu.
- c** Click the **Import** button. The selected transitions are imported into the current time segment.

**11** Repeat step 5 through step 10 for each time segment of interest.

**12** (*optional*) Save the method.

[www.agilent.com](http://www.agilent.com)

## In this guide

This Quick Start Guide describes how to use MassHunter Optimizer Automated MS Method Development software.

This guide is valid for the B.06.00 revision or higher of the Agilent MassHunter Optimizer Automated MS Method Development Software, until superseded.

Windows® is a registered trademark of Microsoft Corporation in the US and/or other countries.

© Agilent Technologies, Inc. 2012

Printed in USA  
Revision A, November 2012



G3793-90007



**Agilent Technologies**