



Refrigerants

Application Note

Environmental

Authors

Agilent Technologies, Inc.

Introduction

Fast GC analysis of halogenated refrigerant mixtures is achieved in less than 50 seconds using an Agilent PoraPLOT U column and Agilent 490 Micro GC.



Agilent Technologies


Conditions

Technique : Micro GC
Column : Agilent PoraPLOT U, 0.25 mm x 10 m fused silica
Temperature : 119 °C
Carrier Gas : Helium
Pressure program : 100 kPa, 5 sec, → 250 kPa, 200 kPa/min
Injector : chip injector heated, 100 °C
Injection time : 5 ms
Detector : chip TCD, Sensitivity: med
Concentration Range : % level

Courtesy : Mario Voglino
Stuart Wallman,
A-GAS (UK) Ltd

Peak identification

1. air
2. CFC 32
3. CFC 143A
4. CFC 125
5. CFC 134A
6. CFC 22

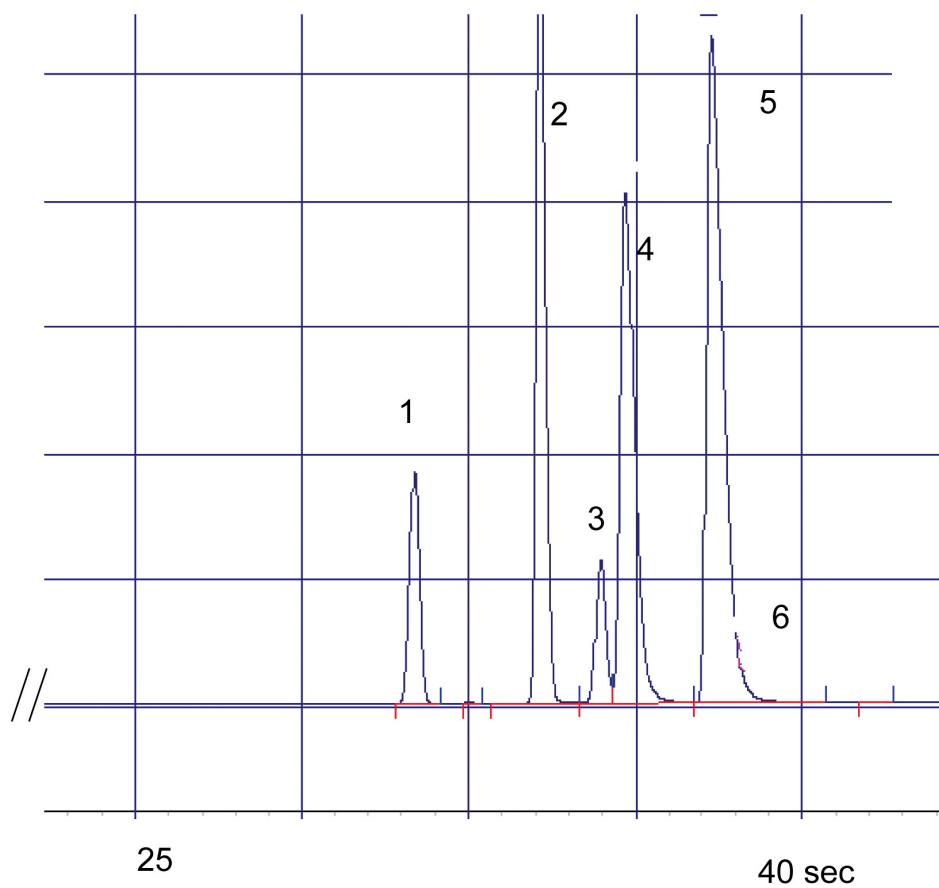

www.agilent.com/chem

This information is subject to change without notice.

© Agilent Technologies, Inc. 2011

Printed in the USA

31 October, 2011

First published prior to 11 May, 2010

A01957



Agilent Technologies