
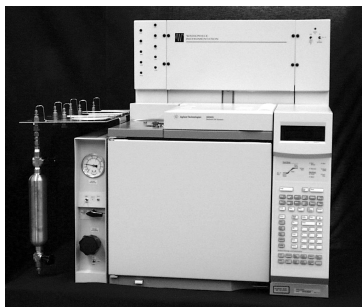


Application 383D-00

Agilent Refinery Gas Analyzer

Technical Overview

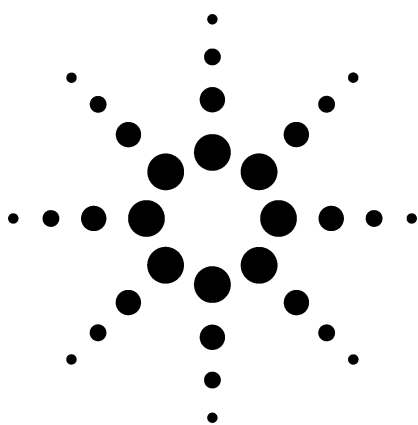



Application Highlights

- A Flame Ionization Detector (FID) is used to detect the C1 through C7 paraffins and olefins to a lower detection limit of 20 ppm, except for trace peaks eluting on the tail of a major component.
- A Thermal Conductivity Detector (TCD) is used to detect hydrogen in a nitrogen carrier to a lower detection limit of 100 ppm.
- A second TCD is used to detect carbon dioxide, ethane, ethylene, acetylene, hydrogen sulfide, oxygen/argon composite, nitrogen, methane, and carbon monoxide to a lower detection limit of 200 ppm except for carbon monoxide (400 ppm), and hydrogen sulfide (500 ppm).
- Analysis time is approximately 15 minutes.



Optional Configurations

- Refinery gas analysis with trace sulfurs by FPD or SCD
- Additional boiling point column for the analysis of heavy hydrocarbons (C1–C30)
- Standard analysis with the addition of trace CO by methanizer
- Custom analyzer for performing ASTM D2163, ASTM D2712, ISO 7941, and ASTM D1945
- High temperature injection for heavy fractions
- High temperature reactor effluent with percent level water
- TCD/TCD/MSD for the analysis of reactor effluent gases
- Liquid sample valves for the injection of pressurized liquid samples.

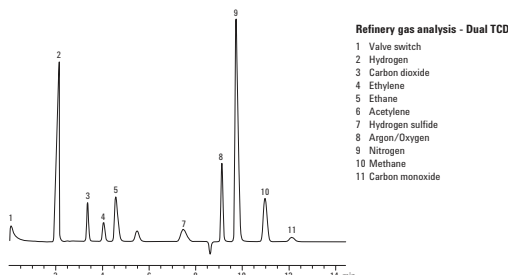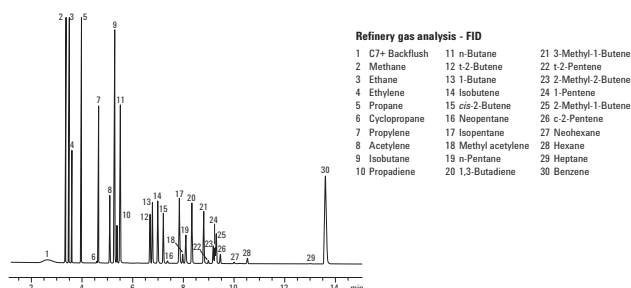

For More Information


For more information on our products and services, visit our Web site at www.agilent.com/chem.



Agilent Technologies




WASSON·ECE
INSTRUMENTATION

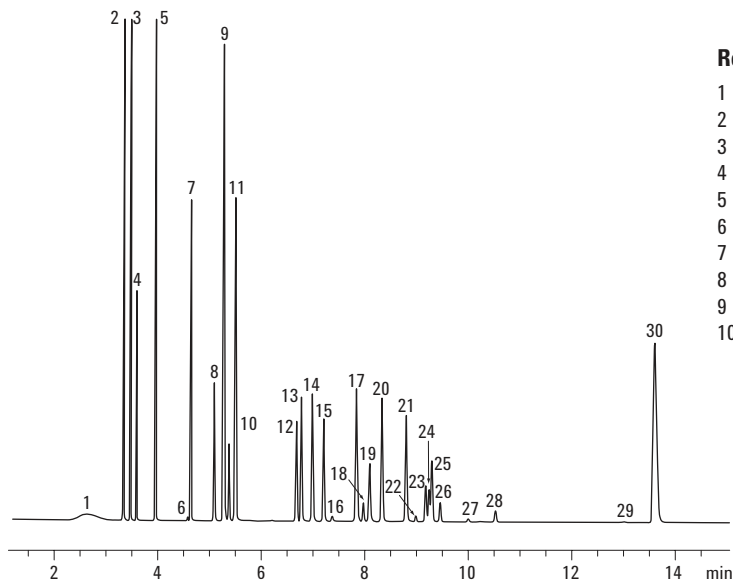

Refinery gas analysis - FID

1 C7+ Backflush	11 n-Butane	21 3-Methyl-1-Butene
2 Methane	12 t-2-Butene	22 t-2-Pentene
3 Ethane	13 1-Butane	23 2-Methyl-2-Butene
4 Ethylene	14 Isobutene	24 1-Pentene
5 Propane	15 cis-2-Butene	25 2-Methyl-1-Butene
6 Cyclopropane	16 Neopentane	26 c-2-Pentene
7 Propylene	17 Isopentane	27 Neohexane
8 Acetylene	18 Methyl acetylene	28 Hexane
9 Isobutane	19 n-Pentane	29 Heptane
10 Propadiene	20 1,3-Butadiene	30 Benzene

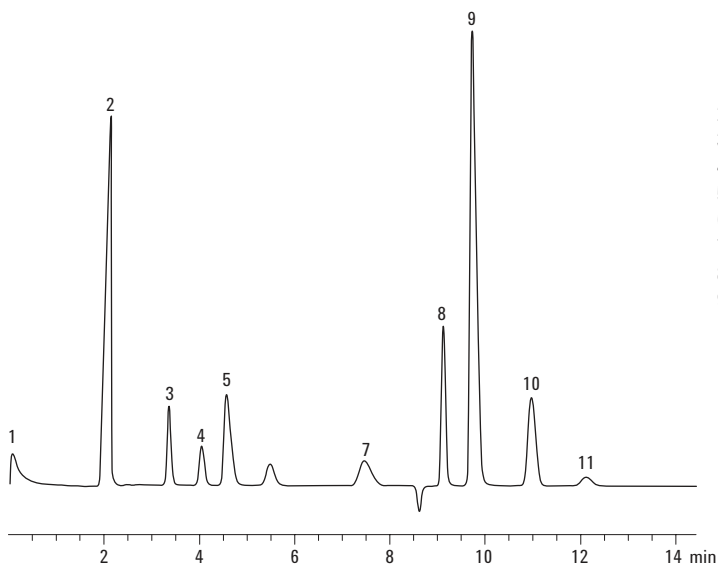

Refinery gas analysis - Dual TCD

1 Valve switch
2 Hydrogen
3 Carbon dioxide
4 Ethylene
5 Ethane
6 Acetylene
7 Hydrogen sulfide
8 Argon/Oxygen
9 Nitrogen
10 Methane
11 Carbon monoxide

FID and TCD output from the Agilent Refinery Gas Analyzer.

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice.

© Agilent Technologies, Inc. 2002

Printed in the USA
December 9, 2002
5988-6738EN