Agilent 8860 Gas Chromatograph

Safety Information

Agilent Technologies

English

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Safety Information

Dear Customer,

Thank you for purchasing a product from Agilent Technologies. Before you start to use your product, please read the corresponding documentation provided on the DVD/USB flash drive. It provides an extensive collection of user manuals, safety information and technical reference for individual instruments.

General Safety

The following general safety precautions must be observed during all phases of operation of this instrument. Failure to comply with these precautions or with specific warnings or operating instructions in the user manuals violates safety standards of design, manufacture, and intended use of the instrument. Agilent Technologies assumes no liability for the customer's failure to comply with these requirements. User manuals are provided with your instrument on this DVD/USB flash drive and/or in printed form. Manuals may also be available on the Web. Go to www.agilent.com and type in your product number in the Search field at the top of the page.

General

Do not use this product in any manner not specified by the manufacturer. The protective features of this product may be impaired if it is used in a manner not specified in the operation instructions.

Before Applying Power

Verify that all safety precautions are taken according to the user manual of the individual instrument. Verify that the voltage range and frequency of your power distribution matches to the power specification of the individual instrument. Never use cables other than the ones supplied by Agilent Technologies to ensure proper functionality and compliance with safety or EMC regulations. Make all connections to the unit before applying power. Note the instrument's external markings described under "Safety Symbols" in the Appendix of the corresponding user manual.

Ground the Instrument

If your product is provided with a grounding type power plug, the instrument chassis and cover must be connected to an electrical ground to minimize shock hazard. The ground pin must be firmly connected to an electrical ground (safety ground) terminal at the power

outlet. Any interruption of the protective (grounding) conductor or disconnection of the protective earth terminal will cause a potential shock hazard that could result in personal injury.

Do Not Operate in an Explosive Atmosphere

Do not operate the instrument in the presence of flammable gases or fumes.

Do Not Remove the Instrument Cover

Only Agilent authorized personnel are allowed to remove instrument covers. Always disconnect the power cables and any external circuits before removing the instrument cover.

Do Not Modify the Instrument

Do not install substitute parts or perform any unauthorized modification to the product. Return the product to an Agilent Sales and Service Office for service and repair to ensure that safety features are maintained.

In Case of Damage

Instruments that appear damaged or defective should be made inoperative and secured against unintended operation until they can be repaired by qualified service personnel.

Solvents

Toxic and hazardous solvents and flammable liquids.

The handling of solvents and reagents can hold health risks.

When working with solvents observe appropriate safety procedures (for example, goggles, safety gloves and protective clothing) as described in the material safety data sheet supplied by the solvent vendor, especially when toxic or hazardous solvents and flammable liquids are used.

Do not use solvents with an auto-ignition temperature below 200° C.

Fume Hood

During normal operation of the GC with many detectors and inlets, some of the carrier gas and sample vents outside the instrument through the split vent, septum purge vent, and detector exhaust. If any sample components are toxic or noxious, or if hydrogen is used as the carrier gas, these exhausts must be vented to a fume hood. Place the GC in the hood for proper ventilation.

Safety Notices

CAUTION

A CAUTION notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a CAUTION notice until the indicated conditions are fully understood and met.

WARNING

A WARNING notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a WARNING notice until the indicated conditions are fully understood and met.

Safety Symbols

The apparatus is marked with this symbol when the user should refer to the instruction manual in order to protect risk of harm to the operator and to protect the apparatus against damage.

Confirms that a manufactured product complies with all applicable European Community directives. The European Declaration of Conformity is available at: <http://regulations.corporate.agilent.com/DoC/search.htm>

Manufacturing date

Power symbol indicates On/Off. The apparatus is not completely disconnected from the main supply when the power switch is in the Off position.

To prevent personal injury, two person lift recommended.

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